



LPG valve block installation: pump must be switched off

LPG block variants

The AZURA® Pump P2.1L and the BlueShadow Pump 80 P can be equipped with a low-pressure gradient (LPG) valve block to deliver binary or ternary LPG gradients. The following variants of the LPG valve block are available:

Pump	Gradient type	Article no.
AZURA® Pump P 2.1L	Binary	AZZ00AA
	Ternary	AZZ00AB
Blueshadow 80P	Binary	AZZ00AA-1
	Ternary	AZZ00AB-1

Pre-installation procedure

NOTICE

Component defect

If the LPG valve block is installed while the pump is switched on, severe damage to components such as the pump's circuit board may occur.

→ Make sure that the pump is switched off before installation.

Procedure

1. Ensure that the pump is switched off and pull the power plug.
2. Remove the yellow protection cap (see Fig. 1).
3. Follow the steps given in the pump's instructions:
 - [V6840](#) for AZURA Pump P2.1L
 - [V7680](#) for BlueShadow Pump 80P

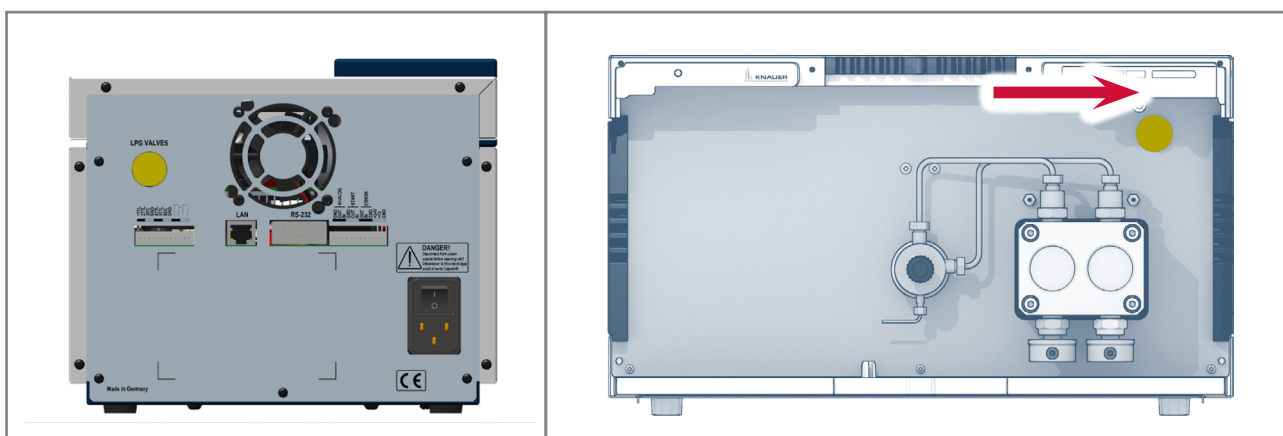


Fig. 1 LPG valve block protection cap on Blueshadow 80 P (left) and AZURA® Pump P 2.1L (right)

More information

For more information, contact support@knauer.net or visit www.knauer.net.