Produktinformation

AZURA Organizer

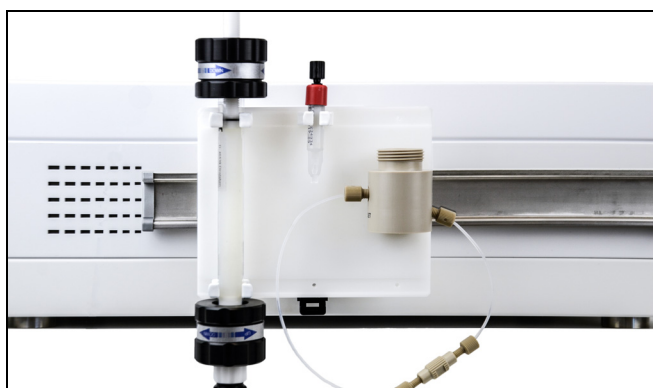
Abb. 1 AZURA Organizer

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AZURA Organizer

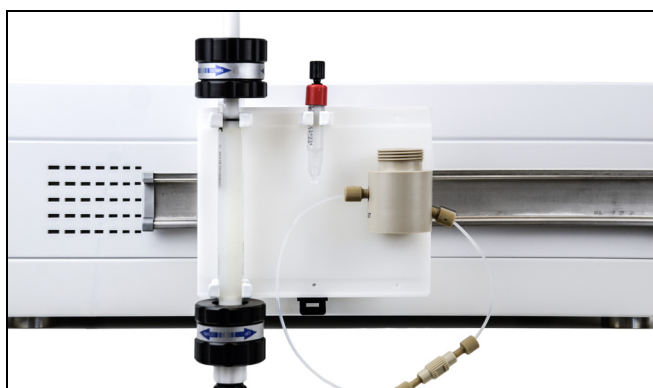
Abb. 1 AZURA Organizer

Mount different columns with diameters between 5 mm to 26 mm, Falcon tubes, back pressure regulators or a pH electrode onto AZURA Organizer. AZURA Organizer is attached to AZURA devices with L footprint by means of an accessory AZURA Click. Attach with clamps in various sizes that are part of the delivery.

12 mm clamp:

- for columns up to 13 mm outer diameter (example: His Trap 1 ml, ECO 10 mm OD, ECO Plus 10 mm OD)

16 mm clamp:

- for 15 ml Falcon tubes
- for columns up to 16 mm outer diameter (example: GSTrap 5 ml)
- for back pressure regulator

25 mm clamp:

- for 50 ml Falcon tubes
- for columns up to 26 mm outer diameter (example: HiLoad 16/60 Superdex)

Attaching the AZURA Organizer onto AZURA Click

Note: In case you are mounting columns to the AZURA Organizer, pay attention that there is enough space between the organizer and the bottom surface.

Prerequisites

- AZURA Click has been attached to the AZURA device.

Process

1. Hook the AZURA Organizer into the upper rail of the railing.
2. Press the AZURA Organizer against the lower rail of the railing.

Removing the AZURA Organizer

Process

1. Pull the handle of the support downward.
2. Remove the organizer upward.

Mounting the Clamps of the AZURA Organizer

Note: We recommend to use to organizers for SEC columns.

Note: Use two clamps to mount longer columns.

Note: Use the Allen head screws supplied with the delivery.

Tools

- Allen wrench, size 2.5 mm

Process

1. Pull approx. 1 cm of silicone tubing over both sides of the clamp.
2. Hold the clamp against the upper bore hole of the AZURA Organizer.
3. Using the wrench, fasten the screw from behind.
4. Adjust the clamp opening with the screw.

Mounting a pH Electrode onto the AZURA Organizer

Tools

- Allen wrench, size 2.5 mm

Process

1. Hold the pH electrode against the middle, diagonally arranged bore holes of the AZURA Organizer.
2. Using the Allen wrench, fasten the pH electrode from behind.

Repeat Orders

Name	Order number
AZURA Click	A70089
AZURA Organizer	A70085
12 mm clamp	A70085-1
16 mm clamp	A70085-2
25 mm clamp	A70085-3