Azura

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Table of Contents

Note: For your own safety, **read** the manual and **always** observe the warnings and safety information on the device and in the manual!

Safety	17
Symbols and Signs	18
Disassembling the Housing	19
Device Overview	20
Removing the Side Panels	21
Removing the Hood	23
Replacing the Deuterium Lamp	24
Removing the Lamp	24
Installing the Lamp	26
Repeat Orders for Lamp Components	28
Certificate for KNAUER Service Training	29

Safety

Note

It is absolutely necessary to retain this service supplement and to have it ready if service is required!

Professional group

This service supplement is exclusively addressed to persons who fulfill the following conditions:

1. Professional qualification as electrician or comparable vocational training has been successfully completed.
2. A written certificate confirms the successful participation in the KNAUER service training, which includes the lamp replacement in this device (see this supplement).

If you do not belong to this or a comparable professional group and do not fulfill both requirements, under no circumstances may you perform the work described in this service supplement. In this case, please contact your superior.

NOTE! If a person is to perform service tasks on the device and does not fulfill any of the professional group conditions, the contractual liability exclusion applies and KNAUER can not be held liable.

Symbols and Signs

Symbol	Meaning
Warning signs in the manual	
	High-voltage hazard
	Danger caused by hot components
Warning signs on the device	
	Electrostatic-discharge hazard

Disassembling the Housing

WARNING!

Burn hazard from hot lamp!
Switch off the lamp meanwhile the device remains switched on for further 15 minutes, otherwise the lamp can not cool down.
Afterwards, switch off the device and pull the power plug.

DANGER!

High voltages pose a life-threatening risk.
Before opening the device, turn off the device and pull the power plug.

Disassembling the housing is necessary when replacing the lamp.

Device Overview

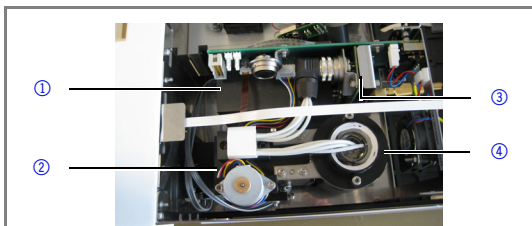
Abb. 1 Interior view

Legend

- ① D2 power supply
- ② Optical bench
- ③ Connection board with lamp plug
- ④ Lamp socket

Removing the Side Panels

Prerequisite

- Device is switched off.
- Power plug is pulled.

Tools

Allen screwdriver, size 2

Process and Figure

1. Open the locking device ①. The locking device can only be opened with the power plug being pulled.

Fig. 2 Lock for side panel

Process and Figure

- Loosen 2 screws ① for each side panel from the rear panel.
- Pull out the side panel towards the back.

Fig. 3 Screws side panel

Next steps

Remove the hood.

Removing the Hood

Prerequisite

- Device is switched off.
- Power plug is pulled.
- Side panels are removed.

Tools

Allen screwdriver, size 2

Process and Figure

1. Loosen 3 screws ① from the rear panel.
2. Pull out the hood towards the back.

Fig. 4 Screws hood

Result

The device has been opened.

Replacing the Deuterium Lamp

The recommended operating time for D2 lamps is 2000 hours. Replace the lamp if it malfunctions or its intensity is low.

Removing the Lamp

Prerequisite

- Device is switched off.
- Power plug is pulled.
- Side panels and hood are removed.
- Lamp has cooled down.

Tools

Allen screwdriver, size 2.5


WARNING!


Burn hazard from hot lamp!
Switch off the lamp meanwhile the device remains switched on for further 15 minutes, otherwise the lamp can not cool down.
Afterwards, switch off the device and pull the power plug.

Process and Figure

1. To remove the lamp connector, turn the locking ring ① .
2. Remove the connector.

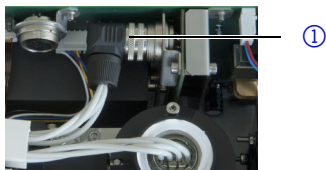**Fig. 5** Lamp connector

3. Loosen both screws ② .
4. Remove the lamp socket ③ with lamp from the lamp pod.

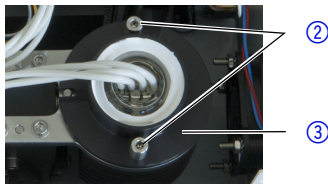**Fig. 6** Screws on the lamp socket

Installing the Lamp

Note

Do not touch the glass body with the hands. Deposits can falsify measuring results and can damage the lamp.

Prerequisite

- Housing is removed.
- Power plug is pulled.

Tools

Allen screwdriver, size 2.5

Process and Figure

1. Hold the lamp at the lamp socket and guide the glass body into the lamp pod ①.

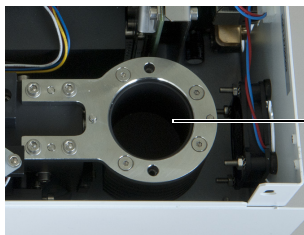
Fig. 7 Lamp pod

Process and Figure

2. Using the screwdriver, tighten the screws ② of the lamp socket ③ .

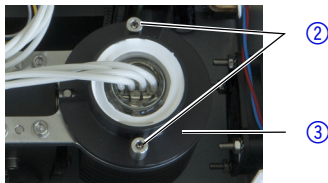
Fig. 8 Lamp socket

3. Plug in the connector and turn the locking ring ④ until the connector is secured.

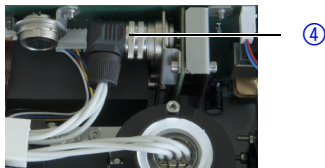
Fig. 9 Lamp connector

DANGER!

High voltages pose a life-threatening risk.

Before opening the device, turn off the device and pull the power plug.

Next steps

1. Close the housing.
2. Plug in the power plug.
3. Switch on the device.
4. Check the functionality of the lamp.

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Fax +49 30 8015010

E-Mail support@knauer.net

Certificate for KNAUER Service Training

Lamps shall only be replaced by persons, who participated in the certified service training for lamp replacement. The instructor hands out the certificate. For questions, contact our Technical Support:

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© Wissenschaftliche Gerätebau
Dr. Ing. Herbert Knauer GmbH
All rights reserved.
The information in this document is subject
to change without prior notice. Translation
of the original German edition of this sup-
plement, version 1.0.
2013-01-16
Printed in Germany on environmentally
friendly paper from sustainable forests.

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® AZURA are registered trademarks of