



Detector ECD 2.1 Instructions







Note: For your own safety, read the instructions and observe the warnings and safety information on the device and in the instructions. Keep the instructions for future reference.



Note: In case you require this instruction in another language, please submit your request including the corresponding document number via e-mail or fax to KNAUER.

Customer Support:	Phone :	+49 30 809727-111 (9-17h Central European Time)
	Fax:	+49 30 8015010
	E-Mail:	<u>support@knauer.net</u>
	Languages:	German, English
Publisher:	KNAUER Wiss	enschaftliche Geräte GmbH
	Hegauer Weg	138
	14163 Berlin	
	Germany	
	Phone:	+49 30 809727-0
	Fax:	+49 30 8015010
	Internet:	www.knauer.net
	E-Mail:	<u>info@knauer.net</u>
Version information:	Document nu	mber: V6685
	Version numb	ber: 1.0
	Release date:	2019/10/23
	Original edition	on
	without prior	on in this document is subject to change notice. For the latest version of the instructions, ite: www.knauer.net/library
Sustainability:	•	d versions of our instructions, we use environmentally r from sustainable forests.
Copyright:		nt contains confidential information and may not be repro- t written consent of KNAUER Wissenschaftliche Geräte
	© KNAUER W All rights rese	issenschaftliche Geräte GmbH 2019 rved.
		egistered trademark of

KNAUER Wissenschaftliche Geräte GmbH.

Table of contents

1.	General inf	ormation
	1.1	Intended use
	1.2	WEEE directive
	1.3	Warning symbols
	1.4	Safety instructions
2.	Introductio	n7
	2.1	Instrument description7
3.	Installation	
	3.1	Site preparation requirements9
	3.2	Unpacking
	3.3	Mains connection
	3.4	PC connection
	3.5	Software
	3.6	HPLC fluid connections
4.	Maintenand	e & shutdown
	4.1	Maintenance
	4.2	Shutting down the system
5.	ECD 2.1 cor	ntroller
	5.1	Introduction
	5.2	Overview of ECD 2.1 screens
	5.3	Parameters
	5.4	ClarityChrom [®] ECD 2.1 control module 34
	5.5	Dialogue Elite
6.	Detection a	nd Parameters
	6.1	Introduction
	6.2	Three-electrode configuration
	6.3	Internal organization
	6.4	Dual flow cell control
	6.5	Parameters
7.	Noise supp	ression: ADF™
	7.1	Introduction
	7.2	Frequency
	7.3	Frequency of signal and noise
	7.4	Amplitude response plot
	7.5	Applying ADF in chromatography

8.	Pulsed Ampe	erometric Detection (PAD)50
	8.1	Introduction
9.	Optimization	n of the working potential
	9.1	Introduction
	9.2	Electrochemical reactions54
	9.3	Hydrodynamic and scanning voltammogram
	9.4	Optimization using a voltammogram57
	9.5	Construction of a hydrodynamic voltammogram
	9.6	Construction of a scanning voltammogram60
10.	. Technical dat	ta64
	10.1	Environmental, dimensions, weight & power requirements
	10.2	General
	10.3	DC mode
	10.4	PULSE mode
	10.5	SCAN mode
11.	Rear panel I/	′O67
	11.1	USB B connector
	11.2	LAN connector
	11.3	VALVE connector
	11.4	ANALOG DATA connector
	11.5	Digital I/O connector
	11.6	Chassis grounding stud
12.	. Functionality	y tests
	12.1	Installation Qualification (IQ)84
	12.2	Operation Qualification (OQ)84
13.	. Troubleshoo	ting
	13.1	Instrument errors
	13.2	Analytical troubleshooting
	13.3	Dummy cell test
	13.4	Stop flow test
14.	. Reorders	
15.	. Index	

1. General information

1.1 Intended use

The ECD 2.1 is used in combination with (Ultra) High Performance Liquid Chromatography for the electrochemical detection of suitable analytes in liquid samples. With this technique the amount of electroactive substances in mobile phase can be quantified. The instrument can be used for the chromatographic analysis of a wide range of electroactive analytes in the fields of for example:

- Bioanalytical analyses
- Food analyses
- Environmental analyses



For research purposes only. While clinical applications may be shown, this instrument is not tested by the manufacturer to comply with the In Vitro Diagnostics Directive.

Operation of an electrochemical detector can involve the use of hazardous materials including corrosive fluids and flammable liquids. The instrument should only be operated by users with the following expertise:

- Completed degree as chemical laboratory technician or comparable vocational training
- Fundamental knowledge of liquid chromatography
- Participation in an installation of the system performed by the manufacturer or a company authorized by the manufacturer and suitable training on the system and chromatography software.
- Knowledge and experience in the safe handling of toxic and corrosive chemicals and knowledge of the application of fire prevention measures prescribed for laboratories.

Information on safety practices is provided with your instrument and operation instructions. Before using your instrument or accessories, you must thoroughly read these safety practices. These instructions are written for laboratory technicians who use the detector for (U)HPLC analysis.



Unskilled, improper, or careless use of this instrument can create fire hazards, or other hazards which can cause death, serious injury to personnel, or severe damage to equipment and property. Observe all relevant safety practices at all times. Only use the device for applications that fall within the scope of the specified intended use. Else the protective and safety equipment of the device could fail.

1.2 WEEE directive

All equipment of KNAUER which are subjected to the WEEE directive shipped after August 13, 2005 are compliant with the WEEE marking requirements. Such products are labelled with the "crossed out wheelie", depicted on the left site.

The symbol on the product indicates that the product must not be disposed as unsorted municipality waste.



Collection & recycling information

Please ship the instrument back to the manufacturer (KNAUER Wissenschaftliche Geräte GmbH, Germany) at the end-of-life time of the product. The manufacturer will take care of the proper disposal and recycling of the instrument at its facilities.

Shipping address for the end-of-life products:

KNAUER Wissenschaftliche Geräte GmbH

Hegauer Weg 38

14163 Berlin

Germany

In case of questions, or if further information is required about the collection & recycling procedure, please contact your local distributor.

1.3 Warning symbols

The following symbols are used in this guide:

ELECTRICAL SHOCK	This sign warns about the risk of electric shock. It calls attention to a procedure or practice which, if not adhered to, could result in loss of life by electrocution. Do not pro- ceed beyond a danger sign until the indicated conditions are fully understood and met.
WARNING	The warning sign denotes a hazard. It calls attention to a procedure or practice which, if not adhered to, could re- sult in severe injury, loss of life or damage or destruction of parts or all of the equipment. Do not proceed beyond a warning sign until the indicated conditions are fully understood and met.
	The caution sign denotes a hazard. It calls attention to a procedure or practice which, if not adhered to, could result in damage or destruction of parts or all of the equipment and/or erratic results. Do not proceed beyond a cautions sign until the indicated conditions are fully understood and met.
BIOHAZARD	The biohazard sign draws attention to the fact that use of biological materials, viral samples may carry a significant health risk.
TOXIC HAZARD	The toxic hazard sign draws attention to the fact that use of toxic solvents or samples may carry a significant health risk.
	The attention sign signals relevant information. Read this information.
0	The note sign signals additional information. It provides advice or a suggestion that may support you in using the equipment.

1.4 Safety instructions

Adhere to the following standard quality control procedures and the following equipment guidelines when using the ECD 2.1 detector. The following safety practices are intended to ensure safe operation of the instrument.

Working environment & safety



The intended use of the instrument is to detect electroactive substances in liquid samples in combination with a (U) HPLC system in a GLP-approved environment. Operators using the system should have the appropriate education an extensive understanding of GLP rules and be skilled in the art. Use this system ONLY for the intended use. Use of the system for any other purpose will cause unsafe situations.



System operation

To assure optimal performance of the detector we recommend that the instrument is checked regularly and maintenance procedures are carried out. Preventive maintenance contracts are available for that purpose. Please contact your local distributor or the nearest sales office for more information.

WARNING - RISK OF ELECTRIC SHOCK DISCONNECT POWER BEFORE SERVICING

AVERTISSEMENT - RISQUE DE CHOC ELECTRIQUE COUPER L'ALIMENTATION AVANT LA MAINTENANCE




Electrical safety

The removal of protective panels on the instrument can result in exposure to potentially dangerous voltages. Therefore, disconnect the instrument from all power sources before disassembly.



Untrained personnel should not open the instrument, **this may only be done by authorized service engineers.** Replace or repair faulty insulation on power cords immediately after discovery of the fault. Check that the actual power voltage is the same as the voltage for which the instruments are wired. Make sure power cords are connected to correct voltage sources: grounded AC power source, line voltage 100 - 240 VAC. The instrument should be connected to a protective earth via a ground socket. The ECD 2.1 must only be used with appliances and power sources with proper protective grounding to prevent damage through build-up of static electricity. The power source should exhibit minimal power transients and fluctuations. If necessary connect to a filtered mains socket.

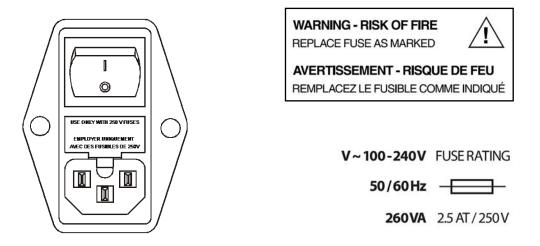

Replace blown fuses with fuses of proper type and rating as indicated on the rear panel and as listed in the list of accessories and spares (appendix D). The fuse holder is integrated in the mains connector. Ensure that the instrument is never put in operation with fuses of a different type. This could cause fire. Only use manufacturer-supplied I/O cable(s) to connect with other devices. Thoroughly connect shielding to common. Manufacturer will not accept any liability for damage, direct or indirect, caused by connecting this instrument to devices and with cables which do not meet relevant safety standards.



Solvents

The solvents used may be flammable, toxic or corrosive. The room in which the system is installed should be well ventilated to prevent that solvent vapors cause poisoning or ignite and cause a fire. Use of open fire in the vicinity of this system must be strictly prohibited. Do not install the system in the same room with any other equipment that emits or could potentially emit sparks. Provide protective equipment near the instrument, when solvent gets into the eyes or on the skin, it must be flushed away immediately. Provide equipment, such eye wash stations and safety showers, as close to system as possible. Use proper eye and skin protection when working with solvents. Additional safety requirements or protection may be necessary depending on the chemicals used in combination with this equipment. Make sure that you understand the hazards associated with the chemicals used and take appropriate measures with regards to safety and protection.

Sample containers (vials) should be sealed to minimize any risks related to solvent vapor.

Biological hazard



When you analyze biological fluids you need possible precautions and treat all specimens as potentially infectious. Always wear protective gloves when handling toxic or biologically infectious samples to prevent bio hazards or hazards while working with the ECD 2.1. If necessary the instrument must be decontaminated before decommissioning or shipment of the instrument for repair to KNAUER or its representatives. When shipped to KNAUER every instrument has to be accompanied with a decontamination form which should be completely filled in and signed by the customer. Without this decontamination form the instrument will not be processed by KNAUER (either repaired or disposed).

Waste disposal



Perform periodic leak checks on LC tubing and connections. Do not close or block the drain in the oven compartment. Do not allow flammable and/ or toxic solvents to accumulate. Follow a regulated, approved waste disposal program. Never dispose of flammable and/ toxic solvents through the municipal sewage system.

Applications: quality control



It is recommended that you routinely run several quality control samples. Quality control samples should represent low, average and high levels of a compound. Make sure that quality control sample results are within an acceptable range, and evaluate precision from day to day and run to run. Data collected when quality control samples are out of range may not be valid. Do not report this data until you are certain that system performance is acceptable. Apart from use of quality control samples, we recommend that you use blanks. The blanks will help you assess whether carry-over is within an acceptable range and monitor the integrity of your data.



Using the ECD 2.1 in other ways than indicated in the instructions or defined by good laboratory practice may result in erratic or unsafe operation.

2. Introduction

Congratulations on your purchase of the AZURA® Detector ECD 2.1. This detector enables you to perform all (U)HPLC applications using electrochemical detection. The ECD 2.1 includes a highly stable Faraday-shielded oven compartment accommodating column and flow cell. This flow cell has surprised researchers for its unsurpassed S/N ratio and therefore you now possess the best possible combination for extremely sensitive EC analyses. The ECD 2.1 has 3 operational measurement modes: DC, SCAN and PULSE mode. Furthermore, a Service, Diag(nostics) and Config(uration) mode are available. In addition, crucial parameters can be controlled by either relays or TTL. The ECD 2.1 can support up to 4 flow cells (optional), which makes it possible to perform 4 independent measurements with one detector.



These instructions cover the installation, set-up and operation of the AZURA® Detector ECD 2.1 only. Detailed operation instructions for other peripheral LC equipment and parts such as flow cells, pumps, auto samplers, valves, column heaters etc. are given in the instructions accompanying those accessories. From this point onwards the abbreviation ECD 2.1 will be used.

2.1 Instrument description

N ECD 2.1 Azura 2 8 3 (2 3 0 3 4 5 6 KNAUER Fig. 1 Front view

Legend

- Instrument housing
- ② LC tubing inlet/outlet
- ③ Instrument door panel
- 4 x 40 Ch LCD display
- 5 Function keys
- 6 <Enter> key
- ,+' and ,-' value keys
- ⑧ Cursor keys
- Door handle (for opening door)


Legend

- Instrument rear panel
- ② Type label (pn, sn etc.)
- ③ Digital I/O connector (25-pins sub-D fem)
- Analog data (9-pins sub-D fem)
- 5 Valve connector (9-pins sub-D male)
- 6 LAN connector (RJ45 jack)
- ⑦ USB connector (USB B)
- 8 Fuse & power rating
- 9 Mains switch/inlet
- Grounding stud
- I Fuse compartment
- Ventilation holes



Legend

- Cell cabinet
- 2 Cell connector (9-pins sub-D fem)
- ③ Top fan heater (intake)
- ④ Door sensor
- (5) Mounting plate (for cells & columns)
- 6 Doorlock
- ⑦ Flow cell clamp (for SenCell)
- ⑧ Column clamp
- In Mounting hole for cell clamp (M4)
- Image: Bottom fan heater (exhaust)
- ① Mounting hole for column clamp (M3)
- I Fuse compartment
- 1 Door panel, rear
- 🕼 Type label

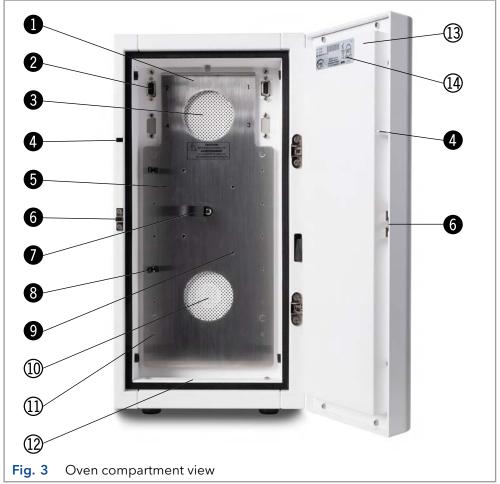

AZURA® Detector ECD 2.1 Instructions, V6685

3. Installation

3.1 Site preparation requirements

For a successful onsite installation of the instrument, please arrange the following requirements at your location in advance:

Personal Computer

In case the instrument is used via remote control by PC software (Dialogue Elite, ClarityChrom®) or firmware (FW) update needs to be performed a desktop computer is required with the following requirements:

- Free LAN port (onboard, PCI, PCI express or PCI-X)
- Free USB port (required for FW updates)
- Microsoft Windows 7 Operating System or higher

Detailed information about the use of the ECD 2.1 in combination with ClarityChrom[®] chromatography software are listed in the ClarityChrom[®] instructions (document no. V1670, Website: www.knauer.net/cc-instructions) and in the list of supported drivers and

<u>www.knauer.net/cc-instructions</u>) and in the list of supported drivers and software packages for all KNAUER devices (document no. V1667, Website: <u>www.knauer.net/softwarecontrol</u>).

Laboratory requirements

Your instrument is intended for indoor use only in an industrial or commercial environment (EN55011 group 1 class A ISM equipment). It is suitable for the following categories: Installation category II, Pollution degree 2, equipment class I.

Table I. Environmental specifications

Parameter	Requirement
Operating temperature	10 - 40 °C (50 - 104 °F)
Operating humidity	20 - 80%, non-condensing



For optimum analytical performance it is recommended that the ambient temperature of the laboratory is between 20-25 °C and is held constant to within ± 2 °C throughout the entire working day. Note: for optimal temperature stability of the cell cabinet the oven temperature should be set to a temperature at least 7 degrees higher as ambient temperature. Requirements for the laboratory bench on which the instrument will be installed:

- Stabile, clean, flat and smooth surface.
- Enough mechanical strength to hold at least the weight of the detector: a full-dressed detector with flow cells, columns and valves can weigh up to 20 kg.
- A detector has the following dimensions 44 (D) x 22 (W) x 43 (H) cm = 17.3" (D) x 8.7" (W) x 16.9" (H). Take into account that additional space is necessary on all sides (at least 10 cm) to prevent obstruction of ventilation holes and allow sufficient heat dissipation.

Chemicals

Mobile phase and flush/storage solutions must be clean as it is in direct contact with the working electrode in EC detection. High purity of chemicals including water is a prerequisite. So all chemicals should be electrochemically clean, HPLC grade or better. For water used for the preparation of mobile phases a water purification apparatus is advised which is able to supply high purity deionized water with resistivity of >18 MOhm.cm and low TOC level (<10 ppb).


3.2 Unpacking

Inspect the transport box for possible damage as it arrives. Immediately inform the transport company in case of damage, otherwise it may not accept any responsibility. Keep the transport box as it is designed for optimum protection during transport and it may be needed again. Carefully unpack the system and inspect it for completeness and for possible damage. Contact your supplier in case of damage or if not all marked items on the checklist are included. Prior to shipment, your detector has been thoroughly inspected and tested to meet the highest possible demands. The results of all tests are included.


1	Delivery is in accordance with order	
2	Delivery is undamaged	
3	All items on checklist(s) are included	
	Certificates of performance are included:	
4	- detector	
4	- detector - flow cell(s)*	

*) Note that flow cells are not part of the detector and have to be ordered separately.

To unpack the detector, lift it from its box by both hands (Fig. 4). **Never lift the detector at its front door**, but at its sides.

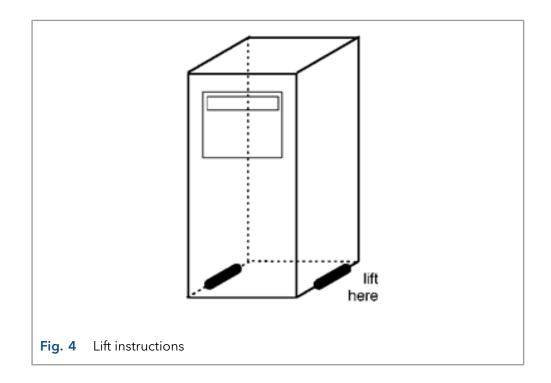

With both hands under the instrument lift the detector to its operation location. Install the detector in an area which meets the environmental conditions.

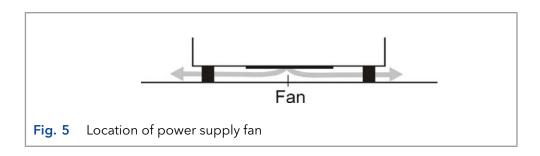

Remove the protective tape from the detector LCD screen. Leave the instrument to adopt ambient temperature for at least half an hour in the place of installation.

Use the detector indoors only. Place the detector upright (on its instrument feet) on a stable, flat and smooth surface. Do not place the instrument in an area subject to excessive dust or shocks.



Do not place it near a source of heat or in direct sun light, as this may influence the heating capabilities of the instrument. Make sure the detector is placed in such a way that the mains power connection can be reached easily to disconnect it from the mains power by removing the mains power cable. Do not block the ventilation holes at the back of the instrument and do not block the fan located at the bottom of the detector (Fig. 5). Blocking the fan will impair the cooling capability of the power supply. Do not place object/instruments on top of the detector. Objects can be placed on any side of the detector; however, make sure these objects are placed at a distance of 5 cm from the detector, in case that objects are placed at only one side of the instrument. and 10 cm from the instrument, if objects are placed on both sides of the instrument.

3.3 Mains connection

Check that fuses and voltage range on the rear side of the instrument match that for the power outlet to be used.



This device complies with the requirements of EN 55011 group 1 class A ISM equipment. In a domestic environment this product may cause radio interference in which case the user may be required to take adequate measures on his own expenses. Understanding power surges. Power surges, line spikes, and transient energy sources can impede instrument operation. Ensure that the instrument's electrical supply is adequately protected from these conditions and properly grounded. If necessary connect the instrument to a filtered mains inlet.



Power supply and protective earth: The system must be connected to a suitable mains power supply with a correctly installed protective earth conductor. Never use the system without a properly connected protective earth conductor.

Leave the instrument powered off until specifically mentioned in the procedure below.

3.4 PC connection

Follow the instructions in this paragraph when the instrument is used with PC control over LAN using the Dialogue Elite or ClarityChrom® Chromatography Software. This section can be skipped if the instrument is used stand-alone. To be able to communicate over LAN a computer is required with a free (PCI, PCI Express or PCI-X) LAN port.

The ECD 2.1 has a fixed IP address: 192.168.5.1, with subnet mask: 255.255.255.0. Gateway and DNS are not filled in.

The instrument is standard delivered with a special crossover LAN (UTP) cable which is part of the ECD 2.1 accessory kit that is shipped with the detector.




To insure stable and error-free communication only use the manufacturer-supplied LAN connectivity kit to connect the ECD 2.1 to LAN. Create a small dedicated local area network to connect the ECD 2.1 to the PC. Do not connect the ECD 2.1 over a company Local Area Network. If needed a second network adapter with a different (unique) IP address range can be applied. In the following section the procedure to connect the instrument to the PC using the crossover UTP cable is described. Configure the IP address of the PC LAN network card by executing the following steps (in this example Windows 7 OS is used, in Windows 8 the example screens may look different):

 Right click on the Network icon in the bottom right of the Windows taskbar and open the menu ,Open Network and Sharing Center'. Alternative: open Windows start menu, open control panel and go to ,Open Network and Sharing Center'):



- **2.** Open the menu ,Change adapter settings' on the left panel of the ,Open Network and Sharing Center' window.
- **3.** Right click on the Local Area Connection icon of the LAN card in your PC and click on properties to open the Network card setting.

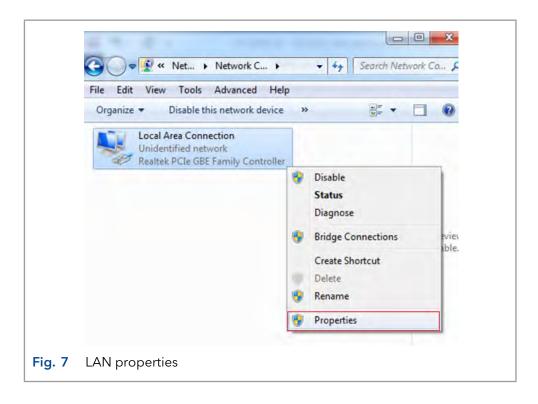


Networking	General	
Connect using:		automatically if your network supports eed to ask your network administrator
Configure	Obtain an IP address autom	atically
This connection uses the following items:	Use the following IP address	s:
Client for Microsoft Networks	IP address:	192.168.5.10
File and Printer Sharing for Microsoft Networks	Subnet mask:	255 . 255 . 255 . 0
Internet Protocol Version 6 (TCP/IPv6) Internet Protocol Version 4 (TCP/IPv4)	Default gateway:	
 ✓ Link-Layer Topology Discovery Mapper I/O Driver ✓ Link-Layer Topology Discovery Responder 	 Obtain DNS server address Use the following DNS server 	
Install Uninstall Properties	Preferred DNS server:	
Description	Alternate DNS server:	le se e
Transmission Control Protocol/Internet Protocol. The default wide area network protocol that provides communication across diverse interconnected networks.	Validate settings upon exit	Ad <u>v</u> anced
OK Cancel		OK Cance

- **4.** Open the menu ,Internet Protocol Version 4 (TCP/IPv4)' menu (double click).
- Configure the network IP address and subnet mask as depicted in the screen dump below (IP 192.168.5.10, Subnet mask: 255.255.255.0). Gateway and DNS fields are not filled.
- **6.** Close the menu(s) by clicking the ,OK' buttons. The network IP address of the LAN network card is now set up for communication with the ECD 2.1.
- **7.** Connect the crossover (UTP) cable to the RJ45 Jack of the LAN card of your PC (typically located on the backside of a desktop PC).
- **8.** Connect the other end of the crossover LAN (UTP) cable to the LAN port on the rear panel of the ECD 2.1 as depicted on the photograph.
- **9.** Switch on the ECD 2.1. Set the detector temperature to 35 °C if a Performance Verification (PV) will be executed, or set it to the temperature at which your application is running. Allow the instrument to stabilize for at least one hour before starting analysis.



Functionality tests like OQ and PV are described in separate instructions. Please contact the Customer Support for more information.

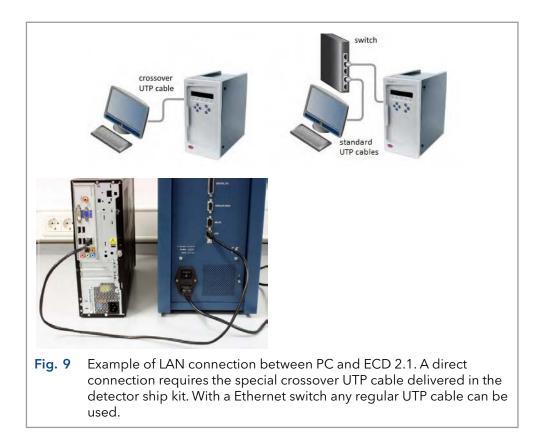

3.5 Software

The ECD 2.1 can be used remote in combination with PC control software. Currently, there are two software packages available for control of the ECD 2.1 electrochemical detector:

- Elite Dialogue software from Antec, The Netherlands.
- ClarityChrom[®] Chromatography software (Version 6.x and higher) from DataApex, Czech

The Elite Dialogue software is also required for the upload of new firmware (= embedded software controlling the ECD 2.1 electronics). In this section the installation and configuration of the ECD 2.1 is shortly described. Note: this is by no means a replacement of the installation documentation available for the software packages. Please refer to this documentation for details.



Ensure you have Administrator access rights in your system before you start with the installation of the software packages. Elite Dialogue and ClarityChrom[®] users must have read/write access to all software folders and subfolders.

Dialogue Elite

To install the software:

- Download the latest version of the Dialogue Elite from the Antec website <u>www.myantec.com</u> (register to get access).
- Double click on the setup.exe file to start the installation wizard.
- Follow the instructions of the installation wizard for successful installation of the software.
- Insert the Dialogue Elite license dongle to get full access to the software (without dongle it will operate in demo mode).
- Make sure that the LAN connection is configured and the LAN cable is connected.
- Power up the detector by means of the mains switch on the rear panel.
- Start the program Dialogue Elite from the Windows start menu.
- During start-up the 'Select devices' menu will pop-up as shown in Fig. 10.

Select devices	Devices	Port settings port scan
DECADE Elite	DECADE Elite	192.168.5.1 👻 🛄
DECADE II DECADE II SDC ROXY ROXY HC Syringe pump		192.168.5.1 / 00 80 A3 9C 85 40 192.168.5.129 / 00 80 A3 A3 64 79 192.168.5.221 / 00 80 A3 A3 69 5F
Connect Bite via USB	Cancel	ОК

- When a ECD 2.1 detector is available it is automatically detected and the IP address is shown in the port settings box. If not press port scan or type in the default IP address 192.168.5.1.
- The pull down field shows all responding devices with their IP and MAC address. In case of doubt check the device unique MAC address on the rear panel IO connector.
- Type OK. The instrument will connect and is ready for use.

ClarityChrom[®] chromatography software

You find a detailed description on chromatography software in the corresponding software instructions (document no. V1670) which can be found on the KNAUER website: <u>www.knauer.net/claritychrom</u>

You find more information about supported software packages and firmware versions in the list "KNAUER software device support" (document no. V1663)which can be found on the KNAUER website: <u>www.knauer.net/softwarecontrol</u>

3.6 HPLC fluid connections

In this section the installation & priming of all relevant fluidic connections are described to be able to use the ECD 2.1 for the analysis of substances with (U)HPLC-ECD 2.1. When working with HPLC solutions and mobile phases take the following precautions:



Use proper eye and skin protection when working with solvents. The solvents used may be flammable, toxic or corrosive. Organic solvents are toxic above a certain concentration. Ensure that work areas are always well-ventilated! Use of open fire in the vicinity of this system must be strictly prohibited. Do not install the system in the same room with any other equipment that emits or could potentially emit sparks. Wear protective gloves, safety glasses and other relevant protective clothing when working on the device!

With respect to third-party (U)HPLC equipment, such as LC pumps, auto samplers, injection valves, column heaters etc. used in combination with the ECD 2.1 detector: The equipment connected to the system should be specifically designed for use in (Ultra) High Performance Liquid Chromatography and capable of delivering flow rates typically in the range between 1 μ L/min up to 5 mL/min.



The manufacturer will not accept any liability for damage, direct or indirect, caused by connecting this instrument to devices that do not meet the relevant safety standards.

Tubing connectors

For optimal operation it is of the utmost importance that all tubing connections on the injection valve, columns and flow cells of the injection valves are made without introducing dead volumes to minimize peak broadening, carry-over etc.




Use only the original polymeric fingertights supplied by the manufacturer with the purchased flow cell to make LC connections on the flow cells inlet and outlet. Do not use metal tubing on the flow cell because it may lead to damage or incorrect operation of the flow cell. Use PEEK, PEEKsil or Fused Silica tubing (with FEP sleeves). For columns and injection valves etc. use only nuts, ferrules or finger tights recommend by the respective manufacturer of the parts. The use of unsuitable connectors may lead to damage of the parts or the introduction of dead volumes.

Please note that the tubing length (length L below) required to make a good connection differs for each brand of connection. If length L is not correct, this will result in faulty peaks and carry-over. Essentially, when you create a connection the ferrule on the tubing is compressed into the valve to make sure that the connection is leak-tight. Take the following into account when creating the connection:

If L is too long, the ferrule cannot form a seal in the connection. This may cause irreparable damage port of the valve, column or other part, as:


Part of the tubing may end up in the connection port internals

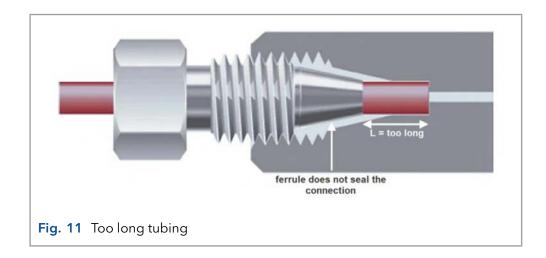

Internals of the port may be damaged.

If L is too short, this may result in:

- Leakage
- Dead volume at the end of the ferrule (a 'mixing chamber').

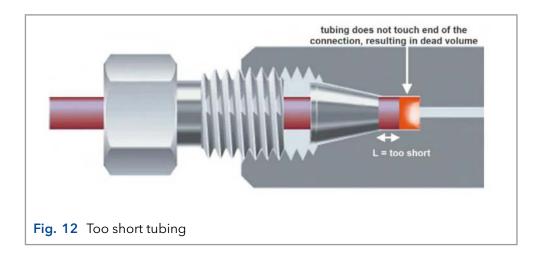

Every ferrule type needs an appropriate length of tubing for connecting it to the type of connection port, depending on the depth of the connection port. Refer to the information provided by the manufacturer for specific information.

18

Mobile Phase

Electrochemical detection is a sensitive detection technique characterized by extremely low detection limits. A typical detection limit of 100 pmol/L or lower for catecholamines is no exception. Improving detection limits will always be limited by the weakest link in an LC-EC system. In daily practice a couple of 'rules' must be obeyed to fully exploit the incredible linear dynamic range and low detection limits of an EC detector. These are not only hardware related, but also refer to mobile phase composition, degassing, temperature and pH stability and several other issues.

Mobile phase requirements:

- Electrochemically clean, HPLC grade or better
- Ion strength 20 200 mmol/L
- Buffer pH near pKa
- In-line 0.2 μm filter & degassing of mobile phase
- EDTA for trapping of metal ions

Mobile phase must be clean as it is in direct contact with the working electrode in EC detection. High purity chemicals including water is a pre requisite. In some applications EDTA is added to the mobile phase to traps electrochemically metals such as Fe2+ by forming an inactive complex.

However at higher working potentials (typically > 1.2 V vs. salt bridge AgCl ref) also EDTA can become electrochemically active and is not recommended. In that case only a passivation step with 15% HNO3 is recommended (see page 20 and following).

Electrolytes assure contact between 3 electrodes in an electrochemical flow cell. Low ion strength destabilises an EC system and noise will increase. Extremely high buffer concentrations cause problems of salt formation. Therefore concentrations between 20 and 200 mmol/L are recommended.

Also constant pH is important for baseline stability and reproducible results. Stability of pH is best when close to pKa of a buffer ion. Often used buffers are phosphate, carbonate, acetate and citrate. Modifiers such as methanol, propanol and acetonitril can be used without problems in DC amperometry, but not in pulsed amperometric detection as peaks are strongly attenuated. In our experience the quality and expiration date of organic modifiers can be a problem resulting in increased noise levels.

Metal inlet frits in mobile phase bottles are not advised as metal ions are introduced and baseline noise is increased. Instead 0.2 μ m nylon inlet filters are recommended.

Vacuum degassing: Considerable amounts of the gases N2, O2 and CO2 may be dissolved in HPLC mobile phases. Whenever the temperature changes, solvents are mixed or a pressure reduction occurs, these gases may show up as very small air bubbles. To avoid noisy baselines an in-line vacuum degasser with a low dead volume (less than 0.5 mL) and high degassing efficiency should be added to the system. The advantage over helium degassing is that the degasser does not change the mobile phase composition.

Helium degassing: Degassing using helium is an effective and universally applicable method but only recommended when working in reductive electrochemical detection and pulsed electrochemical detection (analysis of carbohydrates using anion-exchange chromatography with NaOH as mobile phase). All gases except helium are removed completely. Helium is not EC active and does not change the mobile phase properties significantly. To prevent mobile phase contamination, only high-purity helium must be used.

Installation & start-up

Please carefully follow the next steps for a successful installation and start-up:

- 1. The installation of the flow cell and column is shown in Fig. 13.
- If a manual injector is applied with position sensor, the sensor cable must be connected to the VALVE connector on the rear panel to enable INJECT/LOAD functions. In the CONFIG screen the valve option should be set to ,Valve = present' manually.
- 3. An electrically actuated valve of VICI Valco (type E2CA, EHCA) can be connected & controlled by the ECD 2.1. For that purpose a serial valve cable is required. The serial cable should be connected to the VALVE connector on the rear panel of the instrument to enable INJECT/LOAD functions. This presence of this valve is automatically detected and no further configuration has to be performed.



4. Prior to connection of the HPLC system to the detector all metal parts should preferably be passivated with 15% nitric acid during 20 min. The acid is flushed through the pump, the pump tubing, the dampener, the injector (in load and inject position) and to waste.



Make sure that all parts that are not acid-resistant such as: nylon inlet filters, column and flow cell are not connected during this step.

5. After flushing with nitric acid, the system must be thoroughly flushed with demi water. Make sure that no traces of nitric acid are left in the tubing or pulse dampener (check with pH paper). Flush the system with HPLC buffer.



If an ISAAC[™] reference electrode is used, make sure that the mobile phase contains at least 2 mmol/l chloride (KCl or NaCl) ions.

- 6. Before connecting a new column read the manufacturer's instructions. Our experience is that thorough pre-conditioning of a column is always required. Only a pre-conditioned column is electrochemically clean. If not, the background current may be unacceptably high and substantial fouling of the working electrode occurs. For reversed phase columns flushing with 50% methanol in water for 3 days at a low flow rate is highly recommended. Before switching to mobile phase, flushing with water (10 column volumes) is recommended to prevent precipitation of buffer salts.
- 7. Passage of air bubbles through the flow cell will lead to unacceptable noise levels and 'spikes'. Therefore, the use of an in-line degasser is strongly recommended. In our experience, a one-time degassing step of the HPLC buffer is almost never sufficient. If the ECD 2.1 is used for reductive ECD (at a negative working potential) additional steps should be taken to remove oxygen from the mobile phase. These include degassing with Helium and the use of stainless steel tubing (impermeable for oxygen).
- 8. Consult your flow cell instructions for detailed information about the installation of the flow cell. See Fig. 14 on the next page for reference. Connect the flow cell to the corresponding cell connector in the oven compartment. All cell connectors are marked with a label for identification. For the ECD 2.1, connect the flow cell to the cell connector on the left side marked "Cell 1".




The cell connector inside the oven compartment is ESD sensitive. Make sure that the flow cell is OFF when removing or connecting the cell cable.

Never switch ON the flow cell when: (1) the cell cable is not correctly connected, (2) the cell is only partly (or not at all) filled with mobile phase, (3) the outside of the flow cell is wet, particularly the part between the auxiliary and working electrode connection, because substantial damage to the working electrode or electronics may occur.

Legend

- Cell clamp
- Cell Inlet (tubing connection from column-to-cell)
- Cell outlet (tubing connection from cell-towaste)
- ④ WE contact (red)
- 6 AUX contact (blue)
- 6 REF contact (black)

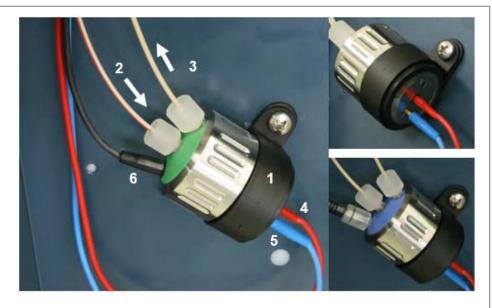

- Fig. 14 Left: SenCell with ISAAC reference mounted under an angle of approximately 45° in the detector. Make sure that the outlet [3] is positioned on the top side to prevent entrapped air bubbles. <u>Top-right</u>: electrical connections of WE (red connector) and AUX electrode (blue connector). <u>Bottom-right</u>: SenCell with SaltBridge reference electrode.
- **9.** Before switching ON the flow cell, make sure that the mobile phase contains sufficient electrolyte (buffer ions). A stable baseline will never be obtained if the cell is switched ON with only water or another non-conducting mobile phase. Also be sure that no air bubbles are trapped in the flow cell.
- **10.** The outlet tubing from the flow cell should lead to a reservoir that is at a higher level than the flow cell. This ensures a small back pressure which prevents air-bubble entrapment. The outlet tubing should be under the liquid level, to avoid electrical noise induced by 'dripping' of mobile phase.
- **11.** Set the cell potential (see page 54 for optimization of the potential), switch ON the flow cell and allow the system to stabilize for approximately 30 min. A 'good' stabilization curve shows a mono-exponential decline without jumps and/or spikes.
- 12. Connect the ECD 2.1 to the PC control software (ClarityChrom® or Elite Dialogue). In the case the instrument is used stand-alone connect an external A/D converter or recorder to the 'Analog data output' (1 V full scale) on the rear panel. Refer to the documentation of the third-party software controlling the A/D converter for detailed instruction how to set up a measurement with the A/D device.

Your system is now ready for use. The ECD 2.1 has been developed for continuous operation. For maximum stability it is advised to leave the system ON continuously. If preferred, the flow cell may be switched OFF at night.

4. Maintenance & shutdown

4.1 Maintenance

In this paragraph all maintenance is described which can be performed by the end-user, all other maintenance & service procedures may be performed by authorized service engineers only.

Periodic check for leakage

Perform leak checks on LC tubing, flow cells and connections on a daily basis and check if the drain on the bottom of the oven compartment is not blocked or closed. Do not allow flammable and/or toxic solvents to accumulate. Follow a regulated, approved waste disposal program. Empty and clean waste container regularly. Never dispose of such products through the municipal sewage system. Check daily that the mobile phase bottles contain enough mobile phase for the number of analyses planned to be executed.

Periodic check of the oven temperature

The operator should perform regular checks to verify if the actual oven temperature is in accordance with the set temperature of the ECD 2.1.



In case the actual temperature exceeds 70 °C switch off the detector and contact the manufacturer or its representatives for service.

Flow cell



Check the performance of the detector & flow cell on a daily basis by evaluating background current, noise and signal. An increase in background current, noise and/or loss of sensitivity may be a sign of contamination of the working electrode (WE) and/or a sign that maintenance is required on the reference electrode (REF) of the flow cell. If necessary perform maintenance on the flow cell. Detailed information about flow cell maintenance instructions are described in the instructions supplied with the specific flow cell (SenCell, VT-03 or FlexCell) purchased in combination with your detector.

Cleaning

In general, the ECD 2.1 needs very little maintenance. The outside of the detector may be cleaned with a non-aggressive cleaning liquid.




Do not use any organic solvents to clean the exterior of the detector, because this may lead to damage of the paint layer.

In case of leakage in the cell cabinet (tubing, connectors, cell, column etc.) remove the spilled mobile phase or other solutions as soon as possible because this may damage the paint layer, or result in the deposition

of salt crusts (in case of buffered mobile phases), which could block the drain in the bottom of the cell cabinet. Remove any dust on the protective screens that cover the fans in the oven compartment.

Replacement of fuses

Replace blown fuses with fuses of proper type and rating as stipulated on the rear panel and specified in the installation section of these instructions. The fuse holder is integrated in the mains connector. Ensure that the instrument is never put in operation with fuses of a different type. This could cause fire.

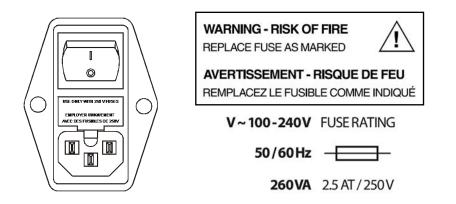

In case the fuses blow out repetitively contact KNAUER or its representatives for instructions and/or service of the instrument.

4.2 Shutting down the system

- There are a couple of steps to take to switch off an LC system with electrochemical detector for a longer period of time. Shutting down is not different from most other HPLC systems. Perform the following procedure:
- Switch off the flow cell using the keyboard (standalone) or via the software (ClarityChrom[®] or Dialogue Elite).
- Check the column(s) documentation for the appropriate storage liquid, apply this and make sure the column is properly flushed. A reversed phase C18 column is usually stored with 50% Acetonitrile/water.
- Take out the column, mount the corresponding end-caps and store the column in an appropriate place.



Avoid precipitation of high salt concentrations in organic solvent, first wash out salts with water if necessary.

- Flush and store the system with 50% water/acetonitrile (or methanol). Switch the injector valve between load and inject a few times. Make sure all tubing, filters etc. are flushed so no traces of salt are left that could precipitate and clog the system.
- Remove the flow cell from the system by disconnecting the inlet and outlet capillary.

- Open the cell, flush with water, use some tissues to carefully dry the cell. Be careful not to damage the spacer in the case of a VT-03 or Flex-Cell (the SenCell does not have a spacer).
- Close the cell and store dry. In case of salt bridge REF, store the REF separately with a cap on it to prevent drying out. Alternatively, put the sb-REF in a 10 mL vial under a KCl solution and close the vial with a cap.
- Switch off the detector (and other LC equipment) via the mains switch (switch to position '0') on the rear panel.


5. ECD 2.1 controller

5.1 Introduction

The ECD 2.1 has been designed for maximum functionality and ease of use. The control of ECD 2.1 parameters via the keyboard & LCD display is designed in a way that without reading this chapter, it should be possible to operate the detector. This chapter is intended as a reference guide in case questions arise during operation.


The information shown in the numerous screens is presented in alphabetical order. For each item an explanation is given, together with the item's nature and the screen(s) of appearance. The nature of an item can be:

- 1. Control: parameters with a cursor box ('□') can be attained via cursor buttons and changed by the 'value' button.
- 2. Status: without a cursor box a parameter reflects the current status.
- **3.** Functions: parameters in CAPITALS are commands accessible via function buttons F1 F5.
- **4.** The ,Enter' button is only used to accept changes in cell potential. In the top right corner of each screen the name of the present screen is displayed. If available, the bottom left function button displays a previous screen, and the bottom right one the next screen.

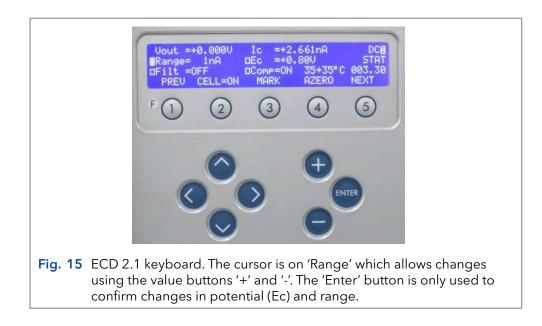


5.2 Overview of ECD 2.1 screens


DC mode

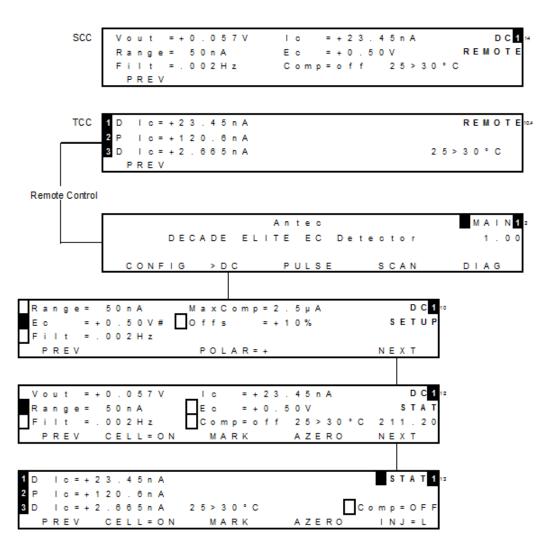


Pulse mode

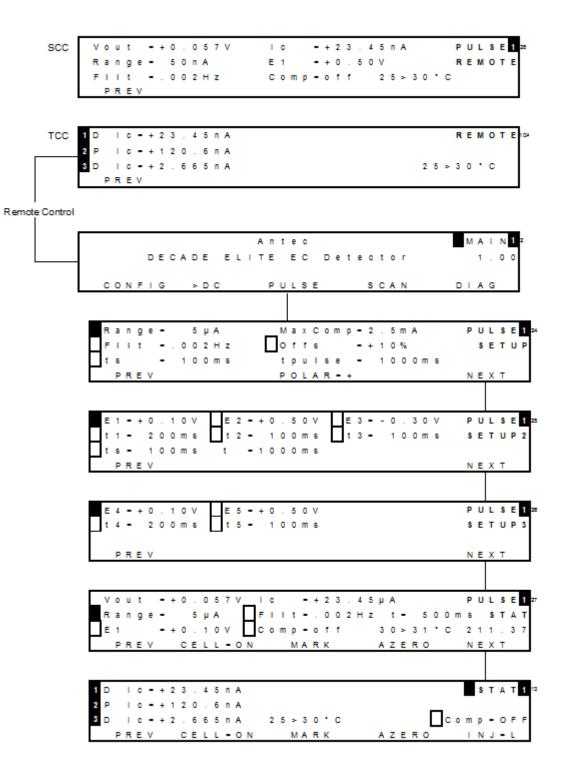


Scan mode

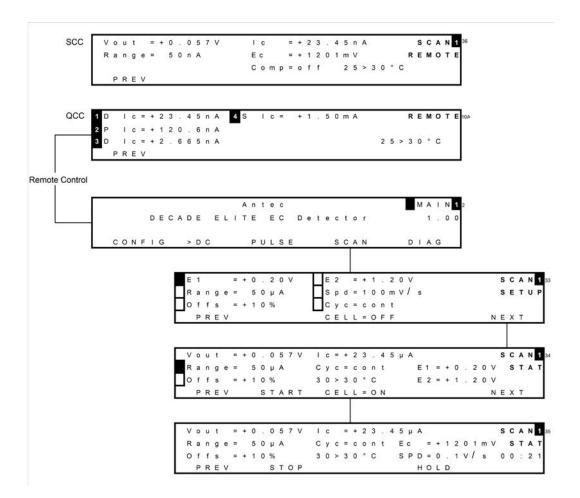

CONFIC menu

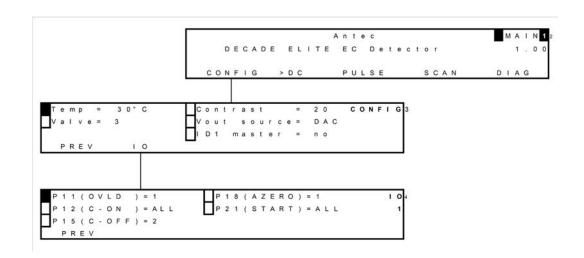

DIAG menu

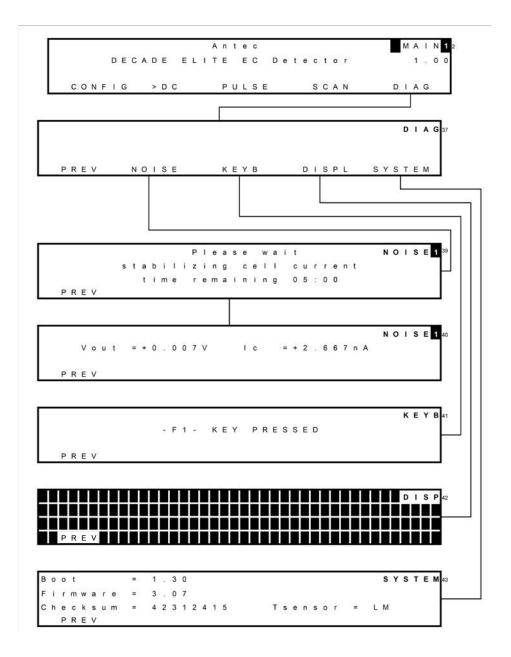

5.3 Parameters

Explanation: Type S is status, F is function and C is control.

Parameter	Screen	Description	Туре
28 > 30°C	dc stat pulse stat scan stat run	Displays the actual (left value) and the pre-set oven tempe- rature (right value).	S
AZERO	dc stat, run, pulse stat, scan stat	Sets the output voltage to 0 V, or to the offset voltage. Control Comp = off changes to Comp = on. If cell current exceeds the max. compensation a message "cell current exceeds max. compensation" appears. In that case max. compensation will be applied, which may not be the 0 Volt level but higher.	F
Boot	system	Displays boot firmware version	S
CELL=ON/ OFF	dc stat, pulse stat, scan set- up, scan stat	Toggles between cell 'ON' and 'OFF'. Confirmation is required "Switch cell on (off)?" Switching on resets the clock to 0.00. Pulse mode: pulsation occurs as long as the cell is on, irrespective which screen is selected. Scan mode: potential E1 is applied.	F
Checksum	system	Displays checksum	S
Comp	dc stat, pulse stat	Toggles between 'ON' and 'OFF', releases auto zero offset. Switches ON if AZERO is pressed. Affects auto zero com- pensation only, not the % offset!	С
CONFIG	main	Enters config screen	F
Contrast	config	Sets the contrast of display	С
Сус	scan setup	Controls the nature of the cycle: half, full and continuous. 'Half' means that the cell potential runs from E1 to E2 and stops at E2 (/). 'Full' means that the cell potential runs from E1 to E2, and back to E1, and then stops (/\). 'Cont' means that the cell potential runs from E1 to E2 and back to E1 continuously (/\/\/\/). Pressing "STOP" or finishing the cycle, sets the potential to E1.	С
DIAG	main	Enters Diag screen	F
DISPL	test	Enters DISP screen for display test.	F
E1, E2, E3, E4, E5	pulse setup2 pulse setup3	Controls the cell potential settings of the pulse. Can be set between +2.50 and -2.50 V with 10 mV steps. Can only be set or changed after confirmation with the 'enter' button.	С
Ec	dc setup	Controls the cell potential, can be set between +2.50 and -2.50 V with 10 mV steps. Can only be set or changed after confirmation with the 'enter' button.	С
Ec	run (dc only), scan stat (du- ring scanning)	Reflects the set cell potential. Displays the actual cell poten- tial in the scan mode.	S
Filt (DC mode)	dc setup, dc stat	Filter settings: RAW (100 Hz), Off (10 Hz) and 1 Hz to 0.001 Hz cut off frequency, in 1, 2, 5 steps.	С

Parameter	Screen	Description	Туре
Filt (PULSE mode)	pulse setup, pulse stat	Filter settings: Off and 0.5 Hz to 0.001 Hz cut off frequency, in 1, 2, 5 steps. (Fcut-off / filter coefficients based on 1 Hz input frequency in pulse mode)	С
Filt	run	Reflects the actual filter setting.	S
Firmware	system	Displays firmware version	S
Hold resume	run, scan stat	Toggle, holds or resumes execution of scan.	F
HOLD=0,1	run, scan stat	Holds or continues execution of scan. Toggles between 1 and 0. Pressing hold again continues scan were it has been hold.	F
lc	stat (dc, pulse, scan), , run, noise	Displays the true, non-compensated cell current, unaffected by auto zero or offset.	S
ID1 master	config	Sets sensor board 1 as master. When this setting is set to 'yes' all parameter settings from sensor board 1 are automa- tically copied/transferred to all other sensor boards present.	С
IO	Config	Enter IO menu	F
INJ=I/L	dc stat, pulse stat	Displays or switches the position of the injection valve, toggles between inject (I) and load (L). If a manual injector with position sensor is applied, it echoes the position of the injector. If an electrically actuated injector is used (optional) it is possible to switch the injector with this function button.	F/S
KEYB	test	Enters 'KEYB' screen, for keyboard test. Press 2x F1 to leave.	F
MARK	dc stat, pulse stat	Triggers a marker signal on output.	F
MaxComp	dc setup, pul- se setup1	Maximum cell current that can be compensated for using auto zero.	S
Next	several screens	Enter next screen	F
NOISE	test	Enters NOISE screen for performance test.	F
Offs	dc setup, dc stat, prog, pulse setup1, pulse stat, scan setup, scan stat	Percentage offset, can be set between -50 and +50%.	С
POLAR	dc setup, pul- se setup2	Inverts output polarity, toggle between + and Requires confirmation.	F
PREV	several screens	Return to previous screen	F
P11(OVLD)	IO	Programmable output: can be configured that the overload (OVLD) signal of cell 1, 2 or 3 only is present on pin 11 when active or ALL cells.	С
P12(C-ON)	IO	Programmable input: can be configured that only cell 1, 2 or 3 is switched ON when active, or ALL cells.	С
P15(C-OFF)	IO	Programmable input: can be configured that only cell 1, 2 or 3 is switched OFF when active, or ALL cells.	С

Parameter	Screen	Description	Туре
P18(AZERO)	IO	Programmable input: can be configured that the signal of cell 1, 2 or 3 is zeroed when active, or ALL cells.	С
P21(START)	IO	Programmable input: can be configured that the data-ac- quisition on sensor board 1, 2 or 3 is started when active, or on ALL sensor boards.	С
Range	dc setup, dc stat, prog, pulse setup1, pulse stat, scan setup, scan stat	Range setting, varying from 10 pA to 200 μA full scale, in 1, 2 and 5 steps. In the pulse and scan mode 10 nA to 200 μA full scale can be used.	С
S	scan setup	Scan speed, can be set from 1 - 100 mV/s in 1, 2, 5 steps.	С
SPD	scan stat	Scan speed, can be set from 1 - 100 mV/s in 1, 2, 5 steps.	С
START	run, scan stat	In DC and pulse mode: toggle between STOP and START execution of a time file. Starts a scan in scan mode.	F
STOP	run, scan stat	Scan mode: STOP aborts scan and resets cell potential to E1.	F
		DC and pulse mode: toggle between STOP and START to control execution of a time file. Pressing 'STOP' aborts this run, cycle counter (Cy) is reset to 1. STOP also deactivates the outputs Aux 1 and 2, and Relays 1 and 2 (status: 0000) and sets the electric valve to load (if present).	
SYSTEM	diag	Enter SYSTEM menu	F
t	pulse setup2, pulse stat	Displays the total duration of one pulse (t1 + t2 + t3 + t4 + t5).	S
t1, t2, t3, t4, t5	pulse setup2 pulse setup3	Duration of potential step E1, E2, E3, E4 and E5. Time can be set between 0 (t2 - t5) or 100 (t1) and 2000 ms in 10 ms increments. Maximum pulse duration is 9999 ms.	С
Temp	config	Controls the temperature of the oven. Range: off, 15 - 60°C, selectable in 1°C steps. The oven is stable from 7°C above ambient oven temperature.	С
Toven	dc setup, pul- se setup1	Controls the temperature of the oven. Range: off, 15 - 60°C, selectable in 1°C steps. The oven is stable from 7 °C above ambient.	С
ts	pulse setup2	Controls the duration of the sampling time in the pulse mode. The time can be set between 20 ms and maximum t1-60ms with 20 ms increments.	С
Tsensor	system	Displays active temperature sensor	S
Valve	config	User confirmation whether a manual valve is connected to phone jack C on rear panel. If present: INJ=I or INJ=L appears in DC/Pulse Status screen	S
Vout	stat (dc, pulse, scan), run, noise	Displays output signal.	S
Vout source	config	Sets the output source from the analog data output: DAC (processed digital signal after 16-bit AD conversion) or I/E (true analog signal from the I/E converter)	S

5.4 ClarityChrom[®] ECD 2.1 control module

Full control and data acquisition of the DC and Pulse measurement mode for both the ECD 2.1 is supported in the ClarityChrom[®] chromatography software.

lew Oper			• as	Report se	tup A	udit trail	Send method by e-mail	() Help	About 🖂 💷
ct Detector			Detec	tor 1		•] [✓ Enabled		
				DECADE	lite Detec	or Methor	d	Det Status	
Main Ou	put	Time Table Mode	Pulse		*				
Sensor			Puse		-				ClarityChrom®
	Rate	1/pulse du	ration	• 1,59 [H2]				
	E1	0,1	[V]	t1	0,4	[2]			a second s
	E2	-2	M	12	0,02	[2]			
	63	0,6	M	13	0,01	[2]	ts 200 - (ms)		
	E4	-0,1	M	14	0,2	[2]			
	ES	0	M	15	0	[9]			ClarityChrom version 8.1.0.87 Krauer GribH, 2019 Build - Mar 28 2019
									S/N - 015-061340 http://www.knauer.net/klaritychrom
									ClantyChrom What's New Readme System Files
									OK
et Status	1	Not Ready (C	æll off, Þ	lethod has r	ot been s	int)		From Det	
ent Table	AS	LC Gradien	t LC	Manager			Thermostat Integration	Colordation Advanced	
ens rable	AD.	LC Gradien	t te	PROBABILIT	ACC	uisition	inemosiat integration	Cachedon Advanced	20.0
ок	Cane	el						Sgnd Me	thod
		***!**	_	-		-	-		
	-	- ~ ~	~ 4						
16	· F	-(.)	11	me	tho	d w	undow in	(Jarity(hro	m®. On the main tab the m

Within a ClarityChrom[®] method it is possible to execute timed event by programming time lines under the Time table tab. Running a Time table enables time-based, automated parametric control of the electrochemical detection (ECD 2.1) during a run. This is particularly useful when during a run settings have to be changed such as the sensitivity, auto zero or control of external equipment.



The SCAN mode is not supported in ClarityChrom[®]. Scanning Voltammetry is supported in the Elite Dialogue software.

5.5 Dialogue Elite

Dialogue Elite for Windows is a multi-functional program to control the ECD 2.1.

File Tools Monitor Deteo Cell	Options Events Sequence Window or Sequence Events Log 1 V On V Enabled	Help Mode DC v Dev status
Output Range Offset Filter D Data refe Polarity	1 v	Potential E 0.80 M
Temperatu Oven Set Measured	V On 35 (C) 363 (C)	Run time t 15.000 [min] In/Output Relay TTL contact 1 2 1 2 3 4
Read from device	I IP: 192.168.0.135 🔲 send/recieve	Send to device
main meas		Dialogue Elite. On the detector tab the ns can be set/controlled (measurement

Dialogue Elite is a tool for:

- controlling all operational parameters
- programming timed events in one run or sequence
- scanning voltammetry
- firmware upgrades using the FW upgrade wizard
- OQ, PQ and dummy cell measurements
- device calibration

There are no tools for further data analysis, therefore Dialogue cannot be considered as fully featured chromatography software package. The free demo version of this program has already all functionality implemented, functionality is unlocked by a license dongle.

The free version is primarily used for controlling all operational parameters of the detector. The free version cannot be used for starting a run, running a calibration or OQ.

For details about the use and functionality of the Dialogue Elite software please refer to document 175.0015 Dialogue Elite user instructions which can be downloaded from the Antec website (<u>www.myantec.com</u>).

6. Detection and Parameters


6.1 Introduction

One of the characteristics of electrochemical detection is its tremendous dynamic range. In amperometric detection peak heights may vary from micro-amperes down to the pico-ampere range. The ECD 2.1 covers such a wide range from 200 μ A down to 10 pA full scale, without being limited by electronic noise. For this reason the ECD 2.1 is equipped with a 24-bit ADC and 16-bit DAC for analogue data output. One of the key features is that data can be sampled with data collection rates up to 100 Hz (100 pts/ sec) in DC mode which assures that fast peak responses typical for UHPLC can be detected with sufficient resolution.

6.2 Three-electrode configuration

The circuitry of the ECD 2.1 detector is designed for operation with electrochemical flow cells with a three-electrode configuration (see Fig. 18). The working potential is set between the working electrode (WE) and the auxiliary electrode (AUX). The AUX is kept at a precisely defined reference electrode (REF) potential by means of the so-called voltage clamp. This is an electronic feedback circuit that compensates for polarization effects at the electrodes.

At the WE, which is kept at virtual ground, the electrochemical reaction takes place, i.e. electrons are transferred at the WE. This results in an electrical current to the I/E converter, which is a special type of operational amplifier. The output voltage of the I/E converter is digitized in the instrument by means of a 24-bit A/D converter and processed, and the resulting output current Ic can be acquired digitally by PC control software (Elite Dialogue or ClarityChrom®) or analog using the 'Analog Data output' on the rear panel connected to a recorder or an external A/D converter.

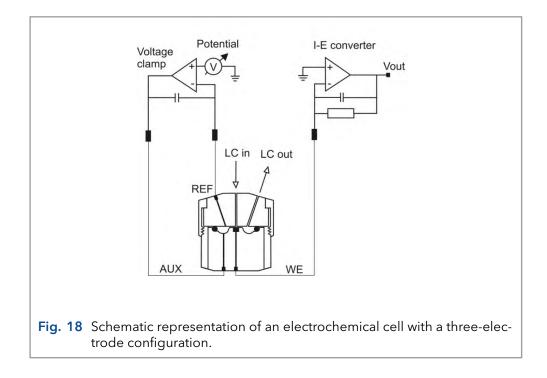


Essentially, for the oxidation or reduction reaction it would be sufficient to use only two electrodes. However, the three-electrode configuration has several advantages over a two-electrode configuration. If the working potential would be applied only over an AUX versus the WE (without REF), the working potential would continuously change due to polarization effects at the electrodes, resulting in highly unstable working conditions.

If the working potential would be applied only over the REF versus the WE (without AUX), the working potential would be very well defined. However, the potential of a REF is only well defined if the current drawn is extremely low (pico-amperes) resulting in a very limited dynamic range.

A three-electrode configuration, combines the best of both electrodes. The REF stabilizes the working potential and the AUX can supply high currents. This results in the tremendous dynamic range of a three-electrode system.

6.3 Internal organization

At the working electrode (WE) in the electrochemical flow cell the electron transfer takes place due to an oxidation or reduction reaction. The resulting electrical current is amplified by the current-potential (I/E) converter (Fig. 19).

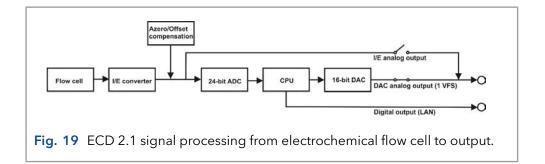


The signal from the I/E converter can be compensated with auto zero or offset, and is digitized using a 24-bits ADC. In the CPU the signal is processed, for example noise filtering, or more complex data processing in PAD. Finally after the 16-bits DAC the signal is set to a 1 V full scale analog data output (by default Output = ADC). Also the true analog signal from the I/E converter (before AD conversion) is available via the 'Analog data output' connector. This output can be selected in the CONFIG menu by setting the parameter Output = I/E.

6.4 Dual flow cell control

The ECD 2.1 electronics are located on 2 different PCB's (printed circuit boards). The control board and the sensor board. The control board is dedicated to communication with PC (LAN) and keyboard & display. It has a processor with a so called ,event handler' that takes care of all user commands and hardware interrupts. The sensor board is fully dedicated to data acquisition and flow cell control. By using this architecture it is possible to extend the functionality of the detector to more than one flow cell by simply adding a sensor board. The control board and other hardware is prepared for more than one sensor board. Typically, a two flow cell configuration can be used in serial or parallel mode detection.


Serial mode detection


In serial mode one LC system is used, with 2 flow cells in series. For data acquisition 2 data channels are applied with the same time base. Serial mode detection is especially suitable for OX-RED or RED-OX applications, examples are analysis of vitamin K and nitro-tyrosine, using micro HPLC. The first flow cell is a FlexCell that converts the analyte of interest in a detectable substance. The second flow cell is a SenCell or VT-03 cell which is used for detection. Note that it is necessary to work with micro HPLC because the conversion rate of the reactor cell is too small when using standard HPLC.

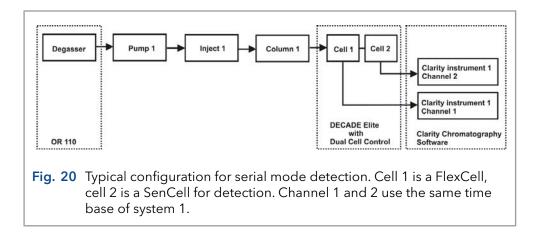

Parallel mode detection


In parallel mode 2 HPLC systems are used with 2 flow cells. In fact, the ECD 2.1 is operated as if 2 independent detectors are in one housing.

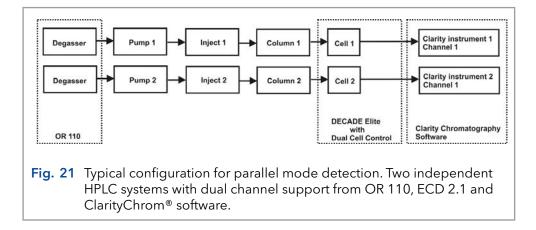

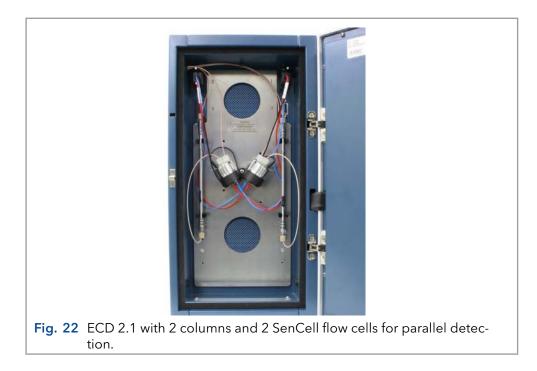


Navigation in dual cell menu

All menus for a dual flow cell system are similar to a single cell system with 2 exceptions. First, in the top right corner a number is visible which indicates the active cell in display. Toggle with the "+" and "-" buttons between sensor boards. If the board number does not change it means that the second sensor board is not installed or not properly recognized. Second, a new status screen is available in dual cell systems which indicates the status of both cells in a single screen. However, for convenience it is advisable to use PC control from the ClarityChrom[®] Chromatography software or Dialogue Elite software when working with 2 flow cells.

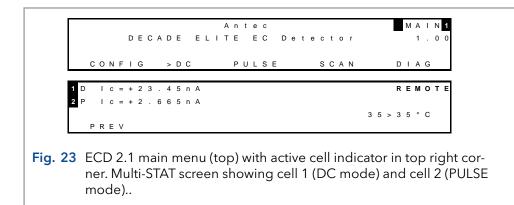

6.5 Parameters

Operational parameters are controlled from the SETUP screens in the ECD 2.1. Parameters are filter, cell potential and offset. Temperature is set in CONFIG menu.

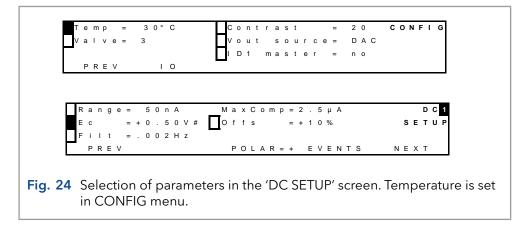

Range

Range selection is done in the 'SETUP' or 'STAT' screen in DC, PULSE and SCAN mode. A number of ranges can be selected; the maximum current that can be compensated for using auto zero and offset differs. The high sensitivity ranges (10 pA - 5 nA) have the best noise specifications. In fact, there is a trade-off between best noise specification at sensitive ranges, and maximum compensation at the less sensitive ranges. This is an inevitable consequence of the tremendous dynamic range that is covered by electrochemical detection.

Range FS	Max comp	Range FS	Max comp
200 µA	2.5 mA	20 nA	2.5 μΑ
100 μA	2.5 mA	10 nA	2.5 μΑ
50 μΑ	2.5 mA	5 nA	250 nA
20 µA	2.5 mA	2 nA	250 nA
10 μA	2.5 mA	1 nA	250 nA
5 μΑ	2.5 mA	500 pA	250 nA
2 μΑ	25 μΑ	200 pA	250 nA
1 μΑ	25 μΑ	100 pA	25 nA
500 nA	25 μΑ	50 pA	25 nA
200 nA	25 μΑ	20 pA	25 nA
100 nA	25 μΑ	10 pA	25 nA
50 nA	2.5 μΑ		

Table II. DC ranges and maximum compensation:

In the PULSE and SCAN mode, current is much higher than in DC mode. Therefore it is not possible to select pA ranges.

Range FS	Max comp	Range FS	Max comp
200 µA	2.5 mA	500 nA	25 μΑ
100 µA	2.5 mA	200 nA	25 μΑ
50 μΑ	2.5 mA	100 nA	25 μΑ
20 μΑ	2.5 mA	50 nA	2.5 μΑ
10 µA	2.5 mA	20 nA	2.5 μΑ
5 μΑ	2.5 mA	10 nA	2.5 μΑ
2 μΑ	25 μΑ		
1 μΑ	25 μΑ		

Table III. PAD ranges and maximum compensation:

Offset


A maximum offset of +50% and - 50% in 5% steps can be set. For example, 20% is a 200 mV offset when the maximum output is 1.0 Volt.

Polarity

The polarity of the output can be inversed. Oxidative and reductive analyses generate opposite currents. For data acquisition, traditionally chromatographic peaks have a positive amplitude. Therefore selection of polarity is useful.

Filter

High frequency noise is efficiently removed and chromatographic peaks can be detected with better signal to noise ratio.

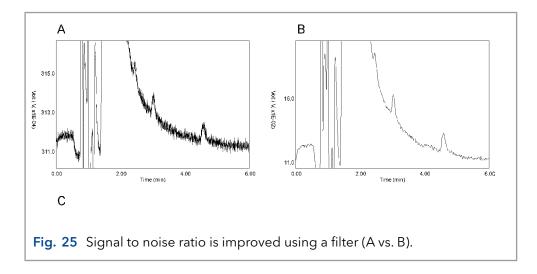

The ECD 2.1 is equipped with ADF (Advanced Digital Filter) as a tool to filter the acquired signal and improve the sensitivity of the analysis (Signal-to-noise ratio). In the next chapter the filter setting is explained including detailed background information about filtering.

DC mode

In the tables below the available filter settings for the DC mode are listed with the corresponding data rate of the output. Data rate is expressed as number of data points per second (Hz). In the DC mode the data rate is not an adjustable parameter but is coupled to the filter setting, except for RAW. RAW is special, the incoming data are not filtered and a data rate between 1 and 100 Hz can be selected.

Table IV. DC mode filter setting and corresponding data rate:

Filter setting DC mode (Hz)	Data rate (Hz)
RAW	100(default), 50, 20, 10, 5, 2, 1
10	100
5	50
2, 1	20
OFF	10
0.5, 0.2, 0.1	10
0.05	5
0.02	2
0.01, 0.005, 0.002, 0.001	1

Filter OFF is also a special case. The data rate is fixed to 10 Hz, and the data is not filtered. Setting OFF is therefore the same as RAW at 10 Hz.

Pulse mode

In the pulse mode the working electrode is dynamically and continuously regenerated by a series of potential steps in a cyclic manner. Data is processed differently, and the data rate is defined by the total duration of the 5 potential steps in a pulse: t1 + t2 + t3 + t4 + t5. Usually the typical pulse duration is between 0.5 and 2s (data rate between 2 – 0.5 Hz). Filter settings in the pulse mode are therefore selectable between 0.5 and 0.001 Hz, and OFF.

The scan mode has no filter at all.

7. Noise suppression: ADF™

7.1 Introduction

Besides for its tremendous linear dynamic range and selectivity, electro-chemical detection is well-known for its very low limits of detection. To further improve these detection limits the manufacturer's engineers have developed ADF (Advanced Digital Filter) and the ECD 2.1 has been equipped with it as a standard. The improvement factor in signal-to-noise (S/N) ratio depends on the frequency relation of signal and baseline noise. S/N improvements from a factor 5 up to more than 100 have been obtained.


To understand how a digital filter works, first the importance of frequencies in chromatographic analysis will be explained. Then we will look at peak width, filter settings, cut off frequency, amplitude response plots and finally at a few chromatograms before and after applying ADF.


7.2 Frequency

A scientific definition of frequency is "the number of completed alterations per unit time". It has two dimensions: count and time. Frequency is usually expressed in Hz, which is counts per second.

The counts themselves can run in a regular, evenly spaced manner, as with sine waves whose curve shapes do not change. Alternatively, the counts can run in an irregular manner within the specified unit of time. If the latter happens, frequencies would vary if broken down into smaller units of time.


In the example of Fig. 26 a signal is shown with a frequency of 12 alterations in 5 minutes. To express its frequency in a more scientific way a full period is precisely determined and expressed in Hertz (or s-1). It is a sine wave with a frequency of 0.04 Hz (Fig. 27).

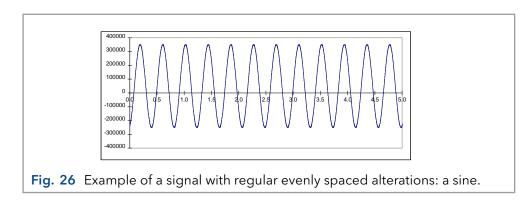

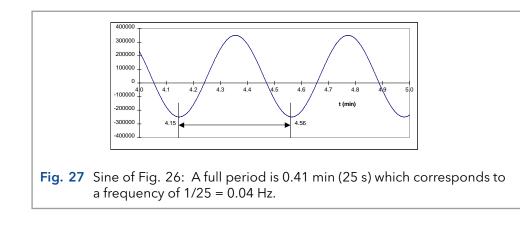


7.3 Frequency of signal and noise


Also a chromatographic peak can be expressed in terms of frequencies. The way to determine this frequency is the same. The duration of the full peak is measured and expressed in Hz.

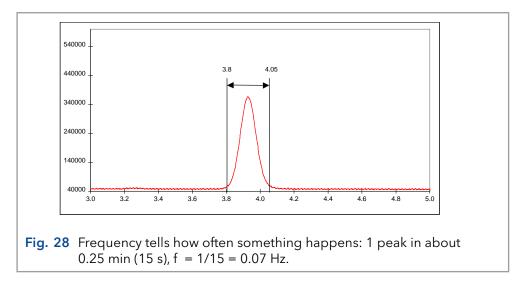


This is further illustrated by an overlay of the same chromatographic peak with a sine of 0.07 Hz (Fig. 29).

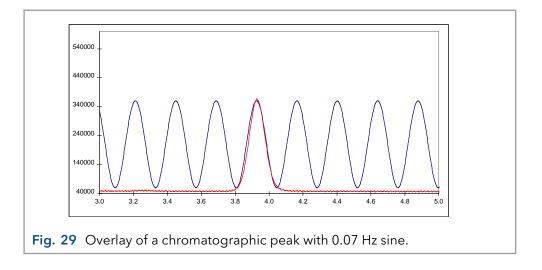

Typically in chromatography narrow peaks are in front of a chromatogram while peaks with longer retention times get wider. As a consequence frequencies are not constant but vary between 0.1 - 0.01 Hz, which corresponds to 10 - 100 s peak width.

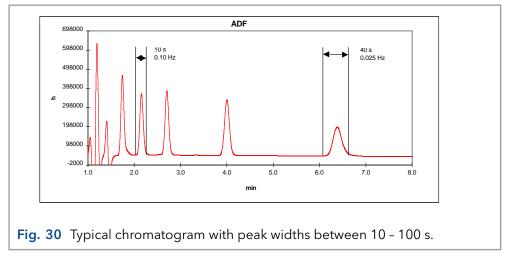

Noise in chromatography can come from different sources. Pump pulsations are typically shown as a very regular noise pattern, while electronic noise has a more random character. This is illustrated in Fig. 31 where a noise trace is shown with an overlay of a 10 and 0.4 Hz sine.

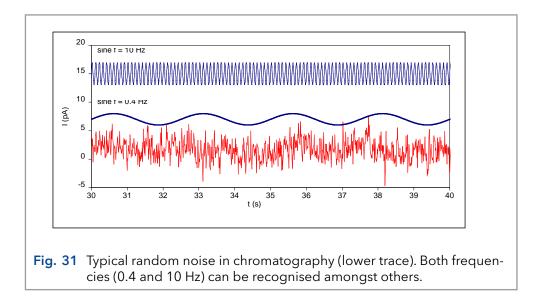

Looking closely to the lower noise trace both frequencies (and others) can be recognised. This is typical for noise in chromatography: a collection of more or less random frequencies.

Low pass noise filters

The way noise filters work is by suppressing certain frequencies in the acquired signal. Typically low pass filters allow chromatographic peaks (low frequency) to pass, while high(er) frequency noise is attenuated. No matter how advanced, it is impossible to use a low pass filter successfully if there is no difference in frequency of signal and noise.

Analogue filters are made of hardware, from capacitors, resistors and amplifiers (opamps). Digital filters are mathematical routines to process an acquired signal.

Traditionally, in many detectors for chromatography an analogue lowpass filter is applied (rise time filter). A 'passive' RC filter consists of resistors and capacitors. An active higher order filter can be considered as a number of these RC filters in series. In a 4th order filter the signal coming from the first filter is filtered again in a second, third and fourth filter. During these steps, loss of signal occurs simply because of all the resistors that are applied. Operational amplifiers, which are 'active' components, are applied in each stage to restore the signal to its original value.

With the availability of powerful processors, digital signal processing has become an excellent alternative for hardware filters. In its most simple form a running average filter takes the average of n data points to create a new data point. For example in a 5-points running average filter output data point y[80] is calculated from measured data points x[80] - x[84] as:

$$y[80] = \frac{x[80] + x[81] + x[82] + x[83] + x[84]}{5}$$

Each input data point has the same weighting factor of 1/5. In more advanced digital signal processing a more complicated equation is used to calculate the output data point y[n]:


$$y[n] = a_0 x[n] + a_1 x[n-1] + a_2 x[n-2] + a_3 x[n-3] + \cdots$$

In contrast to the previous equation, each data point has a different weighting factor a. Sum of these weighting factors a0...n will always be 1.

Characteristic of noise filters is that processing the signal will result in a delay. This is inevitable, as the mathematics of digital signal processing requires a number of previous data points to process a new data point.


The filter characteristic in DSP is often named after the scientist who 'invented' the mathematics behind the signal processing routine. Wellknown names in this field are Bessel, Chebychev, Savitsky, Golay, Hamming and many others.

7.4 Amplitude response plot

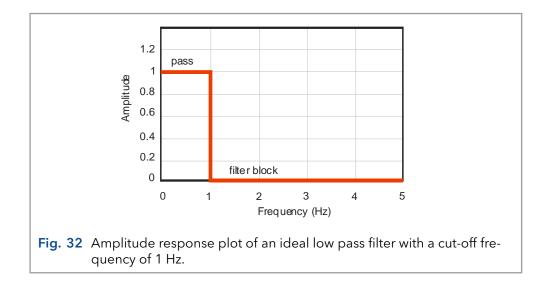


There are several ways to describe the filter characteristics. An amplitude response plot gives important information on filter behavior. Suppose our signal of interest has a frequency between 0 - 1 Hz, and all higher frequencies are noise. An ideal filter is shown in Fig. 32 where signal frequencies between 0 - 1 Hz completely pass while frequencies of higher than 1 Hz are completely blocked.

In practice filters behave a bit different from the ideal situation. Amplitude response plot shows a more gradual attenuation profile at higher frequency. This cut off frequency is where the output signal amplitude is 70% of the input signal, also known as 3 dB point.

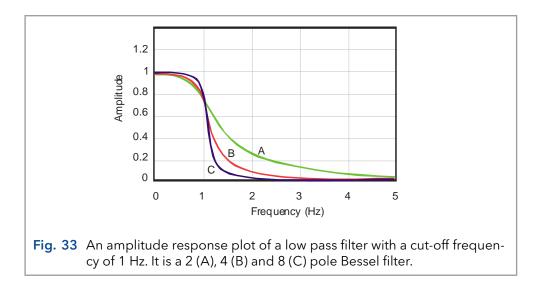

In Fig. 33 it is shown that the number of poles is important, a filter behaves more ideal with increasing number of poles. In a hardware filter the number of poles is the number of filter circuits that are placed in series.

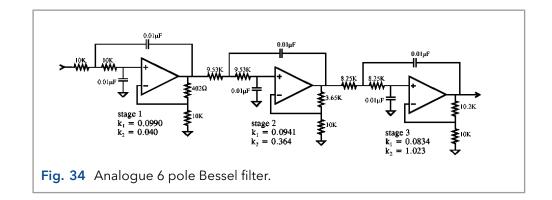


A digital filter does not have poles, but it is characterized by the number of input data points used to calculate a new output data point. For example a 9-point digital filter (Savitzky-Golay) is given as:

```
Y[1] = -0.090909091 X[1] + 0.0606060601 X[2] +
0.168831169 X[3] + 0.233766234 X[4] +
0.255411255 X[5] + 0.233766234 X[6] +
0.168831169 X[7] + 0.060606061 X[8] +
-0.090909091 X[9]
```

Note that the sum of coefficients is exactly 1. Y[n] is the output data point, X[n] are input data points. Generally spoken, the performance of a digital filter improves with more input data points, but also more processor capacity is required for the large number of calculations.

7.5 Applying ADF in chromatography

If noise frequencies in LC-EC differ from the frequency of the signal, noise can be suppressed. Using the right filter setting (cut-off frequency) will specifically attenuate noise and improve the signal-to-noise (S/N) ratio. No matter how 'advanced' a filter is, it is only possible to apply low pass filtering if noise frequencies are higher than the frequency of the signal.

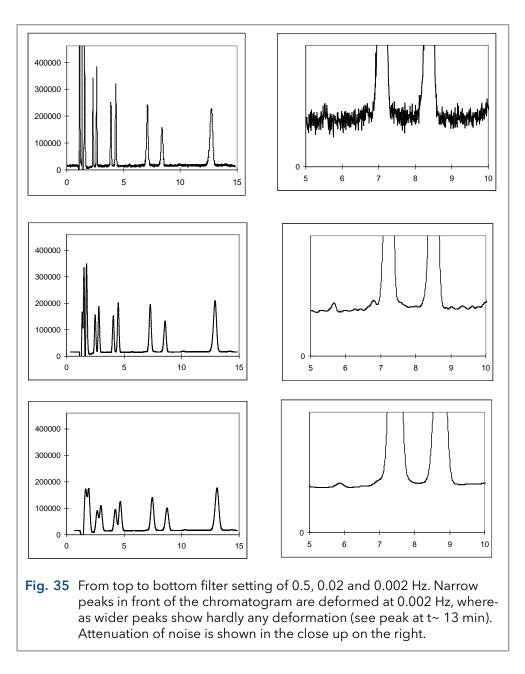


Prerequisite for a 'good' noise filter for data acquisition in liquid chromatography is that it improves the S/N ratio without significant distortion of the signal of interest. This is particularly difficult if the frequency of the signal is close to the frequency of the noise. The ECD 2.1 has a number of filter settings to optimize for best possible signal-to-noise ratio. The width of the peaks of interest is important because wider peaks allow stronger filter settings simply because of the lower frequency of such peaks. Advised filter setting to start further optimization is given as:

Filter setting = 1 / [2 * (peak width)]

So at a 10 s peak width a 0.05 Hz filter setting is advised. If peaks are 50 s a 0.01 Hz filter is advised to start with. Note that if a chromatogram has interesting peaks of 10 s as well as 50 s, it may not be possible to work with one filter setting. In that case it is advisable to switch to a stronger filter setting for the second half of the chromatogram using a timed event. To optimize for the best S/N ratio, use the lowest acceptable cut-off frequency.

After optimization, do not change the cut-off frequency setting during analysis of a calibration sequence. Use the same settings for analysis of samples and calibration standards.

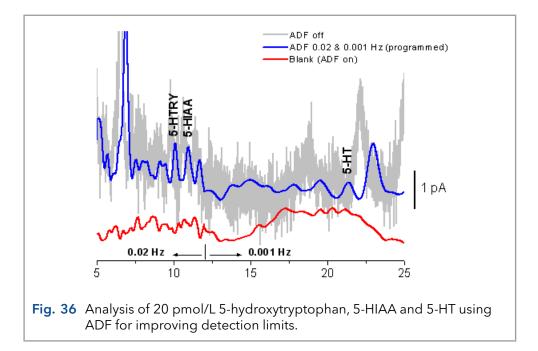


The S/N improvement depends on the composition of the frequency spectrum. Improvement up to a factor 100 may be obtained compared to an unfiltered signal. As high frequency noise is suppressed, remaining noise components will be in the same frequency range as chromatographic peaks.

As suppressing noise will always result in (some) suppression of signal it is advised to switch the ECD 2.1 to the highest acceptable sensitivity.

8. Pulsed Amperometric Detection (PAD)

8.1 Introduction

The ECD 2.1 can operate in PAD mode. In pulsed amperometric detection (PAD) the working electrode (WE) is regenerated at a frequency of 0.5 - 3 Hz by the application of a series of potential changes. This is particularly useful for certain applications where the working electrode is rapidly fouled due to adsorption of insoluble reaction products. A well-known application area of PAD is the analysis of carbohydrates (Fig. 37) [1].

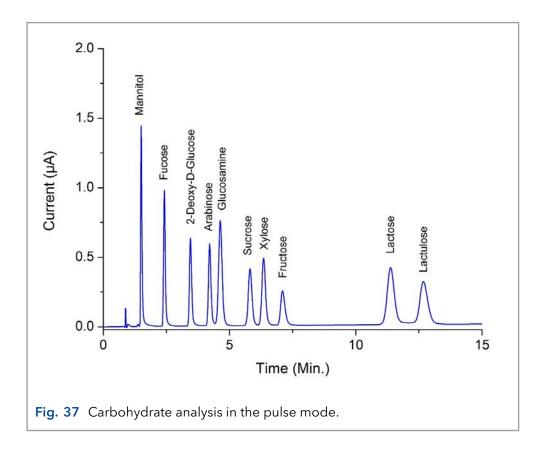

Pulse vs. DC mode

The pulse mode is quite different from the DC mode. Instead of a constant potential, a series of potential steps is applied in a cyclic manner. The signal is sampled during a fraction of the total pulse cycle. During the sampling time (ts) the signal is collected and this value is sent to the detector output. The frequency of data output is determined by the pulse duration: t1 + t2 + t3 + t4 + t5. Usually the duration is between 0.5 and 2s (data rate between 2 - 0.5 Hz).

The background or cell current is usually considerably higher (100 - 1000 nA) than in the DC mode. Therefore, only nano- and micro-ampere ranges are available in the pulse mode. Typically, the background current is between 0.1 - 2 μ A.

When the frequency of the data acquisition system (integrator) is higher than the pulse frequency a typical stepwise pattern may appear in the chromatogram. This is called 'oversampling' and these steps are usually only visible after considerable magnification of the chromatogram. It is recommended to keep the data acquisition at 1 Hz.

High pH of mobile phase

In carbohydrate analysis another 'special' consideration has to be taken into account. Detection is done in the pulse mode using a flow cell with gold electrode. For separation an anion exchange column is used with a mobile phase of around 20-100 mmol/L sodium hydroxide. The pH is between 12 and 14. This puts some demands on the HPLC system, and flow cells.

After prolonged use of the flow cell with a gold working electrode (WE) in the pulse mode, the gold oxide which is generated at the WE, precipitates on the auxiliary electrode (AUX). This gold oxide coating may electrically isolate the AUX and result in an increase of the noise. Cleaning the AUX electrode with metal wool is a way to remove this coating.

The reference electrodes of the Ag/AgCl type are not suitable for carbohydrate analysis. Due to silver oxide formation they require regular (monthly) maintenance. HyREF reference electrodes are maintenance free under these conditions and are therefore particularly suited.


If a mobile phase is used with a high pH (pH>10, carbohydrate analysis), the standard Vespel rotors from the injection valve should be replaced by Tefzel rotors which are pH resistant.

For carbohydrate analysis, only CO2-free sodium hydroxide should be used since carbonate anions may disturb the ion exchange chromatography. It forms CO32- which is a strong modifier. The CO2-free sodium hydroxide is available from several suppliers as a 50% solution (19.2 mol/L). NaOH pellets are not recommended because of their high CO2 content.

Organic modifiers (acetonitrile) strongly attenuate the signal of most carbohydrates in PAD and are therefore not recommended.

Pulse settings

In PAD of carbohydrates a series of potentials is applied in a continuous cyclic manner. During time interval t1 the detection potential is applied. The data collection occurs within t1, during time interval ts (sampling time). The time difference t1 - ts is the stabilization time.

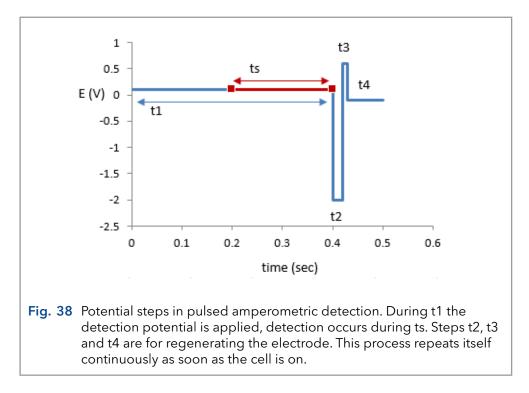

During the next time intervals (t2..t4) the electrode is 'cleaned' by reductive and oxidative potential steps.


Optimization of wave forms

LaCourse and Johnson [2-4] have published several papers on optimization of wave forms in PAD. Several considerations are important for the choice of the pulse duration. Optimization is depending on the working electrode material, the sample constituents and the required detection frequency. The impression may arise that the number of variables, 5 potential steps and 6 time settings, may lead to a time-consuming optimization procedure. In practice, the pulse mode is more straightforward and published in several excellent review papers and application notes.

Output frequency

An important difference between the DC and the pulse mode is the frequency of the output signal. In the DC mode the signal has a 1-100 Hz frequency, in the pulse mode the frequency is determined by the duration of the pulse. Once every cycle, the ts signal is sent to the output.

If a stepwise pattern in the chromatogram is seen when zooming in, it means that the data acquisition has an unnecessarily high sampling frequency. This leads to large data files, but certainly not to a better chromatogram. Usually, data acquisition at 1 Hz is sufficient.

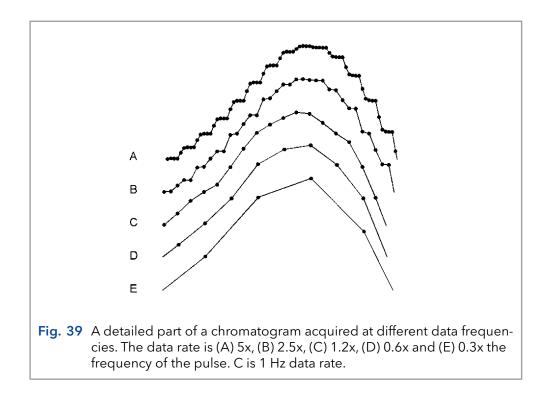

Working electrode material

Gold and platinum are used as working electrodes for PAD. Glassy carbon appears to be unsuitable due to the high electric capacitance of this material. Furthermore, resurfacing of the noble metal working electrode is based on formation and removal of a (metal) oxide layer. This is impossible with glassy carbon.

References

- 1. D.C. Johnson, D. Dobberpuhl, R. Roberts and P. Vandeberg, Review. Pulsed amperometric detection of carbohydrates, amines and sulphur species in ion chromatography - the current state of research, J. Chromatogr. 640 (1993) 79-96
- **2.** D.C. Johnson en W.R. LaCourse, LC with pulsed ECD 2.1 at gold and platinum electrodes, Anal. Chem., 62 (1990) 589A-597A
- **3.** W.R. LaCourse en D.C. Johnson, Optimization of waveforms for pulsed amperometric detection of carbohydrates following separation by LC, Carbohydrate Research, 215 (1991) 159-178
- **4.** W.R. LaCourse en D.C. Johnson, Optimization of waveforms for pulsed amperometric detection of carbohydrates based on pulsed voltammetry, Anal. Chem. 65 (1993) 50-55

9. Optimization of the working potential

9.1 Introduction

A current - voltage (I/E) relationship, or voltammogram, characterizes an analyte. It gives information on the optimum working potential, which can be used to improve detection sensitivity and selectivity.


There are several ways to obtain a voltammogram:

- A hydrodynamic voltammogram is obtained in the DC mode by running several chromatograms at different working potentials. Both peak height and background current are plotted against the working potential. A hydrodynamic voltammogram has as an advantage that the I/E relationship of all analytes of interest can be obtained simultaneously in one set of experiments (boundary condition: all analytes should be sufficiently separated under the applied LC conditions). Furthermore, under real chromatographic conditions reliable information about the S/N ratio is obtained.
- A scanning voltammogram is obtained in the so-called scan mode of the ECD 2.1: the voltage runs between two preset potential values (E1 and E2) and scan speed (in mV/s) and the current is measured.

Optimization of the working potential and the construction of a voltammogram is described.

9.2 Electrochemical reactions

In electrochemical detection (ECD) a reaction of the analyte at an electrode surface is monitored. This distinguishes ECD from most other detection techniques where detection is based on the physical properties of an analyte (i.e. mass spectrometry: molecular mass, absorbance detection: molar absorptivity). For electrochemically active compounds, the potential between reference electrode (REF) and working electrode (WE) determines the reactivity of the analyte at the WE. The potential difference supplies the energy level needed to initiate or enhance the electrochemical reaction. Different analytes may have different oxidation or reduction potentials, which determines the selectivity of ECD.

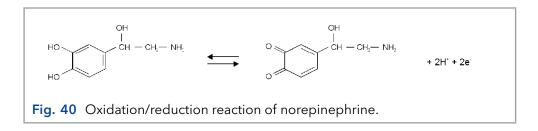


An example of an electrochemical reaction is shown in Fig. 40, norepinephrine is converted into a quinone by oxidation at the WE. Two electrons are transferred at the WE resulting in an electrical current that is amplified by the controller.

9.3 Hydrodynamic and scanning voltammogram


Hydrodynamic voltammogram

A hydrodynamic voltammogram is constructed when the pure analyte is not available and separation over an analytical column is required. Furthermore, under real chromatographic conditions reliable information about the S/N ratio is obtained. The peak heights obtained from the sequence of chromatograms are plotted against the working potential used. Also the background current (Icell) is plotted.

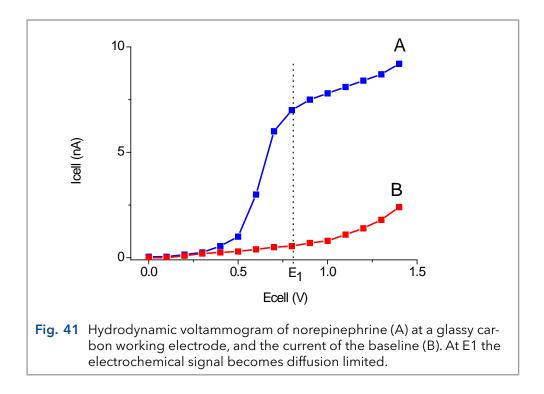

Scanning voltammogram

An alternative for the chromatographic construction of an I/E relationship is the application of scanning voltammetry. In a scanning voltammetry experiment the working electrode potential is ramped up and down between two preset potentials (E1 and E2) and the current is measured while the analyte is continuously flushed through the flow cell. This is repeated as many times as desired. The rate of voltage change over time is defined as the scan rate (mV/s).

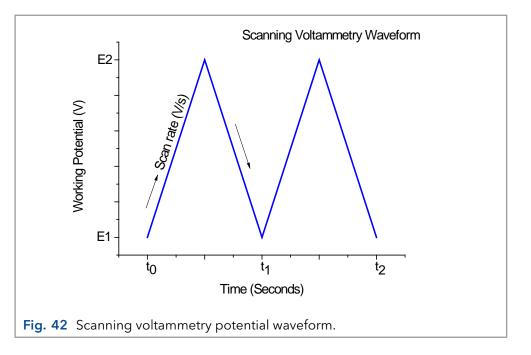


The current is plotted against the working potential to give a voltammogram (I/E curve). An example is shown in Fig. 43. A difference with a hydrodynamic in scanning voltammetry is that no HPLC separation is involved. The signal is the sum of all EC active substances. It takes only a few minutes to construct a scanning voltammogram. This is an advantage, especially when a number of analytes have to be characterized. However, it is a prerequisite to have the pure analyte dissolved in buffer. A scan of the buffer (blank) should be used to distinguish between solvent peaks and analyte peaks.



Any contamination in the buffer may lead to artefacts.

As can be seen in both Fig. 41 and Fig. 43, when the working potential is increased the electrochemical reaction is enhanced hence the signal increases. At a certain potential the I/E curve flattens. All analyte molecules that reach the working electrode are converted at such a high rate that the analyte supply becomes the limiting factor. At the working electrode surface a stagnant double layer exists, where molecular transport takes place by diffusion only. Therefore, the current at (and beyond) this potential is called the diffusion limited current.

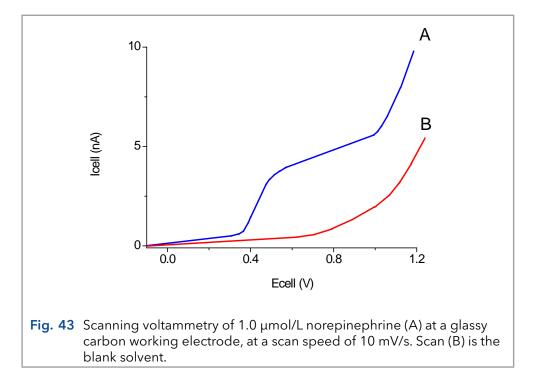


With respect to sensitivity, a high working potential is important. However, at higher working potentials, more analytes are detectable. So, as to selectivity, a low working potential will be favorable.

Working at a potential on the slope of the I/E curve will result in less reproducibility in HPLC. A small fluctuation in the applied potential, or any change in the system (like for instance a pH change) may result in differences in measured peak height. In practice the choice of the working potential is a compromise between sensitivity, selectivity and reproducibility. In the example of Fig. 41 a working potential (E1) of 0.8 V is chosen.


9.4 Optimization using a voltammogram

Sometimes, when interfering peaks appear in the chromatogram, it is possible to optimize the method with regard to selectivity. If the interfering compound has a higher oxidation potential, a working potential is chosen that gives the best selectivity, i.e. the largest difference in peak height. In the example of Fig. 44 the selectivity for compound X is improved considerably by decreasing the potential to E2 or E1. Obviously, if compound Y is the compound of interest, optimization of selectivity in this way is not possible and the chromatography has to be optimized.

Electrochemical detection differs from most other LC detection methods in that a reaction takes place in the detection cell. Due to reaction kinetics an increased temperature speeds up the oxidation/reduction reaction. However, this not only holds for the analyte but also for the background current and possible interferences. An elevated temperature will therefore not automatically lead to a better detection. A constant temperature is of paramount importance for a stable baseline and reproducible detection conditions.

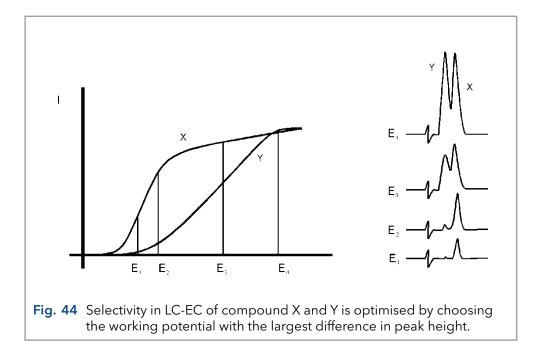

Electrochemical reactions are pH sensitive (Fig. 45). For norepinephrine the I/E curve is shifted to a lower potential at higher pH. When the working potential is high (E2), and the signal is diffusion limited, an increase in pH will result only in a small increase of the peak height. When the working potential is lower (E1), and the signal is not diffusion limited, the signal will strongly increase at higher pH. In both cases the background current increases at a higher pH.

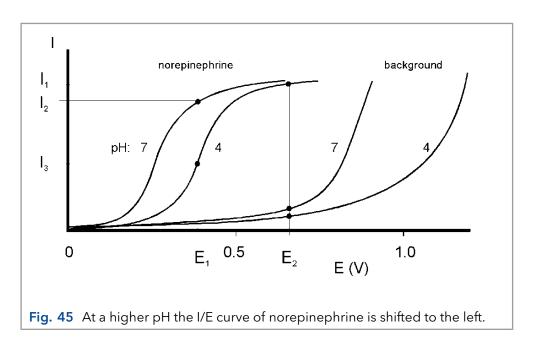


Reaction kinetics predict that electrochemical detection is mass flow dependent. When the LC flow is stopped in LC-EC, the analyte will be oxidized completely and the signal decreases rapidly. This means that the flow rate not only affects temporal peak width and analysis time but also peak height. Also the background signal is sensitive towards fluctuations in the flow rate. Therefore, it is important to use a pulse-free solvent delivery system.

9.5 Construction of a hydrodynamic voltammogram

Before a hydrodynamic voltammogram can be obtained, the chromatographic conditions should be optimized. Then the following steps are taken:


- 1. A solution of the analyte at a concentration between 1 100 $\mu mol/L,$ is prepared in mobile phase.
- **2.** The electrochemical detector is stabilized in the DC mode at a high potential. After stabilization the background current is read from the display of the detector (Icell) and the noise is measured.
- **3.** The run is started by injecting the compound. When at the high working potential no signal is obtained, it may be concluded that the compound is not electrochemically active. In such a case derivatization of the compound may be an option.
- **4.** If a peak is measured, the working potential is decreased by 50 or 100 mV and step 2 to 4 is repeated until the lowest potential setting (Fig. 46).
- **5.** The peak heights and the background currents are plotted against the working potential (Fig. 41).

The working potential which gives the best sensitivity is obtained by plotting the signal-to-noise ratio against the working potential.

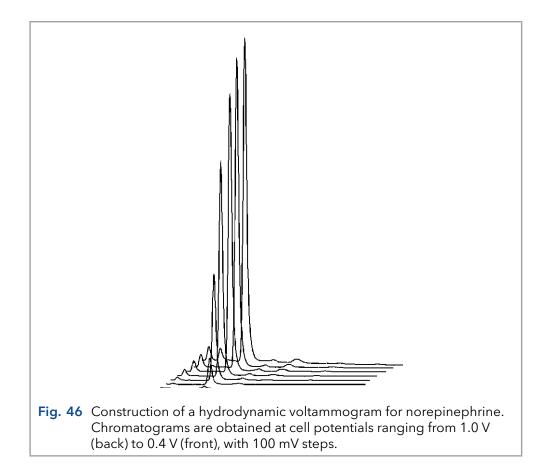

9.6 Construction of a scanning voltammogram

A scanning voltammogram can be recorded using the ECD 2.1 scan mode. The scan mode is programmed in the 'SCAN SETUP' screen of the ECD 2.1. Depending on the data acquisition software that is used and the experimental set-up, a full, half or continuous scan cycle can be chosen.

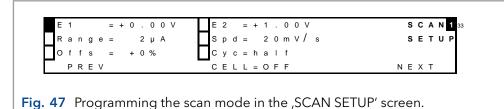


With the Elite Dialogue control software a scanning voltammogram can be programmed under the 'Detector' tab in the main window.

Nonitor Detector	Events	S pump	Log						
Cell	1	•	🔽 On	🔽 Enabled	Mode	Scan	¥		Dev status
Output					Potential				
Range	2 [µA]				E1	0.00	[1]		
Offset	0		[%]		E2	1.00	[V]		
Filter	0.001		[Hz]		scan rate	20	•	[mV/s]	
Data rate	1	· •	[Hz]		sweep	half	•		
Polarity	(i) +	0 -							
Compensation	🔲 On	Autoz	ero						
Temperature					Run time				
Oven	🗸 On				t	0.83	3	[min]	
Set	35 [°C]							
Measured	35.5 [°C]							

In the above example a 'half' scan is programmed, sweeping the potential from 0 V to 1.0 V. A full scan would include the reverse scan, i.e. from 0 V to 1.0 V and back to 0 V. In the continuous mode the voltage is swept up and down between both potentials for a pre-defined run time.


A convenient way to record a scanning voltammogram is by direct infusion of analyte in the flow cell using a syringe pump. In Fig. 49 a scanning voltammetry set-up is shown.

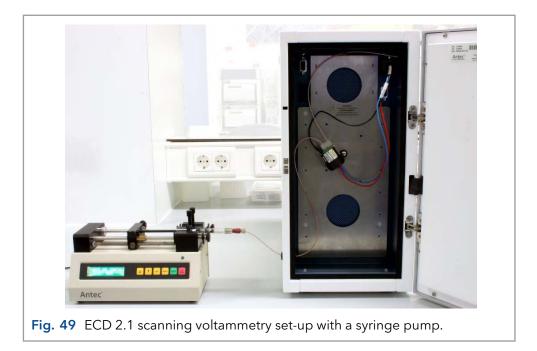

The Antec dual syringe infusion pump which can be obtained as accessory has as an advantage that it can be controlled in Elite Dialogue software as well.

Monitor Detector Events Spump Log	
Monitor Pump started to infuse	Start Stop
Control	Status
Direction 💿 Infuse 🔘 Withdraw	Volume 638.347 [µL] Reset
Flow rate 2.500 [µL/min]	
Target vol. 0.000 [µL] 🔲 on	
Units	Syringe
	Diameter 4.61 [mm]
© mL/min ⊚ mL/hr	Selector 🗸

In the example in Fig. 51 a half scan is shown at a flow rate of 10 μ L/min of a 20 μ M Serotonin (5-HT) standard in mobile phase.

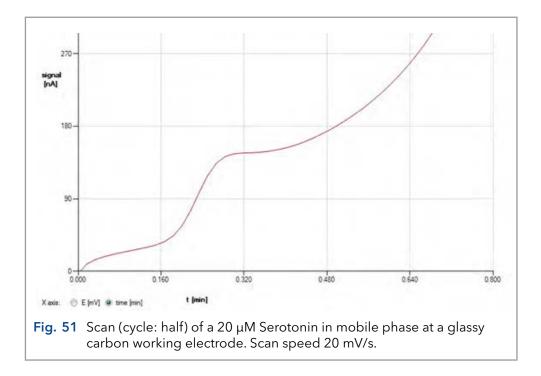

The following procedure is advised to record scanning voltammogram of analytes:

 Use a voltammetry set-up as shown in Fig. 57 (preferably) in combination with the Elite Dialogue software and a syringe pump for direct infusion of analyte in the flow cell.




If Dialogue is not used, connect an A/D converter to the analog output of the detector to record the cell current. Set the A/D converter sampling rate to 1 Hz. This is the same frequency as the voltage steps during the scan. If a higher sampling frequency is chosen a typical stepwise pattern may appear. Note that with such set-up only 'I versus t' curves can be obtained.

- Prepare a solution of the pure compound dissolved in (preferably) the HPLC buffer with a concentration in the range of 10-100 μM.
- Set the lower (E1) and upper potential (E2), the scan rate (Spd), range and scan cycle in the 'SCAN SETUP' menu or under the 'detector' tab in Dialogue.

Typical scan settings to start with: E1 0 mV, E2 1000 mV, scan speed 10 mV/s, range 5 μ A/V. Optimize the settings if required for your specific compounds.

- Prime a 1 mL glass syringe with analyte solution and install it in the syringe holder of the pump.
- Program the syringe diameter and flow rate in the syring pump settings menu. In case a syringe pump is used which can be controlled in Dialogue the settings can be programmed under the 'S pump' tab. A typical flow rate to start with is 10 µL/min. Optimize the flow rate if required during the scan experiments.
- Start the syringe pump and before scanning assure that the flow cell is sufficiently primed with analyte solution.

- A scan can be started (stand-alone) by pressing the 'START' button in the 'SCAN STAT' menu or by starting a single run in Dialogue: press the 'F5' button or click 'start single run' under the 'Options' menu.
- To record a background (blank) scan repeat the experiment with the pure HPLC buffer in which the analyte was dissolved.

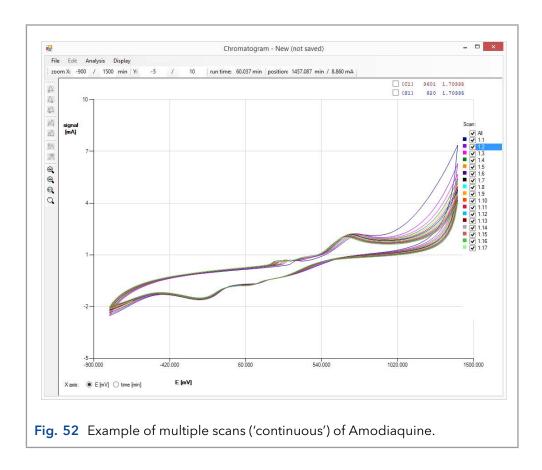

10. Technical data

10.1 Environmental, dimensions, weight and power requirements

Working temperature	10 - 40 °C (indoor use only)
Storage temperature	-25 - +50°C
Humidity	20 - 80% RH
Safety and EMC	According to EC-directives; Emission Group I Class A; cMETus approved
Equipment class	1
Installation category	II
Pollution degree	2
Dimensions	43 (D) x 22 (W) x 44 (H) cm = 16.9" (D) x 8.7" (W) x 17.3" (H)
Weight	max 14.4 kg (32 lbs) without flow cell and column
Installation	Install upright on flat & smooth surface, keep space under the detector free (risk of blo- cking power supply fan).
Power requirements	100-240 VAC, 50/60 Hz, 260 VA, auto-sensing
Mains fuse	2.5 AT / 250V, 5x20 mm, IEC 60127-2
	For safety reasons, any other internal fuse or circuit breaker is not operator accessible, and should be replaced only by authorized personnel by the manufacturer. Only use manufacturer-supplied fuses.



For optimum analytical performance it is recommended that the ambient temperature of the laboratory be between 20-25 °C and be held constant to within ± 2 °C throughout the entire working day. Note: that for optimal performance of the oven the oven temperature should be set at least 7 degrees higher as ambient temperature.

10.2 General

Operating modes	DC, PULSE SCAN
Other mode	CONFIG, DIAG and SERVICE
Sensors	Up to 3 flow cells
Autozero	triggered by keyboard, rear panel TTL, or remote PC control (LAN)
Max. current compen- sation (Autozero)	25 nA - 2.5 mA in DC and PULSE mode de- pendent on range setting
Offset	+50% to - 50% of max. output voltage, 5% steps
PC control	Parametric control and data-acquisition via LAN port (USB service port)
Embedded software	Flash technology, upgradeable via PC (USB)
Oven	+7 °C above ambient to 60 °C, accuracy 0.5 °C, stability 0.1 °C; accommodates column and flow cell(s)
Rear panel connectors	1x IEC inlet (Mains), 1x USB B, 1x RJ45 LAN, 1x 9-pins sub-D Male (Valve), 1x 9-pins sub-D Female (Analog output), 1x 25-pins sub-D Female (Digital I/O)
Analog output (DAC)	-1 to +1 V full scale (via 16-bit D/A converter)
Analog output (I/E)	-2.5 to +2.5 V full scale (unprocessed I/E con- verter signal)
Digital I/O (HW)	2x Relay, 5x TTL outputs (CMOS 3.3V logic), 13 TTL inputs (programmable), 1x GND
Programmable I/O functionality	Cell on, Cell off, Autozero, Start, Overload, Relay, Auxiliary
Valve control	VICI valco 2-pos electrically-actuated valve (E2CA, EHCA) via serial cable, manual valve, 1x inject marker output

10.3 DC mode

Range	10 pA - 200 μA in 1, 2, 5 increments
Filter (ADF)	10 - 0.001 Hz in 1, 2, 5 increments
	RAW and OFF: for unprocessed data
Potential (Ec)	-2.50 V to + 2.50V with 10 mV increments
Data Rate	1 - 100 Hz in 1, 2, 5 increments, dependent on filter setting
Noise	< 2 pA with dummy cell (load of 300 MΩ/ 470 pF) in 1 nA range, filter off, Ec +800mV and temperature of 35 °C.

10.4 PULSE mode


•••••••••••••••••••••••••••••••••••••••	
Range	10 nA - 200 µA in 1, 2, 5 increments
Filter (ADF)	0.5 - 0.001 Hz in 1, 2, 5 increments
	OFF: for unprocessed data
Potential (Ec)	-2.50 V to + 2.50 V with 10 mV increments
Data Rate	1/(pulse duration) Hz
Waveform	Max 5 potential steps
Pulse times (t1-t5)	t1: 100 ms - 2000 ms; t2, t3, t4, t5: 0 - 2000 ms in 10 ms increments
Sampling times (ts)	20 ms - [t1 - 60] ms

10.5 SCAN mode

Range	10 nA - 200 μA in 1, 2, 5 increments
Potential (Ec)	-2.50 V to + 2.50V with 10 mV increments
Data Rate	1 Hz
Scan rate	1 - 100 mV/s in 1, 2, 5 increments
Cycle	Half, Full, Continuous

11. Rear panel I/O

In this chapter all rear panel functionality is described. The ECD 2.1 has besides the mains inlet in total 5 connectors on the rear panel for communication, data output and I/O. A photo of the rear panel connectors is shown below in Fig. 53 for reference.

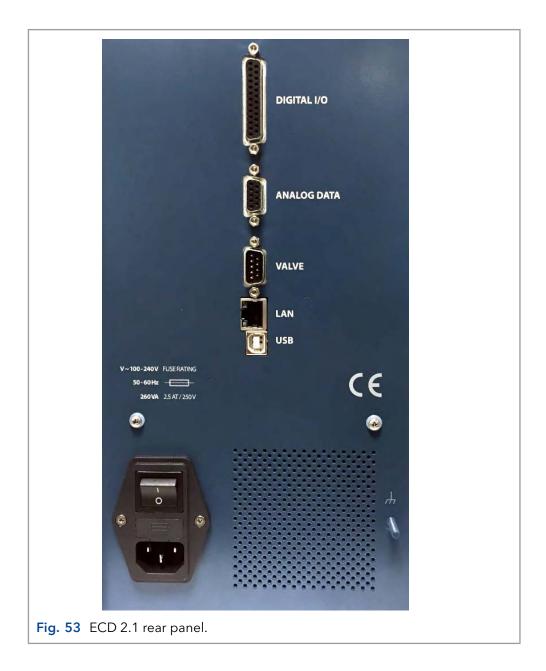

11.1 USB B connector

USB type B connector for serial instrument control over USB, for service use only:

- Based on USB-to-serial UART interface using the FT232R chip from FTDI (Future Technology Devices International Ltd).
- FT232R is fully compliant with USB 2.0 specifications.
- Fixed communication baud rate: 921600 bps.
- Communication over USB is used for software (FW) update of the instrument only.

11.2 LAN connector

RJ-45 bus for serial instrument control over LAN:

- 10Base-T or 100Base-TX (Auto-Sensing) serial-to-ethernet connectivity.
- Network configuration of Xport via Lantronix device installer software utility.
- Fixed communication baud rate: 921600 bps
- Communication over LAN is used for parametric instrument control and data-acquisition.



To establish communication over LAN, the LAN cable must be connected when starting the detector using the mains power switch on the rear panel. If no communication cable (either LAN or USB) is connected during start-up, communication via USB is enabled (default).

Please consult the installation section for details about configuration and set-up of communication over LAN.

11.3 VALVE connector

Serial D9 (9-pins subD male) valve control connector for electrical (VICI) and manual valves:

- Serial (RS232) valve control of VICI 2-position electrically actuated valves via pin 1 - 3 (see VICI technical note <u>tn413.pdf</u> on the official <u>www.vici.com</u> web site). Compatible with the E2CA and EHCA actuator control module.
- Automatic detection of electrically actuated VICI valve, valve control in: STAT screen and Command ID 30 (0=load, 1=inject).
- Manual valve status (load/inj) read-out via pin 5 and 6 (contact closure).
- Inject Marker (pin 9) TTL output. Default: high 3.3V (load), on inject event: low (0 V) for 2000 ms.

Valve connector layout

In the table below the connector layout is shown for the Valve connector.

Table V. Valve connector lay-out.

Pin	Layout
1	GND
2	TRANSMIT(TxD)
3	RECEIVE (RxD)
4	
5	GND
6	SWITCH (Hand valve)
7	-
8	-
9	INJECT MARKER


For a manual valve the status (LOAD/INJ) read-out is established by connecting the LOAD/INJ sensor wires/leads of a manual valve to pin 5 and 6 (contact closure).

Inject marker

A connected valve enables the inject marker (pin 9)|. The contact is high when the valve is in 'load' position (3.3V), and low (0V) in the 'inject' position. On an inject event the contact will go to 0V (low) for 2000 ms, It can for example be used to start the integration software when injection is done.

Electrical valve configuration

A VICI electrically-actuated 2-position valve is automatically detected during start-up of the instrument. Automatically the Valve parameter in the ECD 2.1 IO menu is set to Valve = Present.

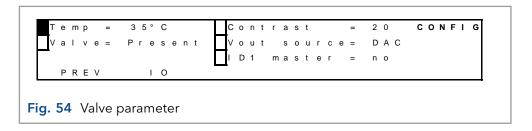



Always restart the detector (by re-powering via the mains switch on the rear panel) when connecting the valve for correct initialization of an electrically-actuated VICI valve.

In the case an electrical VICI valve is detected, the Dialogue Elite software automatically shows the presence of the electrical valve in the device settings window (EC device) and valve control buttons will appear in the detector window.

Settings Acquisition EC device IP switch			
Device info	Serial number	s	
Device: DECADE ELITE	Device:	17530009	
Firmware: 1.00	Control:	15052	
Boot version: 0.08 Checksum: 557841297	Sensor 1:	14012	
User: hendrik-jan.brouwer @ PC-W	Sensor 2:	14010	
IP addr.: 192,168,5,1			
Phys. addr.: 00-80-A3-AC-3B-CF			
comm time-out: 400 [ms]	External valve		
	Manua	I valve 🖂 Electrical	
		i vaive 🕑 Electricat	Jaive
Read from device		Send to device	Close

🚪 Dialogue Elite - I	DECADE ELITE		- 🗆 X
	ptions Events Scripts Window Hel	0	775, 1478
Monitor Detector	Events Scripts Log		
Cell	1 V On V Enabled	Mode DC ~	Dev status
Output		Potential	
Range	50 [nA] 🗸	E -0.01 [V]	
Offset	0 ~ [%]		
Filter Data rate	off ~ [Hz]		
Polarity	• + • •		
Compensation			
Temperature		Run time	
Oven	On	t 15.000 [min]	
Set	off ['C]	In/Output	
Measured	28.6 [°C]	Relay TTL contact 1 2 1 2	3 4
		El. valve position:) Inject sa immediately
Read from device]		Send to device
DECADE ELITE: C	OM3 🔲 send/receive		
56 ECD 2.1 de	tector window wit	th Valve control	

Manual valve configuration

In case of a manual valve the valve present parameter has to be set manually. Configuration of a manual valve can be done via the LCD screen or the Dialogue Elite software.


LCD display: Configuration of a manual valve can be done by setting the Valve parameter in the IO menu manually to Valve = 'Present'.

Dialogue Elite: Go to the device settings and the EC device tab and check the manual valve check box.

	Settings Acquisition EC device IP switt Device: DECADE ELITE Firmware: 1.00 Boot version: 0.08 Checksum: 557841297 User: hendirk-jan.brouwer @ Pi IP addr.: 192.168.5.1 Phys. addr.: 00-80A3AC-3B-CF comm time-out: 400 [ms]	Serial numbers Device: 17530009 Control: 15052 Sensor 1: 14012
Fig. 57 Dia in re		e settings, manual valve check box (encircled

Valve status & control


In stand-alone mode the valve status for both an electrically-actuated valve and manual valve is displayed in the STAT screens on the LCD display. For example see the screen dumps of the STAT screen for a ECD 2.1 SCC (top) and TCC detector. INJ=L represents the LOAD position (position A on the actuator control module of an electrically-actuated valve) and INJ=I represents the INJECT position (B on the actuator control module). The LOAD position is the default position at start-up of the instrument in the case of an electrically-actuated valve.

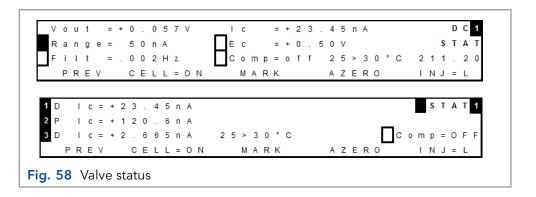

In stand-alone mode the valve position can be controlled via the F5 function button in the STAT screen. In the case of a manual valve the information above the F5 button (INJ=L) in the STAT screen shows the status (position) of the valve only, no valve control is possible with this type of valve.



Make sure that the valve flow path is connected correctly and LOAD corresponds with position A and INJECT with position B. See example in Fig. 59 for a 6-port valve.

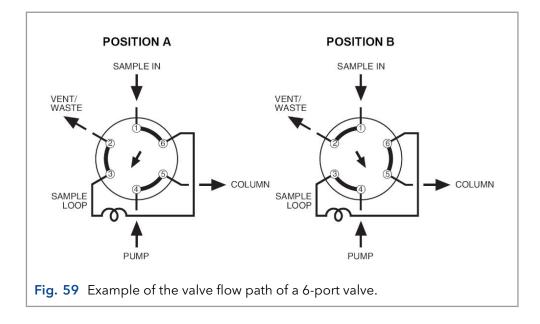

In Dialogue Elite the valve status/position is shown under the monitor tab. To make it visible in the left table, check the valve position checkbox in the item selector (by default the valve position is not shown).

Monitor	Mode	1 Pulse	 ✓ Auto update ✓ Item selector ✓ Mode 	_
	Cell current An. output [V] Potential E [V] Auxiliary E [V] Temperature [*C]	Off 0.000 -0.01 0.00 29.6	Cell current An. Output [V] Potential [V] Auxiliary E [V] Range Filter [Hz]	à
	Valve pos. Autozero Cell on/off Cell enabled	L set	☐ Offset [%] ☐ Offset [%] ☐ Syr flow rate [µL/min] ☐ Syr volume [µL] ☐ Pump 1 flow rate [mL/min] ☐ Pump 1 pressure [bar] ☐ Syr 1%A]	
			AS tray temp. [*C]	

An electrically-actuated VICI valve can also be controlled via Dialogue Elite under the detector tab, see Fig. 56. Note that the valve will switch immediately after a change from Load to Inject (checkboxes) or vice versa, no 'Send to device' action is required.

An electrically-actuated VICI valve can also be controlled via timed events under the events tab in Dialogue Elite. In the Event editor the event 'valve position' is available to switch the valve in the Load/inject position. See example in Fig. 61.


	and the second se	parameter	setting	cell	comment	
	0.00	Start mode	with 'start events'	1		
_	0.01	Valve Position	Inject	1	Valve switching to inject	
	0.05	Valve Position	Load	1	Switching valve back to	o Load position
		Runs	5 1 done: 0	ext start	trigger 🗹 stop run	Run time check
	Edit	Runs		ext start	trigger ⊠ stop run Stop	Run time check
 Eve	12220	Time			Stop	
Ev	ents: C:\Us Read from	Time	e 0.05 [min]		Stop	Start Send to
Eve	ents: C:\Ue	Time	e 0.05 [min]		Stop	Start



For more detailed information about valve control of the electrically-actuated VICI valves in the ClarityChrom® chromatography software or Dialogue Elite please refer to the corresponding instructions.

11.4 ANALOG DATA connector

The ECD 2.1 is equipped with an analog data output connector to provide the measured signals in millivolts (mV) for users who work stand-alone without PC control with the instrument. The ANALOG DATA connector (D9-female) can either be connected to an X-Y recorder or A/D converter. The ANALOG DATA output can supply either a non-manipulated signal directly from the I/E converter, or the processed data signal by the CPU of the ECD 2.1. The type of output can be selected from the CONFIG screen by setting the parameter Vout source to either DAC or I/E.

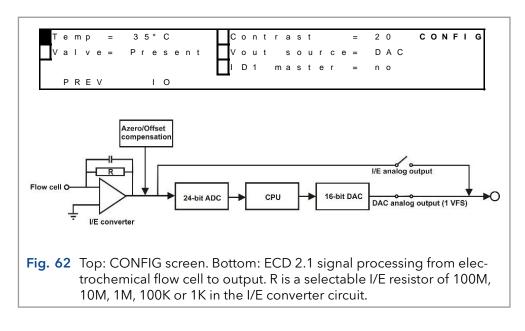

The output parameter Vout source can also be selected/set using Dialogue Elite. Go to the device settings menu and Acquisition tab and set the analogue output to DAC or I/E converter. The instrument is by default set to DAC. For regular measurements it is advised to use the DAC output.

Output Full scale 1 V Polling 2 V Units PA V auto	Plotting in graph: at max. 1 Hz at max. 5 Hz show t=c./f ext. start trigger CRC check	Other channels Actual temperature Cell current Eaux	
Noise monitor	pA after 0 [min] restarts	Analog Output	>
Read from device		Send to Close	

DAC output

The DAC output is the processed signal by the ECD 2.1's CPU and is <u>iden-</u> <u>tical</u> to that of the digital cell current signal obtained via data-acquisition over LAN using the Elite Dialogue or the ClarityChrom[®] chromatography software.

The signal in mV from this output is directly related to the range setting of the ECD 2.1. So for example in the case the 200 μ A measurement range is selected 'Range = 200 μ A' the DAC signal on the analog data connector corresponds with +1000 mV = +200 μ A and -1000 mV = -200 μ A (so +/- 1V full scale). To convert the signal in mV to the actual cell current in nA use the following calculation:

Ic (nA) = Vout (mV) x Range setting (nA) 1000 mV

So for example a signal on the output of 250 mV in the 200 nA range corresponds with an actual cell current of (250/1000)*200 nA = 50 nA.

I/E output

The I/E output is the unprocessed analog signal from the I/E converter circuit. This signal is a true analog signal which is as close as practically possible to the working electrode (WE). The maximum output voltage of the I/E output is +/- 2.5V under all conditions and is independent of the detector range setting. The signal in mV is related to the selectable I/E resistor of 100M, 10M, 1M, 100K or 1K in the I/E converter circuit. The actual cell current can be calculated from the analog I/E output in Volt using the following formula:

 $Ic (Ampere) = \frac{Vout (Volt)}{RI/E (Ohm) \times 10}$

So for example an I/E output signal of 250 mV in the 5 nA range (RI/E = 10 M Ω) corresponds to an actual cell current of 0.25V/(10 x 1E7 Ω) = 2.5E-9A = 2.5 nA. In the table below, the I/E resistor value is listed for every range setting.

Table VI. ECD 2.1 I/E converter resistors:

Current Ranges	I/E Resistor (Ohm)
10рА, 20рА, 50рА, 100рА, 200рА, 500рА and 1nA	100M
2nA, 5nA	10M
10nA, 20nA, 50nA	1M
100nA, 200nA, 500nA, 1μA, 2μA	100K
5μΑ, 10μΑ, 20μΑ, 50μΑ, 100μΑ, 200μΑ	1k

In the detector accessory kit a dedicated Output cable, D9 male - open, 2m (pn 250.0128A) is supplied. It is advised to use this manufacturer-supplied cable for this type of measurements.

For reference, the layout of the analog data out connector is shown in the table below.

Pin	Layout
1	Vout 1
2	Vout 2
3	Vout 3
4	Vout 4
5	Vout 5
6-9	GND

Table VII. Analog data output connector lay-out.

To measure for example the analog signal of cell 1 with an external A/D converter:

- Connect the signal wire, lead of pin 1 (Vout1), to the analog measurement channel of the A/D converter.
- Connect the GND wire, lead of pin 6 (or 7-9V), to the corresponding analog ground connection of the A/D converter.

11.5 Digital I/O connector


The detector has one 25-pins digital I/O connector which enables control of (or by) external equipment. The IO connector contains 18 TTL contacts (5 outputs and 13 inputs, 3.3V CMOS logic), 2 RELAYS (contact closure) and 1 ground (GND connection).

TTL inputs & outputs

Both the TTL inputs and outputs are default = high (3.3 Volt). The TTL inputs are level triggered: the contacts require a minimum TTL-low pulse duration of 100 ms. If multiple activations are required the next pulse should be given after 100 ms TTL high. When the input is kept low, only one activation will occur.

Relays

The ECD 2.1 has 2 free programmable contact closure outputs:

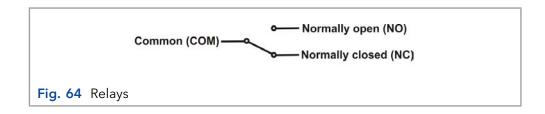

- Relay1: pin 1 normally closed, pin 2 normally open, pin 3 common.
- Relay2: pin 4 normally closed, pin 5 normally open, pin 6 common.


The maximum rating for these contact closure outputs are 24 VDC (switching voltage) and 0.25 A. The relays can be controlled in the Elite Dialogue software and ClarityChrom[®] Chromatography software.

AUX

The ECD 2.1 has 4 free programmable TTL outputs AUX1 - AUX4 (pin 7 - 10). These contacts are default 'high' 3.3V (inactive), when active the status is 'low' 0V.

Overload

The overload output (pin 11) can be used to monitor if the cell current goes out of range during a chromatographic run. An 'Out of range' error appears when the cell current Icell exceeds the limit of the current range at which the measurement is performed (see Fig. 65).

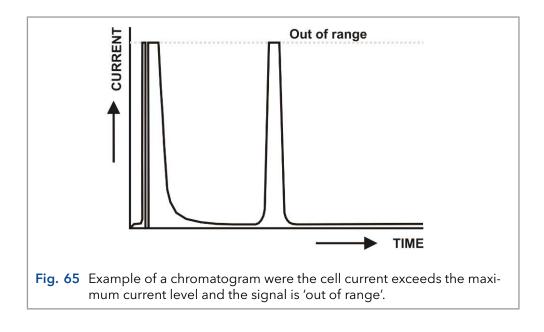




It is important to recognize an 'out of range' (overload) situation, because it may lead to erratic results when quantifying analyte concentrations in samples.

If for instance the cell current goes in out of range during the recording of an analyte peak, it can be (in most cases) easily recognized by a flat top of the peak and a very abrupt transition to the flat top at the edges.

By default the status of the overload output is 'high' 3.3V. When the cell current has the status 'out of range' the overload output will change status to 'low' 0 Volt, until the cell current returns to a value within the measurement range again. The overload output (pin 11) is one of the configurable I/O's.

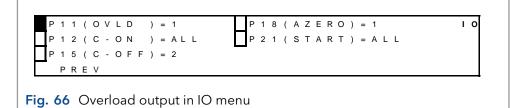

The configurable I/O's can be programmed in the IO menu, which is a submenu of the CONFIG menu. By default, the overload output is assigned to cell 1: 'P11 (OVLD) = 1'. This means that only when the cell current of cell 1 is 'out of range' the status of the overload output will change to 'low' OV. For all other cells present in the ECD 2.1 (in case of a DCC or TCC version of the instrument) an out of range situation will not trigger a response on pin 11.

The following options can be selected for the configuration of pin 11:

P11(OVLD) = 1	Overload output active for cell 1 only
P11(OVLD) = 2	Overload output active for cell 2 only
P11(OVLD) = 3	Overload output active for cell 3 only
P11(OVLD) = 4	Overload output active for cell 4 only
P11(OVLD) = 5	Overload output active for cell 5 only
P11(OVLD) =	Overload output inactive
P11(OVLD) = All*	Overload output active for all cells present

*) When this option is selected, the overload output will be active for all cells present in the ECD 2.1. If the cell current of either one of those cells will go 'Out of range' the overload input pin 11 will become active.

Cell on, Cell off

The ECD 2.1 has 3 TTL inputs to switch on cells (pin 12 - 14) and also 3 inputs to switch off cells (pin 15 -17) AUX1 - AUX4 (pin 7 - 10). This input command can be used for example to switch on and stabilize the flow cell early in the morning by means of a timer. Two of the inputs are configurable (pin 12, pin 15, cell on and off respectively) in the IO menu (see previous chapter about the overload output). The configuration settings of these inputs are: 1, 2, 3, 4, 5,' ', and all. In case 'all' is selected all cells present in the ECD 2.1 will be switched on or off when the corresponding input is triggered.

Autozero

The ECD 2.1 has 3 TTL inputs (pin 18 – 20) available to autozero the cell current of the cell(s). Triggering these inputs enables external activation of the auto zero command. This function is active only when the 'Icell' is displayed. One autozero input is configurable (pin 18) in the IO menu (see previous chapter about the overload output). The configuration settings of this input are: 1, 2, 3, 4, 5,' ', and all. In case 'all' is selected the cell current of all cells present in the ECD 2.1 will be zeroed when the input is triggered.

Start

The ECD 2.1 has 4 TTL inputs (pin 21 – 24) available to start data-acquisition and/or start a scan. One start input is configurable (pin 21) in the IO menu and can be used for example to start the data-acquisition of all cells synchronously using only one trigger input when the setting 'All' is selected.



The manufacturer will not accept any liability for damage, direct or indirect, caused by connecting this instrument to devices that do not meet the relevant safety standards.

Programming outputs

In the Dialogue Elite software the ECD 2.1 outputs (Relays and AUX) can be controlled and programmed. This can be done via the Detector tab on the main window and via timed event tables. By default the I/O is hidden and not shown in the monitor tab. To be able to see the status and control buttons for the I/O go to the Device settings window (settings tab) and check the 'Show status outputs' checkbox under the controls section.

Cell	1 v 🗋 On 🗹 Enabled	Mode DC V Dev status	
Output		Potential	
Range	50 [nA] 🗸	E -0.01 [V]	
Offset	0 ~ [%]	12	
Filter	off 🛛 🗸 [Hz]		
Data rate	10 ~ [Hz]		
Polarity	• • • •		
Compensatio	on On Autozero		
Temperature		Run time	
Oven	🗌 On	t 15.000 [min]	
Set	off [°C]	In/Output	
Measured	28.4 [°C]	Relay TTL contact	
		a. valve position: Load Inject 	
		valve switches immediately	
Read from		Send to	

The status of the Relays and TTL inputs can be set to active (\square) or inactive (\square) . The selected state of the outputs will be activated after clicking the 'Send to device button'.

The outputs can also be switched via timed events under the events tab. By opening a 'New event' under the Events pull-down menu, a new Event table/file can be created. Subsequently, the Event editor will open. Select the event 'Output A' and assign a time at which the output should be switched and select which output should be activated (Relay, Aux) under the value field. See example in Fig. 68. Make sure to program a second step to inactivate the Output again and put it back in its initial inactive state again.

All events	Selected events					
[Common] A Start mode Repeat all events Repeat all events End events End events End events End events End events events Acquistion nold Noise monitor Run time Scaling XY [File actions] Save chromatogram Export to Excel Select Excel templat Excel Worksheet/Cc Load method [EC settings] Cel on/off Range Auto set unit Filter Autozero Offset	Start mode Durput A	Settings time Cell/Dev. ID Comment Value	relative: to p	[min]		
Polarity Y	Up Down Delete	[Auto save	Sort	Close	

	t [min]	parameter	setting	cell	comment
	0.00	Start mode	with 'start events'	1	
	0.01	Output A	relay 1	1	Switching Relay to active
	0.02	Output A	inactive	1	Always add an inactivation step as well!
	Edit	Run Tim		ext start tr	igger ⊻ stop run v Run time check Stop Start
Ev	vents: C:\U	Tim			Stop Start
Ev		Tim	e - 0.02 [min]		

0	1
O	
-	

Table VIII	DIGITAL I/C	connector	lavout
	DIGITALI/C	connector	layout.

Pin	Type I/O	Function (default)	Configu- rable I/O	Description
1,2,3	Relay 1	Relay 1	-	Contact between 3 (common) and 1 (default), activated by time file Outp 0100 or Command 47
4,5,6	Relay 2	Relay 2	-	Contact between 6 (common) and 4 (default), activated by time file Outp 0001 or Command 47
7	TTL OUT	AUX 1	-	Free programmable TTL output, activa- ted by time file Outp 0001, Command 47
8	TTL OUT	AUX 2	-	Free programmable TTL output, activa- ted by time file Outp 0010, Command 47
9	TTL OUT	AUX 3	-	Free programmable TTL output, activated by Command 47
10	TTL OUT	AUX 4	-	Free programmable TTL output, activated by Command 47
11	TTL OUT	Overload	A*, 1-5	Active in case of signal overload ('OUT OF RANGE', 'PAD OVLD')
12	TTL IN	Cell on 1	A, 1-5	Trigger to switch the cell on
13	TTL IN	Cell on 2	-	Trigger to switch the cell on
14	TTL IN	Cell on 3	-	Trigger to switch the cell on
15	TTL IN	Cell off 1	A, 1-5	Trigger to switch the cell off
16	TTL IN	Cell off 2	-	Trigger to switch the cell off
17	TTL IN	Cell off 3	-	Trigger to switch the cell off
18	TTL IN	Autozero 1	A,1-5	Trigger to zero/null the cell current (com- pensation lc)
19	TTL IN	Autozero 2	-	Trigger to zero/null the cell current (com- pensation lc)

Pin	Type I/O	Function (default)	Configu- rable I/O	Description
20	TTL IN	Autozero 3	-	Trigger to zero/null the cell current (com- pensation lc)
21	TTL IN	Start 1	A,1-5	Trigger to start a Scan and Data-acquisition
22	TTL IN	Start 2	-	Trigger to start a Scan and Data-acquisition
23	TTL IN	Start 3	-	Trigger to start a Scan and Data-acquisition
24	TTL IN	Start 4	-	Trigger to start a Scan and Data-acquisition
25	GND	GND	-	Ground connection

*) A = All boards.

In the detector accessory kit, a dedicated I/O cable is supplied with an I/O connector board with screw terminal connections for easy connection of open-ended electrical wiring for I/O:

ECD 2.1 I/O conn. Board ECD 2.1 I/O cable 25M-25M, 1.8m


For use, connect the I/O cable to the 25-pins I/O connector on the detector rear panel, subsequently connect the I/O connector board to the end of the I/O cable.



On the rear side of the I/O connector board a label is adhered for reference with the lay-out of the digital I/O connector as described in Table VIII.

11.6 Chassis grounding stud

On the rear panel a chassis grounding stud is available at the lower right-side, next to the ventilation holes of the power supply compartment. This grounding stud which is connected to the central grounding point of the instrument can be used for shielding purposes. For example, to shield the flow cell outlet from external electrical interferences (troubleshooting noise issues), or to shield the flow cell from other equipment which might be coupled in series with the electrochemical detector such as an MS.

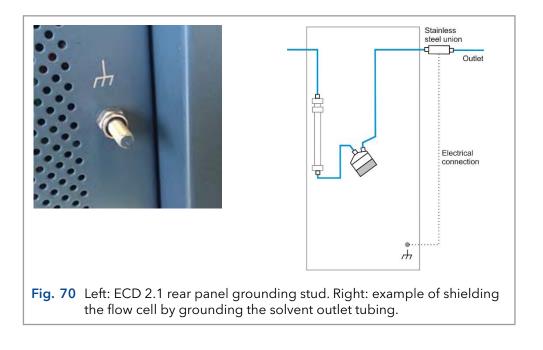



Use the chassis grounding stud for shielding only, not for safety grounding.

12. Functionality tests



Note: Standard processes regarding IQ ans OQ in single devices may be handled differently in individual cases.

12.1 Installation Qualification (IQ)

The customer may request the Installation Qualification, which is free of charge. In case of a request, the Customer Support of KNAUER or from a provider authorized by KNAUER performs this functionality test during the installation.

The Installation Qualification is a standardized document that includes the following:

- confirmation of flawless condition at delivery
- check if the delivery is complete
- certification on the functionality of the device

12.2 Operation Qualification (OQ)

The Operation Qualification includes an extensive functionality test according to KNAUER standard OQ documents. The Operation Qualification is a standardized document and free of charge. It is not part of the delivery. Please contact the Customer Support in case of request.

The Operation Qualification includes the following:

- definition of customer requirements and acceptance terms
- documentation on device specifications
- device functionality check at installation site
- **Test Intervals** To make sure that the device operates within the specified range, you should test the device regularly. The test intervals depend on the usage of the device.
 - **Execution** The test can be carried out either by the Customer Support of KNAUER or from a provider authorized by KNAUER (for a fee).

13. Troubleshooting

Even though great care was taken in the design of the ECD 2.1, problems may occur during operation of the instrument. The information in this chapter may help you to identify and solve the source of the problems. Errors can be categorized in two types:

- Instrument errors
- Analytical problems

In the next sections both types of errors are described. In the event that the problems cannot be solved after following the instructions in this troubleshooting section, contact your local supplier for further assistance.

13.1 Instrument errors

Incidental fault conditions may occur in any instrument. The ECD 2.1 will generate an error message containing an error number with a short description for several hardware fault conditions, which will be shown in the instruments LCD display.

Table IX. Error messages.

11 Checksum error	
12 Temperature sensor 1 error	
13 Disconnect flow cell x	
14 Control board SRAM error	
20 Checksum error	

Please contact your local supplier if one of the above errors occur for further instructions. In the case the instrument does not power up at all check the following remedies.

No detector response

Possible cause	Remedy
No power	Check line voltage setting, plug in power cord
Power switch off	Turn this switch ON (at the rear panel)
Faulty fuse	Replace fuse
Divergent mains voltage	Check line voltage




Make sure the ECD 2.1 is connected to a grounded power source with a line voltage which is within the specified ratings. If the ECD 2.1 does not respond, a fuse in the mains inlet may be blown. Furthermore the following messages can be displayed on the LCD screen or PC control software during a measurement:


Table	эX.	Messages.

Message	Advice
01 Out of range*	Output is either above +1.0V or below -1.0V. Pressing AZERO may give an adequate read-out again. If the message remains after pressing AZERO, the autozero function is unable to compensate the background cell current. Advice: use a less sensitive range in the SETUP menu.
02 PAD overload	Charging current in pulse mode out of range. Pressing AZERO may give an adequate re- ad-out again. If not, it is advisable to change the pulse settings (increase t1) or use a less sensitive range.

Make sure that maintenance is performed on a regular basis.

*) An 'Out of range' error appears when the cell current Icell exceeds the limit of the current range at which the measurement is performed (see Fig. 71).

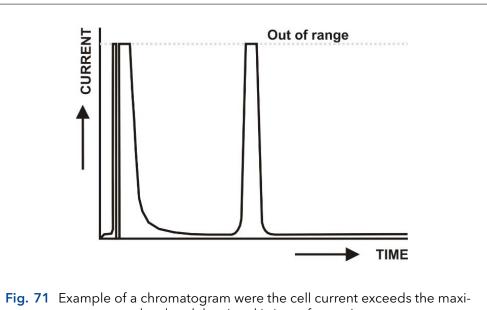

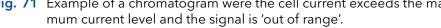



It is important to recognize an 'out of range' (overload) situation, because it may lead to erratic results when quantifying analyte concentrations in samples.

13.2 Analytical troubleshooting

Analytical problems like loss of signal, increase in noise level, high cell current, loss in sensitivity etc. may occur in any (U)HPLC-ECD 2.1 system. It may be hard to find the cause and several checks need to be performed to eliminate the source of the problem. The first step is to determine whether the problem is caused by the ECD 2.1 electrochemical detector or the rest of the (U)HPLC system. For that purpose two basic checks should be performed by the end-user:

- Dummy cell test
- Stop flow test


Dummy cell test: The outcome of the dummy cell test, which is described in the next paragraph will give an indication if the problems are caused by the detector hardware (electronics).

Stop flow test: The stop flow test will eliminate if the problems are caused by the electrochemical flow cell, or is originating from the rest of the (U)HPLC system (pump, autosampler, pulse damper, column, mobile phase etc.).

13.3 Dummy cell test

External dummy cell

An external dummy flow cell is shipped with every ECD 2.1 instrument for troubleshooting purposes and maintenance checks. The dummy cell test can be performed standalone via the LCD display in combination with an A/D converter or via the Dialogue Elite / ClarityChrom® software. A successful dummy cell test confirms that the controller, including the cell cable, functions properly. If the result of the noise measurement with the dummy cell is within specs, the controller is excluded in a troubleshooting procedure.

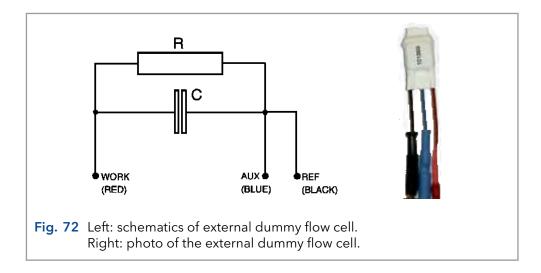

The dummy flow cell consists of a resistor (R) of 300 MOhm and a capacitor (C) of 0.47 μ F in parallel. The current is measured over the resistor according to Ohm's law (V = I x R), hence with a working potential of 800 mV the current drawn will be about 2.67 nA. Slight differences as to this (ideal) value are due to the tolerance of the resistor (± 1%). The capacitor functions as a 'noise generator' and in fact resembles the capacitance of a well-functioning VT-03 flow cell in an ideal (U)HPLC set-up. The noise generated via the dummy should be less than 2 pA if the filter of the controller is set to off, provided that the dummy is within the fully closed Faraday shield at the same position as the flow cell.

Parameter	Setting	
Cell potential	800 mV	
Oven	35 °C, stable	
Filter	Off	
Range	1 nA/V	

Test criteria:

- I cell = 2.67 +/- 0.05 nA
- Noise < 2 pA</p>



The results (cell current and noise) of the dummy cell test should be within the above mentioned test criteria. If the current values for Icell and noise are not within the criteria it is an indication that something could be wrong with the detector hardware. Please consult your local representative.

Internal dummy cell

The ECD 2.1 also has the option to run a so-called internal dummy cell test. This exclusively checks the performance of the electronic circuit boards (amplifier circuitry) only, so it excludes the cell cables and the external dummy flow cell. From the MAIN screen DIAG can be selected to enter the DIAG screen, followed by selecting NOISE. This activates a timer in the NOISE screen, and after 5 min stabilization auto zero is activated and the dummy cell test is ready. Noise of the internal dummy cell can be measured at the output. As with the external dummy cell the noise should be better than 2 pA. Detector settings in the NOISE screen are the same as in the external dummy cell test with the exception of the oven temperature. Temperature is switched off.



Fig. 73 Stabilizing cell current

In the NOISE screen, the cell current is shown and the output voltage.

				NOISE
	V o u t = + 0 . 0 0 7	V I c	= + 2 . 6 6 7 n A	
PR	EV			


To be able to start an internal dummy cell test you need to check the Internal dummy cell checkbox (☑) under the Settings tab (controls section) in the device settings window (see Fig. 75). To run an internal dummy cell test in Dialogue Elite, set all the parameters to the values listed in Table XI and set the cell to OFF. Subsequently, click the 'Send to device' button. Evaluate the cell current in the monitor window. The cell current shown in the monitor window is the current generated with the internal dummy cell. The internal dummy cell current and noise level should fall within the criteria specified on the previous page. Make sure to uncheck the Internal dummy cell option again from the device settings menu after the test.

		1	Auto update
M	ode		litem selector
C	all current	+2.655 [nA]	
A	n. output [V]	0.000	💀 Device settings
Po	otential E [V]	0.80	
A	wiliary E [V]	-0.80	Settings Acquisition EC device IP switch
Te	emperature [°C]	28.2	Parameters Menu's
Δ.	tozero	✓ set	Filter:
	al on/off	lo ser	tem:
	al on/on all enabled		Value:
U.	al eriabled		Value.
			Auto load last used file
			Apply Auto load last used file
			vbbà
			Controls
			Kevlock on Show status outputs Dialogue Elte - service version
			interiock on in interior status outputs 000 - soft dopole
			✓ Logging on ✓ Internal dummy on ✓
			Read from Send to Clos
_			device
	ad from		Jeno to device
Ue	avice		device
Incent	ELITE: COM3	and free days	

13.4 Stop flow test

The stop flow test is a basic test to check if the problems are related to either the flow cell or the rest of our (U)HPLC system. Perform the following steps to execute the stop flow test:


- Switch off the (U)HPLC pump.
- Disconnect the tubing connection from the column outlet (see Fig. 76).

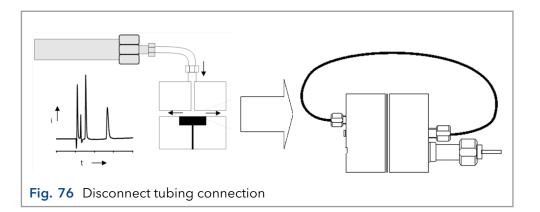


- Disconnect the outlet tubing from the flow cell.
- Connect the other end of the tubing (of the tubing connected to the inlet of the flow cell) to the outlet of the flow cell.
- The fluidic path of the flow cell is now completely isolated from the rest of the LC system.
- Record a run (without injection of sample) to measure/evaluate the background cell current (Icell) and noise.
- Compare the obtained values of the cell current and noise with the values observed before the stop flow with the pump on.

Results:

<u>Icell:</u> In case a significant drop in Icell is observed, for example a drop of more than 50%, it is an indication that the problem is not flow cell related but originates from other parts of the LC system. The most obvious reasons for a high background current are electrochemically active contaminants in the mobile phase, column bleeding, leaking pulse damper. This causes can be systematically eliminated by replacing the mobile phase or disconnecting column or pulse damper etcetera, and reevaluate the cell current.

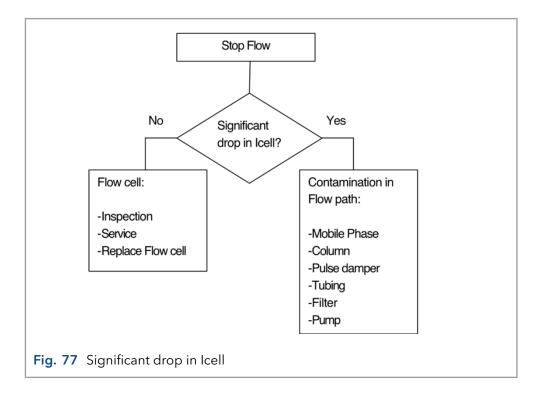

<u>Noise:</u> If a significant drop of the noise only is observed it could be signature of for example pump problem (check valves, air in pump head, compressibility issues or leaking seals).

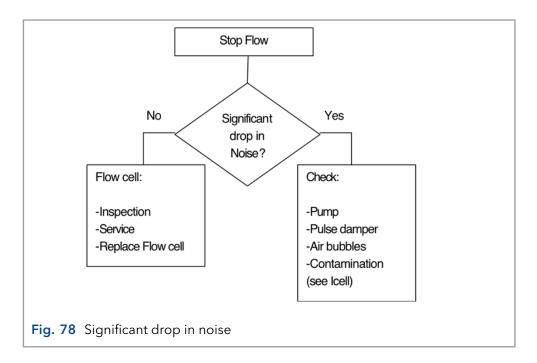

0

In case no significant drop in noise or cell current is observed service or replace the cell. In case you still cannot solve the problem contact your local representative.

Please bear in mind that analytical problems may also be caused by external Influences like temperature, unstable samples. Make sure that the application was running trouble-free before and that no changes have been made to the system. A number of causes and possible solutions for analytical problems is listed below. Contact your local representative if you need further help.

No detector response

Possible cause	Remedy
No power	Check line voltage setting, plug in power cord
Power switch off	Turn this switch ON (at the rear panel)
Faulty fuse	Replace fuse
Divergent mains voltage	Check line voltage
Cell disconnected, or switched off	Check connection
Output disconnected	Check connection
Fouled WE	Clean WE

High cell current

Possible cause	Remedy
Contaminated buffer	Replace buffer, do not recycle the buffer
High WE potential	Optimize potential, if possible: use smaller WE diameter
Salt bridge in REF not saturated	Refill with wetted KCl crystals
Retained peaks from previous runs	Wait for elution of these (very) broad peaks
Column is 'bleeding'	Replace column
High amount of Fe++ in buffer	Add EDTA to buffer, rinse metal parts with 15% HNO ₃

Noisy baseline

Possible cause	Remedy
Salt bridge in REF not saturated	Refill with saturated KCl, add wetted KCl crystals
Air bubble in REF or in cell	Remove air bubble, continuously degas the mobile phase
Slow temperature fluctuations	Isolate detector cell, set oven temperature
Fouled WE	Clean WE
Leaking REF or cell	Tighten connections with care

Decreased sensitivity (low S/N ratio)

Possible cause	Remedy
Fouled WE by dirty samples	Clean WE, if possible: dilute samples
Cell potential too low	Optimize potential
Contaminated buffer (high I _{cell})	Replace buffer, do not recycle the buffer

Saturation of output

Possible cause	Remedy
Damaged REF	Check with spare REF, replace if necessary
Damaged WE	Replace cell block
Cell incorrectly connected	Check connections (REF: black, WE: red, AUX: blue)
Cell potential too high	Optimize cell potential

Baseline oscillations

Possible cause	Remedy
Malfunctioning pump (regular pattern)	Check pump (seals, valves)
Over-tightened cell bolts	Adjust cell bolts, check pump pressure
Air bubbles in cell or REF	Maintenance REF
Temperature oscillations	Set oven temperature
Contaminated buffer (high I _{cell})	Replace buffer, do not recycle the buffer
Fouled WE	Clean WE
Fe++ in buffer	Add EDTA, passivate metal parts with HNO_3

14. Reorders

The electrochemical detector is shipped together with a number of parts. The listing in Table below may not be complete, see check list of delivery for complete listing.

Table XII. Accessory kit A07548-4 (175.0200) for the ECD 2.1 detector

Part number	Component
2.500.040	External dummy flow cell
2.500.107	Column clamp 12 mm
250.0102	Cell clamp kit 40 mm
2.500.113	Fuse 2.5 AT 250 V
2.500.170	LAN (UTP) cable, crossed, 3 meter
2.500.175	USB cable, A-B, 3 meter
250.0131B	I/O conn. board
250.0131C	I/O cable 25M-25M, 1.8 meter
250.0128A	Output cable, D9 male - open, 2 meter
2.500.116	Mains cable (Europe)
2.500.118	Mains cable (USA)
250.0126A	Cell cable

For a VICI Valco electrically-actuated 2-position valve, with an E2CA, EHCA actuator a serial cable is available for control via the ECD 2.1 detector.

For these and other ECD 2.1 parts or flow cells contact your local supplier.

15. Index

Α

ADC 38 ADF 45 Amplitude response 48 ANALOG DATA 8, 75 Applications: quality control 6 Auxiliary electrode 38

В

Biological hazard 6 Buffer pH 21 Buttons on keyboard 28

С

Carbohydrate analysis 52 ClarityChrom® Chromatography software 15 Cleaning of detector 25 Column 23 Compensation 42 Controller 28 Cut off frequency 49

D

DAC output 76 Degassing mobile phase 23 Digital filters 48 Digital I/O connector 8, 78 Door sensor 8 DSP 48 Dual flow cell control 39 navigation 41 Dummy cell external 89 internal 90 test 89

Ε

E1 33 E2 33 EDTA 21 Electrical safety 4 Electrochemical reactions 56 Elite Dialogue software 15 EN 55011 group 1 class A ISM equipment 12 Environmental specifications 9 Errors 87

F

Filter 43 settings 33 Filters low pass noise 47 Flow cell 25 Frequency 45 Fuse replacement 5, 26

Н

Helium degassing 21

I

I/E converter 38, 39 output 77 relationship 56 Inject marker 71 Installation 9 Installation Qualification (IQ) 86 Instrument description 7 Integrator 52 Intended use 1 Ion strength 21 IP address 12

Κ

Keyboard 28

L

LAN connector 8, 70 LC tubing 25 Leakage 25 License dongle ClarityChrom® 17 Dialogue Elite 16, 37

Μ

Maintenance 25 Maximum compensation 42 Messages 87,88 Mobile phase requirements 21

Ν

Noise air bubbles 23 baseline 94 filters 47 increased 24,53 measurement 61 specification 42 suppression 45

0

Offset 43 Operation Qualification (OQ) 86 Output 54 Overload 79

Ρ

Parallel mode detection 40 Passivation with nitric acid 22 pH 21,53 Polarity 43 Potential optimization 56 Power supply fan 11 Pulsed Amperometric Detection (PAD) 52 Pulse mode characteristics 52 controller 30 filter 44 literature references 55 optimization 54 output frequency 54 Pulse settings 53 working electrode 55

R

Range 42 Reference electrode 38 Rotor 53

S

Safety instructions 4 Sampling time 52 Scan mode example 62 scanning voltammogram 57 Serial mode detection 40 Signs 3 S/N ratio decreased sensitivity 94 filter settings 50 Solvents 5 Stop flow test 91 System operation 4

Т

t1 35 Technical data 66 Tefzel rotor 53 Three-electrode configuration 38 Time table 36 Transport box 10 ts (sampling time) 52

U

USB connector 8,69

V

Vacuum degassing 21 Valve configuration electrical 71 manual 72 Valve connector 8 VALVE connector 70 Vespel rotor 53 VICI Valco 22 Voltage clamp 38 Voltammogram construction 61,62 introduction 56

W

Warning symbols 3 Waste disposal 6 WEEE directive 2 Working electrode 38 material 55 Working environment & safety 4

Science Together



Latest KNAUER instructions online: www.knauer.net/library

KNAUER Wissenschaftliche Geräte GmbH Hegauer Weg 38 14163 Berlin Phone: Fax: E-Mail: Internet: +49 30 809727-0 +49 30 8015010 info@knauer.net www.knauer.net

© KNAUER 2019