


Software

► Mobile Control App User Manual

V6851

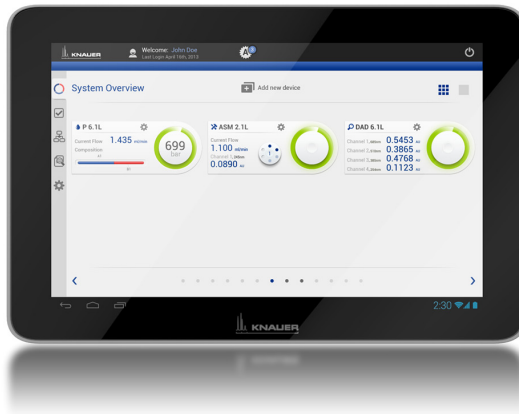

HPLC

Table of Contents

Note: For your own safety, read the manual and always observe the warnings and safety information on the device and in the manual.

Mobile Control	6
Certificate	6
Installing the Mobile Control	7
Demo Mode	7
Full Version	7
Installing the Mobile Control	7
Activating the Full Version	7
License Validity	9
Mobile Control manual	9
Integration into the Network	9
Starting the Mobile Control	9
Exiting the Mobile Control	10
Updating the Mobile Control	10
Changing/Logging-Out User	10
Monitor Mode	10
Operation	11
Prerequisites	11
Entering Values	11
Buttons	11
Network Settings	14
Preferences	15
Instrument Lookup	16
Selecting an device	16
Removing Devices from Selection	16
Device Settings	17
Setting the Leak Sensor	18
Switching the Leak Sensor On/Off	18
Adjusting the Sensitivity of the Leak Sensor	18
Detector	18
Pump	19
Setting the Mixing Chamber	19
Assistant	19
Autosampler	20

Configuration Management	21
Connecting the Router to the Mobile Control	22
Creating the Configuration	23
System Overview	23
Device Status	24
Detail View	24
Events - Setting Short-Circuit Bridges	25
Detector Settings	25
Pump Settings	27
Valve Settings	28
Column thermostat settings	29
Autosampler settings	29
Device Status	29
Programs and Sequences	30
Creating the Program	31
Starting the Program	32
Stopping the Program	32
Creating the Sequence	32
Adding Programs to the Sequence	33
Changing the Program Order in the Sequence	34
Reading Out Program Details in the Sequence	34
Data acquisition	34
Showing the Data Acquisition Window	35
Showing or hiding Traces	36
Zooming by using the Touchpad	38
Zooming by using the Mouse	38
Preview Run	39
Data Viewer	39
Analysing Traces	40
Normalizing Traces	41
Reviewing Program	41
Analysing of Chromatogram	43
Integration Parameters	44
KNAUER Firmware Wizard	45
Installing the Mobile Control	45
Starting the Firmware Wizard	45
RS-232	46
LAN (WIFI)	47

Run Queue	49
Logs and Errors	50
Energy Options	51
Putting Devices into Standby Mode Manually	51
Putting Devices into Standby Mode Automatically	51
Waking Up Devices from Standby Mode Manually	51
Waking Up Devices from Standby Mode Automatically	51
User Management	51
Creating User	52
Changing the Password	52
Changing Own Password	52
Changing the Password of Another User	53
GLP Data	53
Index	56

Mobile Control

The Mobile Control, version V 4.0 or higher, is a program for hand-held devices which operate with Windows 10 or higher. The Mobile Control is used to operate devices of the AZURA product line. As soon as devices are accessed through other software, they cannot be operated via Mobile Control. However, information like the flow rate or the device status continues to be displayed.

Data transfer between the devices and the Mobile Control is usually realized via a WIFI router. Alternatively, the Mobile Control can also be installed on a computer with Windows 10 and access the devices via LAN.

The devices to be operated, the device-specific data as well as the run parameters are recognized and displayed automatically. Different user accounts can be created. It is only possible to switch between users if no run is in progress. Every user profile is password-protected.

Note: This manual only describes the functionalities of the Mobile Control. Note the user manuals of the respective devices.

System requirements	Supported instruments
<ul style="list-style-type: none"> ▪ Windows 10 ▪ Display size: minimum 7", 10" for Data Acquisition ▪ Display resolution: 1024 x 768 ▪ RAM: 1 GB without data collecting, 2 GB with Data collecting ▪ CPU: 1.33 GHz ▪ Processor: Dual-Core ▪ Disc space: Windows: 50 MB without data collecting, 200 MB with data collecting 	Pump P 6.1L Pump P 2.1L Pump P2.1S Pump P4.1S Detector RID 2.1L Detector UVD 2.1L Detector DAD 6.1L Detector DAD 2.1L Detector MWD 2.1L Detector UVD 2.1S Assistant ASM 2.1L Column thermostat CT 2.1 Autosampler AS 3950

Certificate

After a Mobile Control license is purchased, KNAUER provides a certificate containing information about the activation code, serial number of the license and other relevant data relating to your Mobile Control. Please keep the certificate in safe custody.

Installing the Mobile Control

After installing the Mobile Control on your device it can be operated in demo mode or be activated.

Demo Mode

The demo mode offers an overview of the Mobile Control functionalities. You can operate virtual devices with the demo version. After having been downloaded, the Mobile Control can be started by pressing the *TAKE A DEMO TOUR* button.

Full Version

If a tablet PC is delivered, the Mobile Control is installed and activated already. To use your own tablet PC, download the full version from the KNAUER website and install it on your tablet PC.

In order to use the entire range of features of the Mobile Control, a full version has to be purchased and installed.

Please keep in mind that previous license codes (Version 1.x) are not longer valid. For Version 2.x, a new license code must be purchased from KNAUER.

Installing the Mobile Control

Deinstall a previous version of the mobile control, if applicable and delete the data base after storing the previous data at a safe place.

Note: The installer is provided in *DOWNLOADS · SOFTWARE · AZURA MOBILE CONTROL*.

Process

1. Download the installer from the KNAUER website.
2. Execute the installation.

Next steps: Start the Mobile Control and activate the full version.

Activating the Full Version

After the installation is finished, a 20-digit device code is generated. This device code is linked to the MAC address of the network adapter of the hand-held device on which you have installed the Mobile Control.

Note: Before activation, the mobile control will generate a new device code every time the Mobile Control is opened. It is insignificant which of these codes you send to KNAUER together with the serial number of the purchased license, as the Activation Code provided by KNAUER is valid for all Device Codes generated by this end device.




KNAUER generates a license for the device code, which is activated with an activation code. After purchasing a license, KNAUER sends a certificate to you, containing the activation code and all relevant information. Please keep the certificate in safe custody.

If you order a Mobile Control license together with a tablet, KNAUER deliver the configured tablet with activated Mobile Control with certificate. Please keep the certificate in safe custody. If you order a Mobile Control without preconfigured Tablet you will get the Mobile Control certificate with the serial number, but without activation code.

When entering the activation code, include dashes and note the upper and lower case spelling. It might be necessary to switch off the automatic correction of your hand-held device. Store the activation code carefully.

Note: Make sure on your hand-held device that the WLAN is switched on before entering the activation code.

Note: Keep the activation code in safe custody.

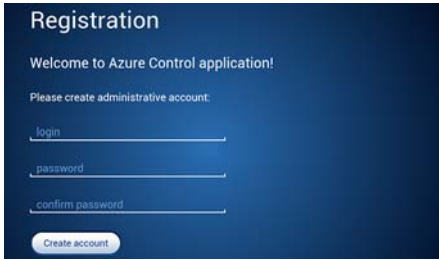Process	Display
<ol style="list-style-type: none"> 1. Tap program symbol. 	 <p data-bbox="564 671 818 699">Fig. 1 Program symbol</p>
<ol style="list-style-type: none"> 2. Submit the 20-digit device code to KNAUER via e-mail or hotline. 3. Enter the activation code sent to you by KNAUER. Pay attention to dashes and upper or lower case spelling. 4. Tap <i>ACTIVATE</i>. 	 <p data-bbox="564 986 962 1013">Fig. 2 Activating the Mobile Control</p>
<ol style="list-style-type: none"> 5. Enter a user name and a password. 6. Confirm the password. 7. Tap <i>CREATE ACCOUNT</i>. An administrator account is created. 	 <p data-bbox="564 1295 972 1323">Fig. 3 Registering the Mobile Control</p>

Next steps: Start the Mobile Control.

License Validity

After the full version has been activated, the license is linked to the MAC address of the hand-held device and cannot be transferred to another device. If the device goes out of operation, one more license may be generated for a new hand-held device. Contact the KNAUER technical support for a new license.

Phone: +49 30 809727-111

Fax: +49 30 8015010

E-mail: mobilecontrol@knauer.net

You can make your requests in English and German.

Re-installing the operating system on the same device has no effect on the validity of the license because the license is linked to the hardware of the device.

Mobile Control manual

The manual is provided on the KNAUER website. When you order a Mobil Control with a tablet, the tablet carries the manual as PDF file.

Integration into the Network

To operate the devices with the Mobile Control, a LAN connection has to be established to the same WLAN router which the Mobile Control communicates with. All devices which are logged into the same network are displayed under *SYSTEM SETUP · INSTRUMENT LOOKUP · DISCONNECTED*.

Starting the Mobile Control

The Mobile Control is started by tapping the icon. The user is prompted to enter the user name and password. If more than one configuration is defined then it is necessary to select the required configuration. For more information please read the chapter on configurations (see "Configuration Management" on page 21).

If you received a tablet with a pre-installed Mobile Control, KNAUER has set up a user account for you already. In this case, the user name is 'Admin' and the password is '12345'. You can find this information on the provided certificate as well. The user name and the password can be changed (see chapter *User Management*).

Exiting the Mobile Control

Process

1. Log out the current user.
2. Tap *CLOSE* ⓘ on the start screen.

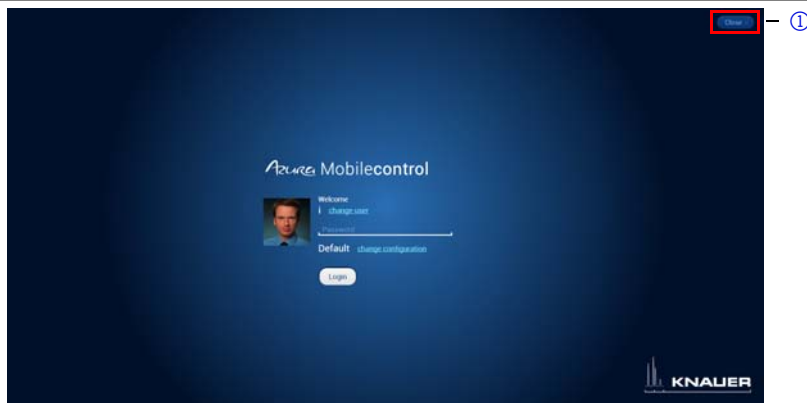


Fig. 4 Closing the Mobile Control

Updating the Mobile Control

Any Mobile Control updates are announced on the website. To update the Mobile Control, download the latest version from the KNAUER website. Mobile Control updates with the same article number are free of charge.

Please deinstall the antecessor version of the mobile control before installing the updated version.

Changing/Logging-Out User

With *LOGOUT*, the user leaves the Mobile Control and all devices are disconnected. The start screen displays the user login. All set data is stored. After that, proceed as described in "Installing the Mobile Control" on page 7.

Monitor Mode

The Mobile Control is automatically set to monitor mode, if a KNAUER chromatography data system (software) accesses the connected HPLC devices. In monitor mode the run parameters of the connected devices can be monitored, but no entry can be made. The Monitor Mode is quit automatically, when OpenLAB is quit.

The latest version of Mobile Control is supported by:

- OpenLAB CDS EZChrom Edition from KNAUER driver package
- ClarityChrom/ClarityChromPrep from KNAUER driver package. Please check release note regarding version compatibility

It is recommended to tap on the *BROWSE* button under *INSTRUMENTS LOOKUP* after changing from Mobile Control mode to monitor mode and after quitting the monitor mode, to refresh the connection of the Mobile Control to the devices.

Operation

APPLY: Numerical input or settings are confirmed with *APPLY*.

CANCEL: Cancels the current process.

The text color indicates the status of the buttons. On active buttons, the text color is blue, on inactive buttons the text color is gray.

Prerequisites

- Hand-held device with Windows 8 (see “System requirements” on page 6)
- WLAN router or LAN router
- The devices are not operated by any other software.
- The full version license has been purchased.
- The Mobile Control has been installed.

Entering Values

Enter values via a slide control ② or via the text input field ① above the slide control. Change the value by tapping the text field/value. If the entered value is above the maximum or below the minimum value, it is corrected to the maximum or minimum value automatically.












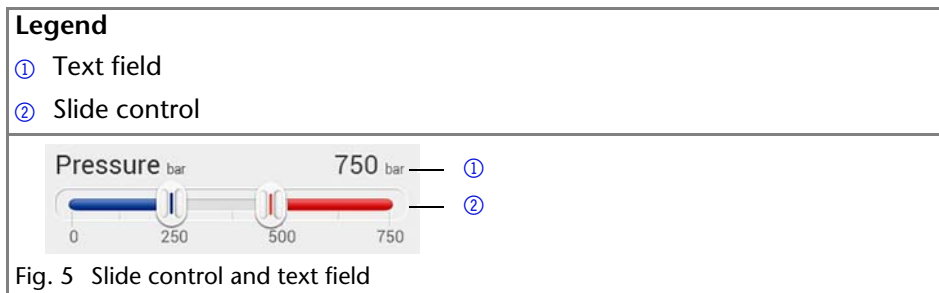


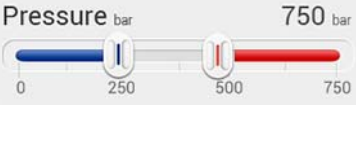


Fig. 5 Slide control and text field

Buttons

The Mobile Control is operated via the touchscreen of a tablet computer or via a personal computer with Windows 8. Submit the set parameters to the respective device with *APPLY*. The following buttons are available:

Control element	Explanation:
	<p><i>SYSTEM OVERVIEW</i></p> <ul style="list-style-type: none"> Shows an overview of all connected devices.
	<p><i>GLP</i></p> <ul style="list-style-type: none"> Displays GLP data for the selected device.
	<p><i>PROGRAMS AND SEQUENCES</i></p> <ul style="list-style-type: none"> List of all saved programs and sequences. Create programs or sequences.
	<p><i>LOGS AND ERRORS</i></p> <ul style="list-style-type: none"> Lists all errors, system logs, and communication logs.
	<p><i>SYSTEM SETUP</i></p> <ul style="list-style-type: none"> <i>NETWORK SETTINGS</i> shows an overview of network settings. <i>PREFERENCES</i> displays basic settings of the devices. <i>ENERGY OPTIONS</i> displays standby mode settings and wakes up devices from standby mode. <i>INSTRUMENT LOOKUP</i> shows an overview of available devices and all devices which are connected. <i>USER MANAGEMENT</i> displays settings and user accounts.
	<p><i>RUN QUEUE</i></p> <ul style="list-style-type: none"> Overview of processable programs and sequences
	<p><i>PROGRAM</i></p> <ul style="list-style-type: none"> Graphical display of the program.

Control element	Explanation:
	<p><i>SYSTEM OVERVIEW, DETAIL VIEW</i></p> <ul style="list-style-type: none"> ▪ Toggles between <i>SYSTEM OVERVIEW</i> and <i>DETAIL VIEW</i>.
	<ul style="list-style-type: none"> ▪ Displays different sets of data depending on the device. ▪ Button displays <i>DETAIL VIEW</i>.
	<p>Error messages</p> <ul style="list-style-type: none"> ▪ Displays error messages.
	<p><i>STOP</i></p> <ul style="list-style-type: none"> ▪ Stops the run.
	<p><i>LOGOUT</i></p> <ul style="list-style-type: none"> ▪ Logs out the user and displays the login screen.
	<p>Button with different functionalities, e.g. <i>RUN</i>, <i>STOP</i>, or <i>STANDBY</i>.</p>
	<p><i>TEXT FIELD AND SLIDE CONTROL</i></p> <ul style="list-style-type: none"> ▪ Slide control sets values. The set value is displayed in the text field. ▪ Enter the numeric value by tapping the text field.

The figure displays an overview of the general buttons:

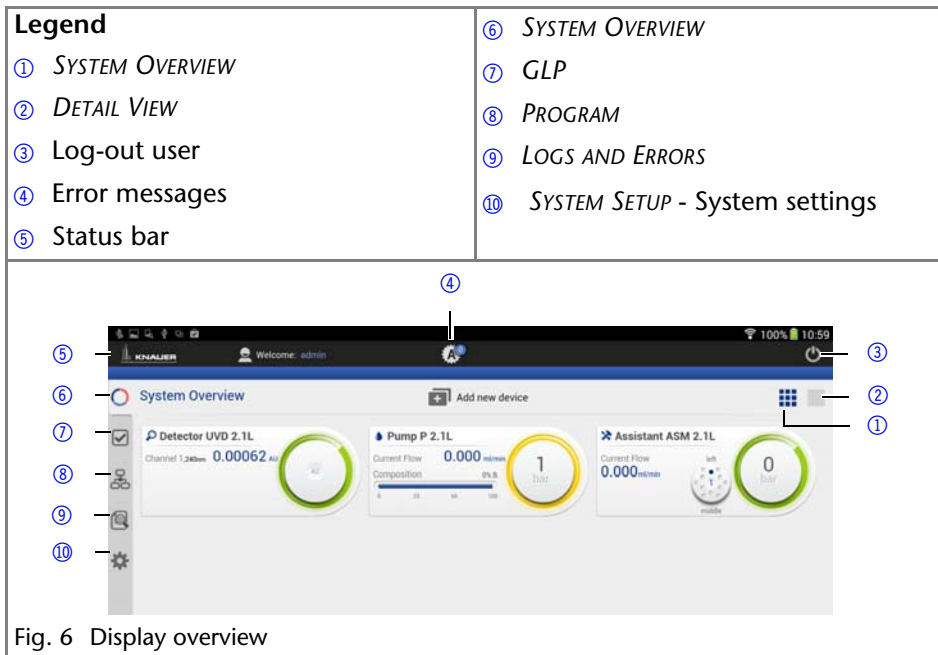


Fig. 6 Display overview

Network Settings

All *INSTRUMENTS LOOKUP BROADCAST PORT*, *IP ADDRESS*, *SUBNET MASK*, and *GATEWAY* data is read-in automatically and does not have to be modified.

NAME: Name of the network.

INSTRUMENTS LOOKUP BROADCAST PORT: Communication in LANs is realized via ports which are part of the network address. If more than one HPLC system is connected to the same LAN and you plan on controlling them separately, you can use different ports to avoid interference. To do so, the port number of every device has to be changed to the same port number in the device configuration of the chromatography software or Mobile Control. We recommend to use the same port number for all devices in the same system.

Note: The port is set to 10001 at the factory. Use identical port numbers in the device configuration of the chromatography software or Mobile Control and in the device, otherwise the connection cannot be established.

ADDRESSING SCHEME: The choice between *DHCP* and *STATIC* is made under *SETUP > GENERAL* for each device. *DHCP* reads-out data like *IP ADDRESS* automatically. If *STATIC* is activated, data is entered manually.

IP ADDRESS: Displays the IP address of the device.

SUBNET MASK: Displays the subnet mask.

GATEWAY: Displays the gateway.

RESET LAN SETTINGS: used to reset communication settings of KNAUER devices for factory defaults (DHCP, 10001 port). A device you want to reset should be on and connected to the same router. To run command you should enter the serial number of the device.

Preferences

SET DATE, SET TIME: The date and time settings refer to the system setups of the hand-held device on which the Mobile Control.

PRESSURE UNITS: The unit for pressure is displayed under *PRESSURE UNITS*. Choose between *bar*, *MPa* and *psi*.

SYSTEM DETECTOR UNITS: The unit for absorption is selected under *SYSTEM DETECTOR UNITS*. Choose between *AU*, *mAU*, μ *AU* (UV detectors), *mRIU*, μ *RIU*, *nRIU* (RI detectors).

TEMPERATURE UNITS: The unit for temperature is selected under *TEMPERATURE UNITS*. Choose between Celsius °C and Fahrenheit °F.

WAIT FOR TEMPERATURE: Choose between *START IMMEDIATELY* and *WAIT*. Here is defined, whether a program starts immediately or after reaching of a programmed target temperature.

SYSTEM LOG: Choose between *SYSTEM* and *COMMUNICATION*. The Mobile Control logs the system protocol/*SYSTEM LOGS* and the communication protocol/*COMMUNICATION LOGS*. The messages are listed under *LOGS AND ERRORS*.

LOG LIFETIME: Choose between 1 week, 2 weeks, and 1 month. *LOG LIFETIME* displays the duration for which the system log or communication log data is stored.

EMERGENCY MODE: This setting specifies how the HPLC devices behave, if the WLAN connection is disrupted. This may happen if the tablet leaved the operating distance of the WLAN router, or the display is switched off. If the setting is set to *CONTINUE*, the devices keep on running. Programs that may run at that moment will continue to the end. If the setting is set to *STOP*, the pumps will stop upon disruption of the WLAN connection, and programs will be canceled.

SHOW VIRTUAL KEYBOARD: This setting is only available with a Windows operating system. The virtual keyboard appears every time a text field is tapped, to make an entry. For tablets without hardware keyboard this setting should always be active. Switching off leads to deactivation of the virtual keyboard not only in the Mobile Control, but for all areas of the tablet. For entries in text fields the virtual keyboard will not appear anymore. If the Mobile Control is installed on a desktop PC with a hardware keyboard, though, it is advisable to switch off the virtual keyboard. Otherwise the virtual keyboard appears each time a text field is clicked on to make an entry, and will overlay the text field, so that the entry is hampered. In addition, all buttons must be clicked on twice to get a reaction.

Instrument Lookup

INSTRUMENT LOOKUP lists all devices which are connected to the router.

BROWSE: Checks all devices which are connected to the router and are available for a configuration.

: Lists all devices under *CONNECTED*.

: Lists all devices under *DISCONNECTED*.

Selecting an device

The Mobile Control is able to connect to one or more devices.

Prerequisite

The Mobile Control has been connected to a router.

All devices which are connected to the selected router, are displayed under *SYSTEM SETUP · INSTRUMENTS LOOKUP*.

1. Tap *SYSTEM OVERVIEW* and *ADD NEW DEVICE* or select *SYSTEM SETUP · INSTRUMENTS LOOKUP · DISCONNECTED* displays all devices which are connected to the router, but not to the Mobile Control. Active devices are displayed in black. Inactive devices, e. g. an device in standby, are displayed in gray.
2. Tapping the button adds the device to *CONNECTED*.
3. Tap *APPLY* to save the settings.

Selected devices are displayed under *SYSTEM SETUP · GENERAL* and under *SYSTEM OVERVIEW*.

Removing Devices from Selection

1. The button removes devices from the configured system.
2. Tap *APPLY* to save the settings.

Device Settings

If a device is connected to the Mobile Control under *INSTRUMENT LOOKUP*, it is displayed under *SYSTEM SETUP · GENERAL*.

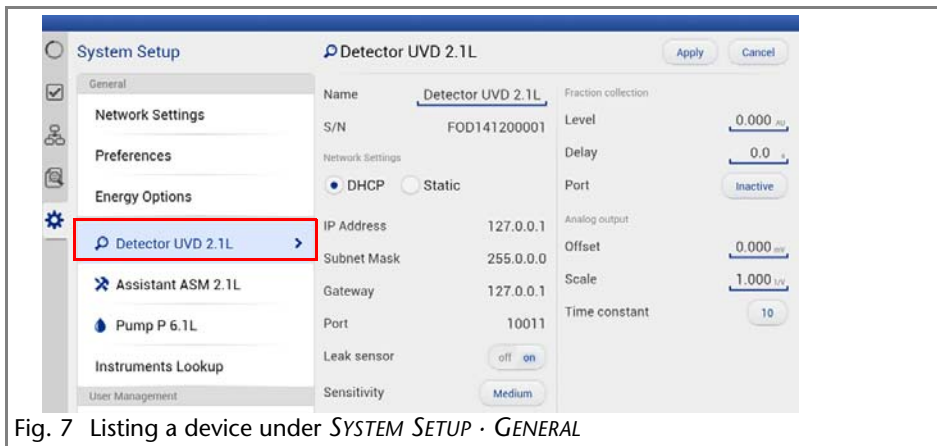


Fig. 7 Listing a device under *SYSTEM SETUP · GENERAL*

NAME: By default, the device designation is used as device name. Tap the name to change it. Change the name if more than one device of the same type is configured in the system.

S/N: The serial number of the device is read out automatically.

PORT: Communication in LANs is realized via ports which are part of the network address. If more than one HPLC system is connected to the same LAN and you plan on controlling them separately, you can use different ports to avoid interference. To do so, the port number of every device has to be changed to the same port number in the device configuration of the chromatography software or Mobile Control. We recommend to use the same port number for all devices in the same system.

Note: The port is set to 10001 at the factory. Use identical port numbers in the device configuration of the chromatography software or Mobile Control and in the device, otherwise the connection cannot be established.

DHCP: DHCP allocates *IP ADDRESS* and *SUBNET MASK* automatically.

STATIC: Static enables you to enter *IP ADDRESS* AND *SUBNET MASK* manually.

IP ADDRESS: Displays the IP address of the device.

SUBNET MASK: An IP address consists of two parts. One part of the IP address designates the network address of the device. The other part designates the distinct address of an device inside of a network. The subnet mask defines which part of the IP address is the network address of an device. It determines which other devices the respective device can communicate with, namely all devices with the same network address. This network is called subnet. This means that all devices of a system and the computer have to operate in the same subnet,

using the same network address. Devices in other networks can only be communicated with via a router.

GATEWAY: If communication has to be established with devices in other networks, a gateway is used. The gateway routes all network requests, which are not directed towards its own network (subnet) to another network (subnet). This task is usually performed by routers which communicate with subnets via IP protocols.

Setting the Leak Sensor

The leak sensor can be switched on and off. Three different settings are available, *LOW* (low sensitivity), *MEDIUM* (medium sensitivity), and *HIGH* (high sensitivity).

Switching the Leak Sensor On/Off

1. Select device in menu *SYSTEM SETUP*.
2. Under *LEAK SENSOR*, activate the button *ON* or *OFF*.
3. The sensitivity of the leak sensor can be adjusted.

Adjusting the Sensitivity of the Leak Sensor

1. Select device in menu *SYSTEM SETUP*.
2. Under *SENSITIVITY*, activate the button *LOW*, *MEDIUM*, or *HIGH*.

Detector

INTEGRATION TIME (DAD): Displays the set integration time in ms.

CELL TYPE (DAD): To enter the installed flow cell choose between *3 mm/2 µl*, *10 mm/2 µl*, *10 mm/10 µl*, *50 mm/6 µl*; *TEST CELL* and *FIBER OPTICS*.

AUTOZERO AT WAVELENGTH CHANGE (DAD): If the function is *ON*, the autozero is carried out automatically at every wavelength change.

ACTIVE CHANNELS (DAD): Activates 1,2,3 or 4 channels.

PORT: Analog output for the signal that triggers e.g. the fraction collection. Choose between *Inactive*, *EVENT 1*, *EVENT 2* and *EVENT 3* depending on the type of detector.

OFFSET: Transfers the analog output signal in mV.

SCALE: Choose between *0 AU/V*, *0.5 AU/V*, *1 AU/V*, *1.5 AU/V*, *2 AU/V*, *2.5 AU/V*, *3 AU/V*, *3.5 AU/V*, *4 AU/V*, *4.5 AU/V* and *5 V*.

TIME CONSTANT (DAD): Smoothes measuring values. Measuring points of a set time interval are combined and the mean value is displayed as a measuring point. A broader interval increases the smoothing proportionally. Choose between *0.00 s*, *0.01 s*, *0.02 s*, *0.05 s* (DAD), *0.1 s*, *0.2 s*, *0.5 s*, *1.0 s*, *2.0 s*, *5.0 s*, and *10 s*.

RESTORE FACTORY SETTINGS: This function enables your to re-set the device to its default settings.

Pump

Availability of options depends on the device type. Unavailable functionalities are grayed out.

PUMP HEAD: Displays the conveying volume of the pump head.

GRADIENT TYPE: Some pumps are pre-configured as isocratic, HPG or LPG pump. In this case, the configuration cannot be changed. For pumps that are not pre-configured, you can select *ISOCRATIC*, *LPG BINARY* or *LPG TERNARY*, *HPG A* or *HPG B*.

START INPUT: Choose *DISABLED*, if you want to deactivate the analogue input. Choose between *START PUMP* and *STOP PUMP*, if the pump should start or stop running upon receiving the trigger signal. Select *ENABLED*, if the trigger signal shall be used to start a programm.

for the analogue input of the pump.

OFFSET (ANALOG OUTPUT): Offsets the analog output signal in V.

FULL SCALE (ANALOG OUTPUT): Choose between 1 V, 2 V, and 5 V.

SIGNAL SOURCE: Choose between *PRESSURE*, *HPG/LPG - A*, *HPG/LPG - B*, *DISABLED*, and *FLOW*.

TIME CONSTANT (Analog output): Smoothes measuring values. Choose between 0.1 s, 0.2 s, 0.5 s, 1.0 s, 2.0 s, 5.0 s, and 10 s.

RESTORE FACTORY SETTINGS: This function enables your to re-set the device to its default settings.

Configuring the Pump for High-Pressure Gradient

For the pumps, which are not configured as HPG automatically, the following settings are made under *SYSTEM SETUP · PUMP*:

1. Set one of the two pumps to *HPG A*.
2. Set the other pump to *HPG B*.

Configuring the Pump for Low-Pressure Gradient

For pumps, which are not configured as LPG automatically, the following settings are made under *SYSTEM SETUP · PUMP*:

Set the gradient to *LPG BINARY* or *LPG TERNARY*.

Setting the Mixing Chamber

The mixing chamber can be set under *SYSTEM SETUP · PUMP · MIXING CHAMBER*. Choose between 50, 100, 200, 350 µl, and a custom size.

Assistant

Depending on the devices built in, the configuration is divided in *LEFT*, *MIDDLE*, and *RIGHT*. Devices are displayed according to device configuration.

D2 LAMP (detector): The lamp can be switched on and off.

PUMP HEAD: Select the size of the pump head. Several sizes are available.

IMIN SENSITIVITY (pump): The motor current is a measure for the current load of the pump and therefore the system pressure for pumps which do not have a pressure sensor. The pump switches off when the current falls below the entered value. Setting for the minimum motor current permitted before the pump switches off (in %).

IMAX SENSITIVITY (pump): The pump switches off when the current exceeds the entered value. Setting for the maximum motor current permitted before the pump switches off (in %).

CONFIGURATION (valve): A list field with different valves is displayed. Choose between *6 PORT MPOS* (multi-position valve), *8 PORT MPOS*, *10 PORT MPOS*, *12 PORT MPOS*, *16 PORT MPOS*, *6 PORT 2 POS*, *8 PORT 2 POS*, and *10 PORT 2 POS*.

Autosampler

Availability of options depends on the device type. Unavailable functionalities are grayed out.

TRAY CONFIGURATION: Opens menu to select used autosampler tray. Choose according to your installed trays

PREP. MODE: Turn on, if autosampler uses Prep Mode

LOOP VOLUME: Enter Volume of installed sample loop volume

TUBING VOLUME: Enter Volume of installed tubing volume

SYRINGE VOLUME: Enter Volume of installed syringe volume

RELAY EVENT: Choose relay event if needed

TRAY POSITION: Choose position of tray. "Home" = backside of tray cabinet, "Front" = frontside of tray cabinet to change vials

SYRINGE POSITION: Choose if syringe needs to be changed

NEEDLE EXCHANGE: Press start to exchange sample needle

Configuration Management

The Configuration Management allows to control and manage different HPLC systems with one tablet. The systems are connected to different routers. Each configuration is linked to the SSID (Service Set Identifier) of the router and contains information of the integrated AZURA devices in the system.

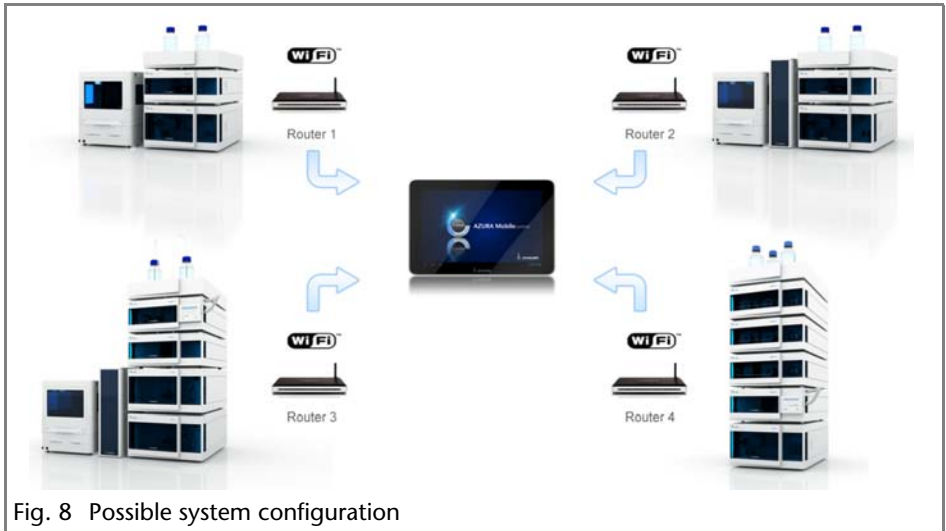
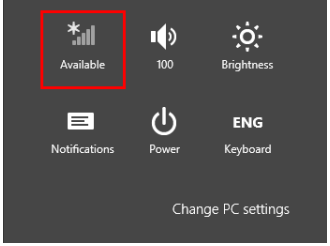



Fig. 8 Possible system configuration

After selecting a configuration the application connects automatically to the corresponding router and enables the communication with the AZURA devices.

Note: Before adding a new configuration the required router must be connected to the application.

Connecting the Router to the Mobile Control

Process	Figure
<ol style="list-style-type: none"> To access the list of available networks from the Start screen, move your cursor to the bottom right corner of the screen and open the charms. <ul style="list-style-type: none"> On a device with touchscreen, swipe from the right side of the screen. Click or tap Settings and then Network. 	
<ol style="list-style-type: none"> Click or tap on the name of the network you wish to connect to. Select "Connect automatically" if you use this network regularly. Windows will automatically connect to the network any time you are within range. 	

Creating the Configuration

A new configuration can be created with the “Create new configuration” button under *SETTINGS > CONFIGURATION MANAGER*. Adding a configuration name and defining the user permission will complete this process.

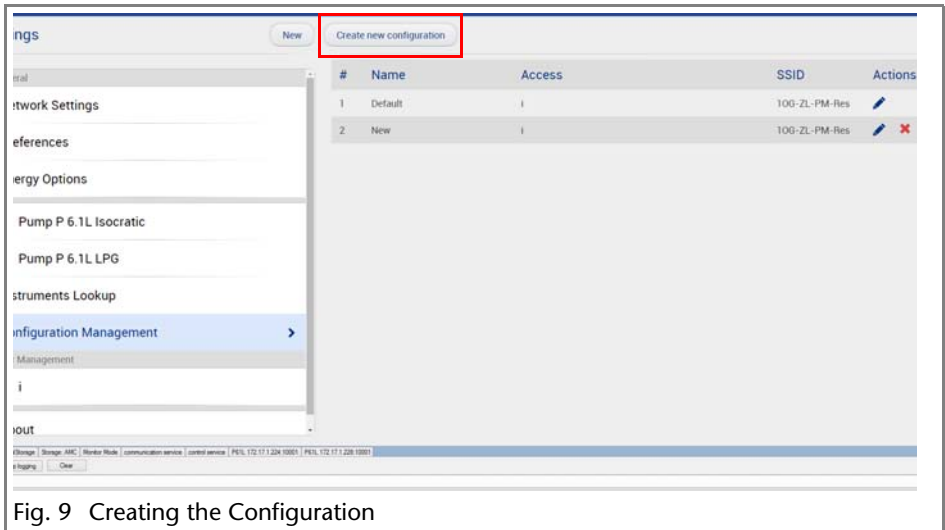


Fig. 9 Creating the Configuration

Name	name of configuration	
Access	users that have permission for controlling the system	
SSID	Service Set IDentifier of connected router	
Actions	edit existing configuration: 	delete new configuration:

System Overview

The system overview offers an overview of all devices which are connected to the Mobile Control at the moment.

Every button displays information about the respective device. The view of an assistant depends on the configuration of the assistant.

Detector	Device name, connection status, wavelength, signal
Pump	Device name, connection status, flow, pressure, the current gradient composition
Valve	Device name, valve position

Column Thermostat Device name, connection status, target temperature, lower and upper temperature safety limit

Autosampler Device name, connection status, vial, temperature

Settings like flow or valve position are made in the *DETAIL VIEW* of the respective device. If a degasser is integrated into an assistant, it is not displayed as a device. The degasser cannot be operated with the Mobile Control.

Device Status

The connection status of the device is indicated in *SYSTEM OVERVIEW* with four different colors:

	green	yellow	red	blue
Detector	The detector is now ready to operate.	The lamp is turned off or the detector starts.	The device transmits an error.	The device is in standby mode.
Pump	The pump conveys.	The pump is ready for operation.	The device transmits an error.	The device is in standby mode.
Assistant with detector and/ or pump	The detector is ready and/ or the pump conveys.	The lamp is turned off or starts or the pump is ready for operation.	The device transmits an error.	The device is in standby mode.
Column thermostat	The column thermostat is heating or cooling or has reached the target temperature.	The column thermostat is ready for operation.	The device transmits an error.	
Autosampler	The Autosampler is ready for operation.		The device transmits an error.	

Detail View

Via *SYSTEM OVERVIEW* · *DETAIL VIEW* or by tapping on the widget in the *SYSTEM OVERVIEW* you can access the *DETAIL VIEW*. Information and settings for the current device are displayed here.

Some devices provide additional settings on further pages. To browse to the next page, tap the < or > (arrow) ④ in the center of the screen.

If more than one device is installed, you can browse through the devices by tapping the arrows < and > ①, at the bottom.

Legend

- | | |
|---|---|
| <ul style="list-style-type: none"> ① Toggles between the devices ② Different buttons are displayed depending on the device.
Run = starts the run
Stop = stops the run,
Autozero = performs an autozero ③ <i>STANDBY</i> = standby mode
<i>POWER UP</i> = wake up from standby mode | <ul style="list-style-type: none"> ④ Further settings for this device ⑤ <i>DEVICE STATUS</i> - window for the status of the device ⑥ Color display device status ⑦ <i>CANCEL</i> = Cancel settings ⑧ <i>APPLY</i> = Apply settings |
|---|---|

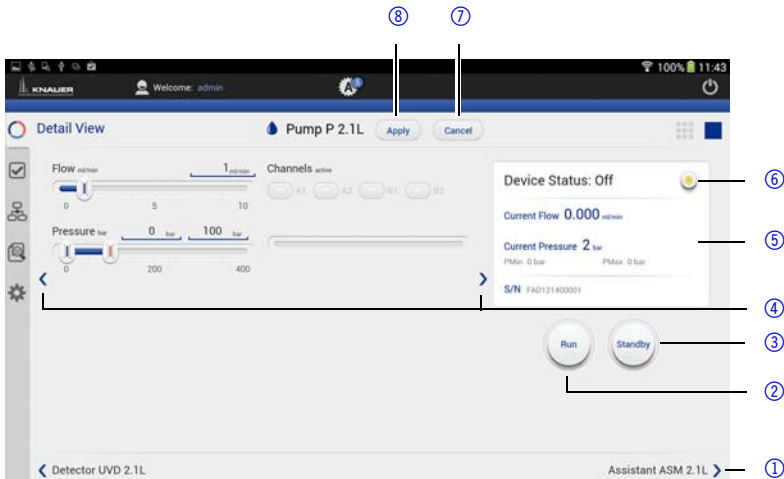


Fig. 10 Detailed view of the pump

Events - Setting Short-Circuit Bridges

Events are short-circuit bridges which can be programmed or manually activated. They operate external devices.

Short-circuit bridges are set under *DETAIL VIEW · EVENTS*.

Note: Note the information about events in the user manual of the respective device.

1. Tap the drop-down list for *ON*, *PULSE*, or *OFF*.

Detector Settings

The following section describes the detector settings.

Entering the Wavelength

The wavelength is entered under *DETAIL VIEW · WAVELENGTH*.

1. Tap the text field and enter the value or adjust the value with the slide control.
2. Save data with *APPLY*.

Entering Reference Correction

A baseline correction can be activated for channels 1 and 2 under *DETAIL VIEW · REFERENCE CORRECTION*.

Signal Mode

Touch *DIRECT* or *INVERTED* to change the signal.

DIRECT SIGNAL(+): Displays the signal without modifications.

INVERTED SIGNAL (-): Inverts the signal.

Switching the Lamp On/Off

The lamp of the detector can be switched on and off.

DEUTERIUM LAMP ON/OFF, HALOGEN LAMP ON/OFF: The lamp is switched on or off using *ON* oder *OFF*. The text color displays the current status of the lamp. On an active button, the text color is blue.

Checking the Wavelength Accuracy

By touching *VALIDATE* or *CALIBRATE* the device checks the wavelength accuracy.

Performing an Autozero

Tap *AUTOZERO*. The *DEVICE STATUS* displays 0.00000 AU after an autozero.

If the function *DEVICE SETTINGS · AUTOZERO AT WAVELENGTH CHANGE* is activated, the device is carried out an autozero at every wavelength change. By default all detectors carry out an autozero at the start of a run.

Flushing the Detector

RID 2.1L only. The flush function activates the reference cell valve enabling this cell to be purged with eluent. The flush valve can be switched on and off immediately either via software or via analog command; alternatively via software a flush time program can be selected, whereby the valve is switched on and after a selected time span (30 s, 60 s, 120 s, 400 s) the valve is automatically switched off. The flush time program can be interrupted at any time with the off command.

Target temperature

RID 2.1L only. It is possible to select the temperature of the optical unit in the range 30-55 °C in 1 °C steps via software. It is recommended to set the temperature 5-10 °C above the ambient conditions, in order to improve and ensure baseline stability.

Pump Settings

The following section describes the pump settings.

Setting the Flow Rate

Set the flow under *FLOW* and confirm with *APPLY*. The pump starts running. The pump can be stopped again with the *STOP*-button in *DETAIL VIEW*.

Setting the Pressure

Set the minimum and maximum pressure under *PRESSURE* and confirm with *APPLY*. Minimum and maximum pressure is set with 2 slide controls or via the text field.

Minimum pressure: The pump switches off after 30 seconds, if the pressure goes below the minimal pressure limit. This may be the case, if a leak is detected or air bubbles are in the system.

Maximum pressure: To protect the column, the pump switches off immediately, if the pressure exceeds the maximum pressure limit, e.g. in case of clogging or an excess flow rate.

Gradients

If a gradient-compatible pump or pump combination has been connected, the gradient can be set under *DETAIL VIEW*. Some pumps have to be configured as gradient pump under *SYSTEM SETUP > PUMP* first (see chapter *Device settings > Pump*).

Setting the High-Pressure Gradient

Under *DETAIL VIEW*, the settings for the composition of the gradient are made.

Some pumps offer a solvent selection option. For these pumps, a selection can be made for every channel between two different solvents.

Pump A: Channel A1 or A2

Pump B: Channel B1 or B2

1. Select the required channels by tapping.
2. Tap the text field to enter the amount of solvent in %. The channel's share of the gradient is indicated by the color of the bar.

Setting the Low-Pressure Gradient

Under *DETAIL VIEW*, the settings for the composition of the gradient are made.

The active channels are shown as colored dots. For pumps, that are pre-configured as LPG pump, channels can be activated or deactivated here separately by tapping. For pumps, that are not pre-configured as LPG pump, this setting can be made under *SYSTEM SETUP > PUMP* by choosing between *LPG BINARY* or *LPG TERNARY*.

Tap the text field to enter the amount of solvent in %. The channel's share of the gradient is indicated by the color of the bar.

Choosing the solvent factor

Use the arrow > in *DETAIL VIEW* to switch to further settings. Select the solvent for each channel from the drop-down list under *SOLVENT A, B, C, or D*. The factors of compressibility for the listed solvents are provided. For other solvents tap the text field in last row of the list and enter the factor of compressibility for a user defined solvent.

Purge

To remove air from the pump head or to change the solvent, purge the pump with *PURGE*.

1. Open the venting screw at the pressure sensor to prevent a pressure surge and damage to the column.
2. Use the arrow > to move to other settings.
3. Enter the flow under *PURGE*.
4. Tap the button *ON* and confirm with *APPLY*.

PMax Mode

The P_{Max} Mode is available for some pumps and enables settings which determine how the pump reacts when maximum pressure is reached.

If the P_{Max} Mode is activated, the pump continues to run with the set pressure. The flow is adjusted in order to keep the pressure. A maximum flow can be set under $P_{Max} \text{ MODE} \cdot \text{MAX. FLOW}$.

At the end of a run, the flow is gradually decreased to prevent a sudden drop in pressure.

1. Use the arrow > to move to other settings.
2. Tap the button $P_{Max} \text{ MODE}$. A window is opened.
3. Tap *ON*, enter a value for *MAX.FLOW* and confirm with *OK*.
4. Confirm with *APPLY*.

Isobar Mode

The Isobar Mode will be available in a later version.

Starting Pump

To start the device, open the *DETAIL VIEW* and tap *RUN*.


Valve Settings

The following section describes the valve settings.

Choosing a Valve

The valve head is recognized automatically and displayed in the menu *SYSTEM SETUP · ASSISTANT · CONFIGURATION*. For 6-port valves, the valve has to be set as a 2-position valve or a multi-position valve under *SYSTEM SETUP · DEVICE SETTINGS*.

Changing the Valve Position

Process	Figure
<ol style="list-style-type: none"> 1. Select <i>DETAIL VIEW</i>. 2. Insert the corresponding valve position ①. The selected valve position is displayed as a number in the center of the graphic. 3. Tap <i>APPLY</i> to save the settings. 	 <p data-bbox="566 363 941 394">Fig. 11 Changing the Valve Position</p>

Column thermostat settings

The following section describes the column thermostat settings.

Setting the target temperature

Set the target temperature under *TARGET TEMPERATURE* and confirm with *APPLY*.

Setting temperature safety limits

Set the minimal and maximal allowed temperature under *TEMPERATURE SAFETY LIMITS* and confirm with *APPLY*.

Autosampler settings

The following section describes the autosampler settings.

Start Wash

The autosampler will start a wash cycle to wash the cycles.

Setting the target temperature

Set the temperature under *TARGET TEMPERATURE* and confirm with *APPLY*.

Move Tray

Move tray to front or to back to enter or remove vials.

Device Status

Under *DEVICE STATUS*, the most important information about the device is displayed. Depending on the device, the type of information may differ.

General

S/N, the serial number of the device. For the assistant, the serial number of the assistant is displayed and not the serial number of the installed devices. The serial number of the installed devices is displayed under *GLP data*.

Detector

Wavelength

Pump

- Current Flow

- Current Pressure (valid for devices with pressure sensor)
- Minimum and maximum pressure (valid for devices with pressure sensor)

Valve

Current valve position

Assistant

Depending on the configuration of the assistant, information about the respective device, as described above, will be displayed.

COLUMN THERMOSTAT

- Target temperature, the entered target temperature
- T Min, the minimum selectable temperature for this thermostat
- T Max, the maximum selectable temperature for this thermostat
- Temperature safety limits
- Measured temperature, the current temperature of the column thermostat

AUTOSAMPLER

- Device Status
- Current Vial
- Serial Number
- Target Temperature
- Current Temperature

Programs and Sequences

Under *PROGRAMS AND SEQUENCES*, a list of saved programs and sequences is displayed. Programs and sequences can be saved, loaded, edited, copied, or deleted.

Tap the program or sequence to display the settings.

Legend

- | | |
|-------------------------------------|-----------------------|
| ① SAVE - Save program | ⑤ Delete program line |
| ② CANCEL - End input without saving | ⑥ Add device or event |
| ③ Name of the program | ⑦ Enter value |
| ④ Add program line | ⑧ Select parameter |
| | ⑨ Select device |
| | ⑩ Enter point of time |

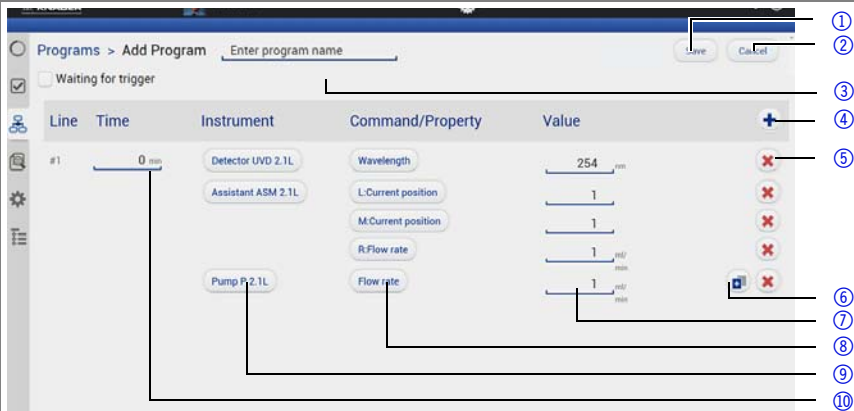


Fig. 12 Programs

+ **ADD**: Adds a new program or sequence.

LOAD: Loads the highlighted program or sequence and starts running the first line of the program to adjust the devices to the starting parameters. Pumps start running.

EDIT: Enables editing of the program.

COPY: Copies the program. This program or sequence can only be saved under a new name.

DELETE: Deletes the highlighted program or sequence.

Creating the Program

Open the window for creating programs under *PROGRAMS* and + **ADD**. Which input can be entered into list fields or text fields depends on the device and is displayed accordingly.

Note: If "Waiting for trigger" is activated, the program can be started with an external trigger.

Process

1. Enter the name of the program into the text field ② .
2. Set the time under *TIME*. To change the time, tap the text field ⑦ and enter the time. The first program line always begins with 0 min.
3. Under *INSTRUMENT*, tap the button ⑥ and select a device.
4. Under *COMMAND/PROPERTY*, tap the button ⑤ and select a parameter.
5. Enter a value into the text field ④ . The displayed value depends on the selected device.
6. Add further program rows ③ .
7. Save the program with *SAVE* ① .

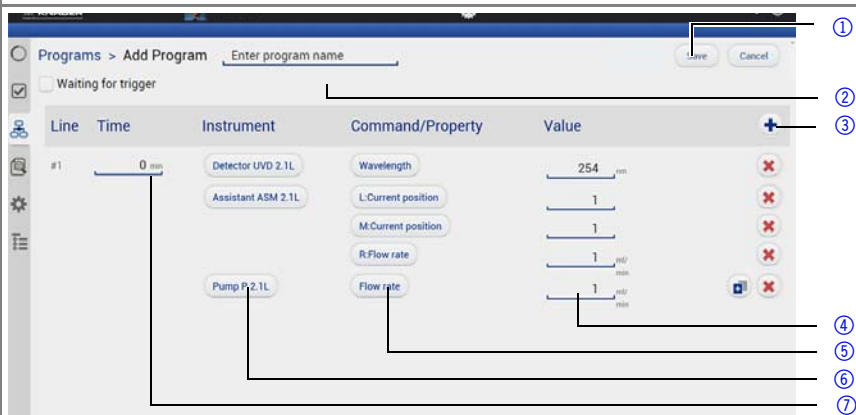


Fig. 13 Creating a program

Starting the Program

1. Select program.
2. Tap *LOAD* to load the program. The first row of the program is executed immediately and settings are kept until the program is started.
3. Tap *START* in the status bar to start the program.

Stopping the Program

Tap *STOP* in the status bar to end the program. Pumps keep running and must be stopped manually under *DETAIL VIEW*.

Creating the Sequence

Open the window under *PROGRAMS* and + *ADD*. Which input can be entered into list fields or text fields depends on the device and is displayed accordingly.

Note: Tapping the button *CANCEL* ends the input without saving.

Process

1. Enter the name of the sequence into the text field ② .
2. Add a program ① .
3. To save the sequence, tap the *SAVE* button ③ .

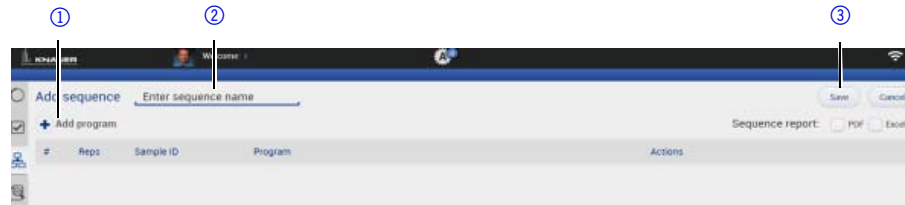



Fig. 14 Add program to a sequence

Adding Programs to the Sequence

Process

1. To edit a sequence, select the sequence line and tap  .
2. Tap the + *ADD* button 14.
3. Tap the *SELECT PROGRAM* ① button.
4. Select the required program and press OK.
5. Enter a Sample ID ② and add the appropriate suffix (e. g. increment number, user name, method name, instrument name, date and time).
6. Enter a Filename ③ and add the appropriate suffix.
7. Enter the number of repetitions ④ .
8. Enter a required comment ⑤ .
9. To save the sequence, tap the *SAVE* button ⑥ .

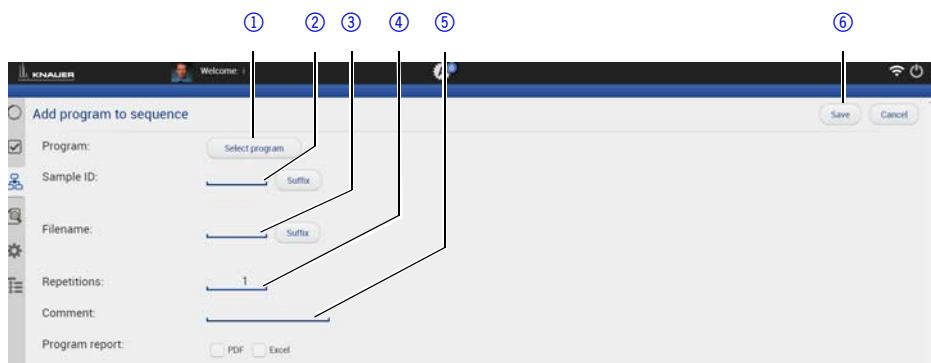




Fig. 15 Adding programs

Changing the Program Order in the Sequence

Process

1. To edit a sequence, select the sequence line and tap .
2. Change the order using the arrows .

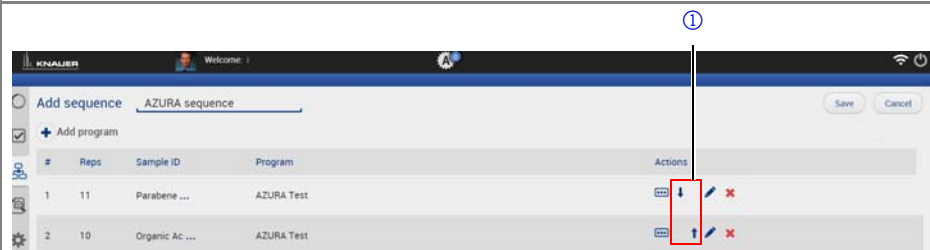


Fig. 16 Changing the program

Reading Out Program Details in the Sequence

Process





1. To edit a sequence, select the sequence line and tap .
2. To read out the program details, tap on the symbol  .
3. To save the sequence, tap the *SAVE* button .



Fig. 17 Reading out program

Data acquisition

The data acquisition tool allows to acquire online data of detectors. Besides detector signals auxiliary traces from AZURA devices can be acquired. Auxiliary traces can be pressure signals from pumps, temperature traces from detectors, autosamplers and column thermostats.



Fig. 18 The data acquisition pool

Showing the Data Acquisition Window

After loading and starting a program the data acquisition window is available. At the end of the program the data file will be stored in the project folder. All data will be stored in **C:\Knauer\Project folder\Result\Data**

Process

1. To select the data acquisition window, tap on symbol ①.


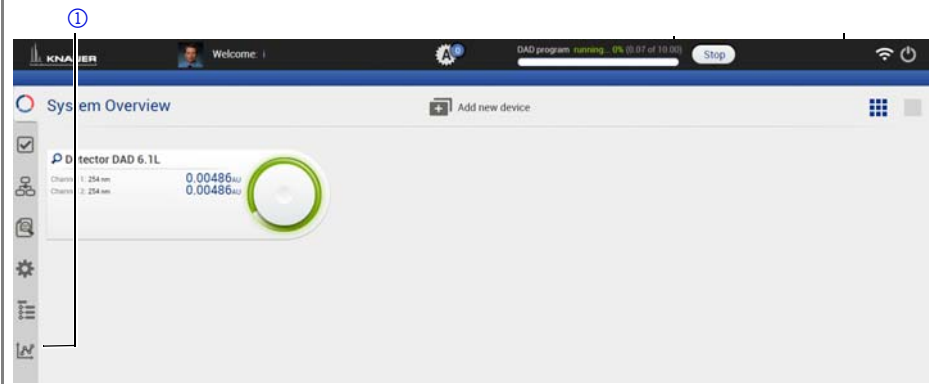


Fig. 19 Selecting the data acquisition window

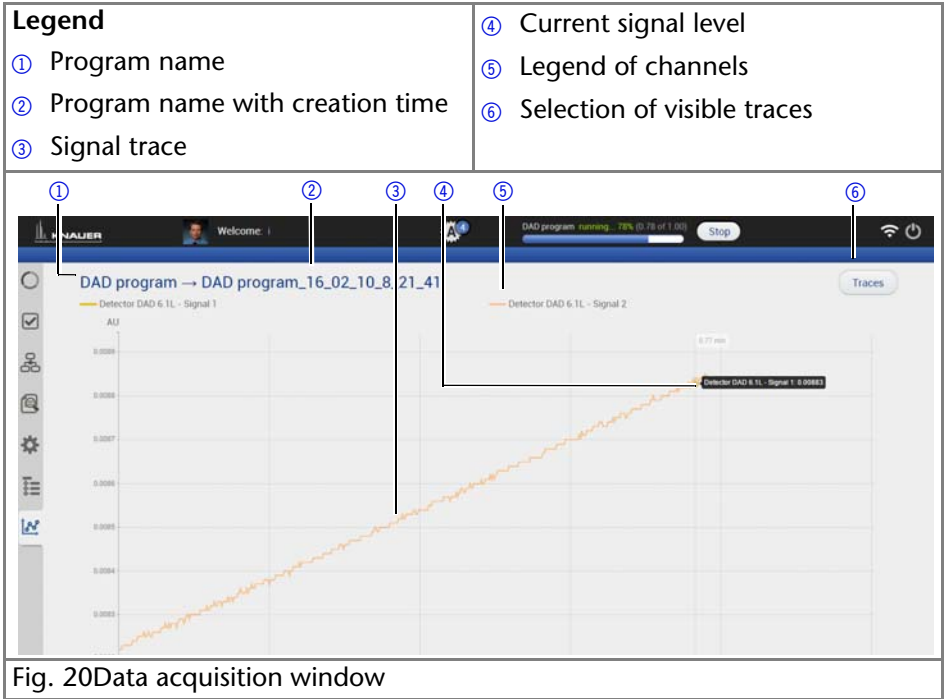

Showing or hiding Traces

During a running program or at the end traces can be shown or hidden. In the list of traces all available data traces, auxiliary traces and program parameters are shown. Program parameters are marked with light grey. This indicates that these traces can be not selected as active trace. The active trace can be selected.

Process

1. To show or hide the traces, tap the *TRACES* button ①.
2. To make this trace visible or hidden, tap the *OFF/ON* button ②.
3. To select the active trace, press the trace name. The active trace is highlighted with bold face ③.

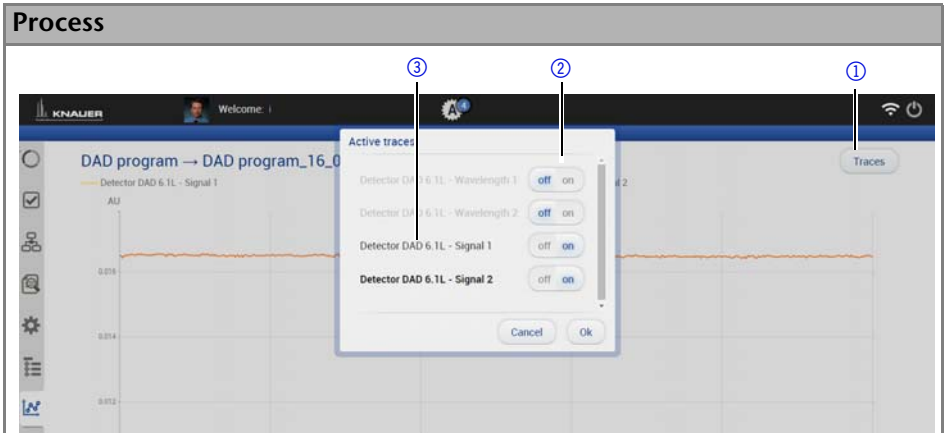


Fig. 21 Showing traces



Fig. 22 Selecting traces

Zooming by using the Touchpad

Process

1. To move the data traces, touch the surface with one finger and moving the finger in the required direction.
2. Use gesture ① to zoom in into the x-axis of the data trace.
3. Use gesture ② to zoom out of the signal trace.
4. By double click on the screen the data trace will be fully unzoomed.



Fig. 23 Zooming

Zooming by using the Mouse

Process

1. The data trace can be moved by moving the mouse with pressed left mouse button.
2. Scroll wheel up/down + ALT: Zoom in/out y-axis
3. Scroll wheel up/down + CNTR: Zoom in/out X-axis
4. By left double click the data trace will be fully unzoomed.

Preview Run

After creating of a new program or modifying of an existing program a preview run can be started. In this run the first program line of all devices will be executed. This allows to check if all devices are running as expected or can be used for monitoring the equilibration of the HPLC system. The preview runs for an infinite time and must be stopped manually. During a preview run signals from detectors or auxiliary traces will be acquired. This data cannot be stored and is only available until a new preview or program/sequence is started.

Process

1. Open the program window, by tapping on symbol ①.
2. To start a Preview run of an saved program, tap the *PREVIEW RUN* button ②.

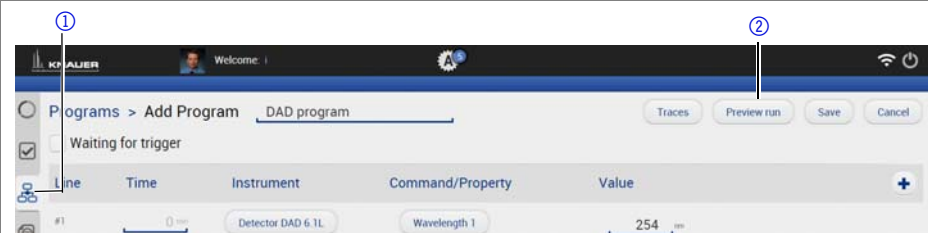


Fig. 24 Program window

Data Viewer

For reviewing acquired data and analysing chromatograms Data Viewer program must be started. After opening all the available traces are overlaid.

Process

1. Start the program "KnauerDataViewer".
2. To open a file, tap on the *CHOOSE FILE* button ①.
3. Select the data file and press the *OK BUTTON* ② to open the data file.



Fig. 25 Choosing file



Fig. 26 Confirming

Analysing Traces

Select the trace you want to analyse. After selecting the active trace this trace will be automatically analysed. The limits for the y-axis for each trace can be adapted.

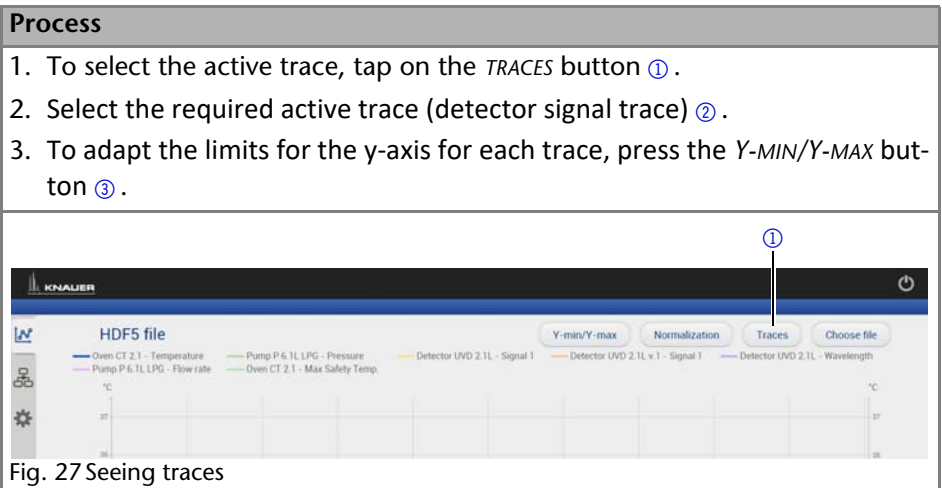


Fig. 27 Seeing traces

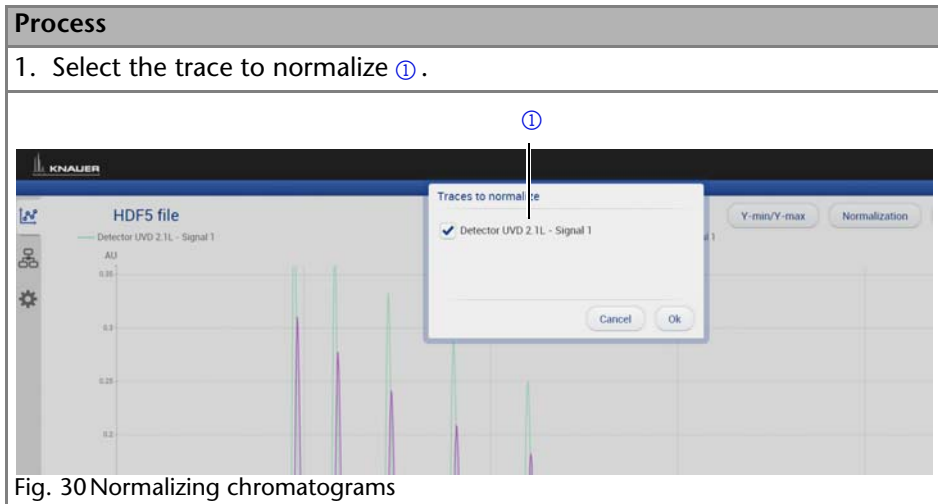


Fig. 28 Selecting traces



Normalizing Traces

Traces can be normalized to other traces. This function allows you to normalize one or more chromatograms to the first chromatogram, adjusting the heights such that the apex height of a selected peak matches that of the peak selected on the first trace.



Reviewing Program

After opening of a data file the method which is related to this program can be reviewed by clicking program symbol.

Process

1. To review a method, tap on symbol ①.

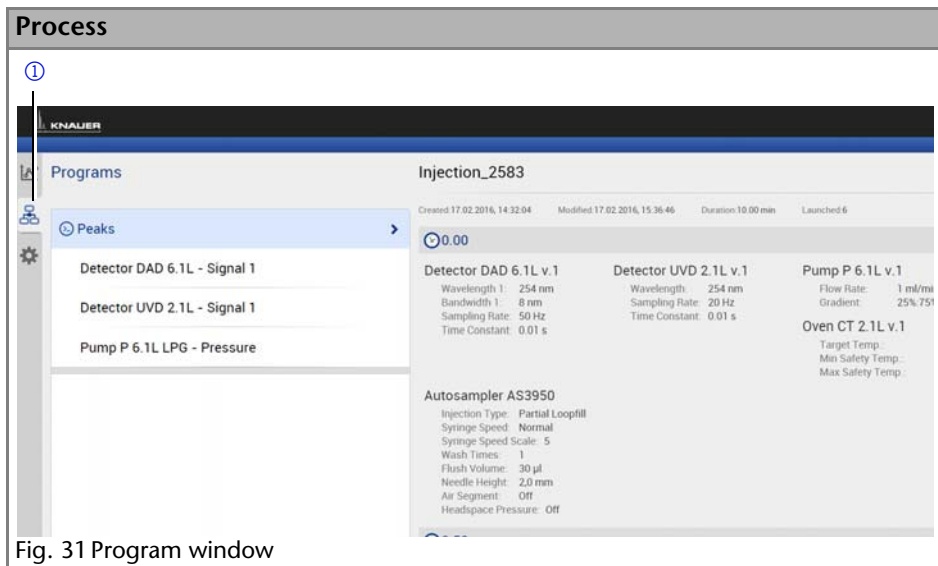


Fig. 31 Program window

Appearance

The appearance of the chromatogram can be adapted in the settings window. The following check boxes are available:

- Grid: Activate or deactivate
- Baseline: Hide or unhide
- Fill peak: The area below the integrated peak can be filled with solid color
- Show second Y-axis: Define an additional axis on the right side of the chromatogram

For each peak two different annotations can be defined. The following annotations are available:

- Peak number
- Retention time
- Peak Area
- Peak Area [%]
- Height
- Height [%]
- Asymmetry
- Platen numbers
- Resolution
- Width

Process

1. To review a method, tap on symbol ① .
2. Choose the appearance by using the check boxes ② .
3. Choose the annotations by using the *WIDTH* and *PEAK AREA* buttons ③ .

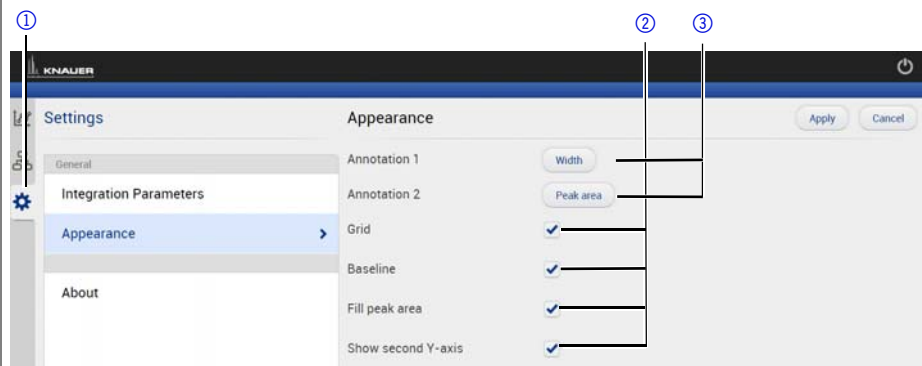


Fig. 32 Setting window

Analysing of Chromatogram

Each detector signal can be analysed. It is possible to define separate integration parameters for each trace. The Data Viewer will determine values for Peak number, Retention time, Peak Area, Peak Area, [%], Height, Height [%], Asymmetry, Plate numbers, Resolution, Width.

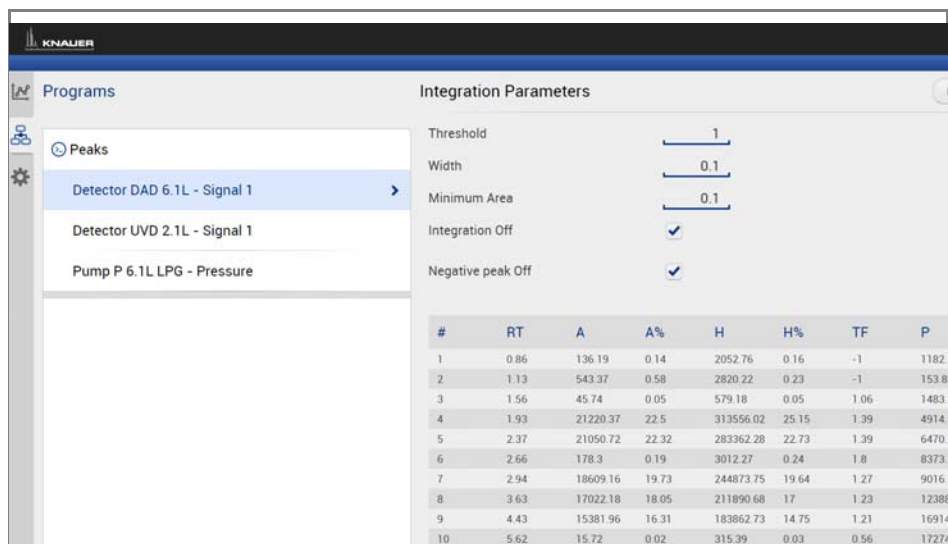


Fig. 33 Analysing of chromatogram

Integration Parameters

Two Integration events are required for each run: Width, and Threshold. These parameters are used to detect peak start, stop, and apex, and to distinguish true peaks from noise.

Width

The Width is used to calculate a value for smoothing, the data points before the integration algorithm is applied. In most circumstances, an initial Width value based on the narrowest peak in the chromatogram will be adequate for proper integration of all peaks.

Threshold

This parameter is the first derivative, used to allow the integration algorithm to distinguish the start and stop of peaks from baseline noise and drift. The recommended Threshold value is based on the highest first derivative value determined in that section of the chromatogram.

Minimum Area

The Minimum Area parameter is used to reject unwanted peaks in the chromatogram. A value of e.g. 1000 will omit all peaks with an area of smaller than 1000.

After entering of integration parameters and pressing the apply button the chromatogram will be re-processed and analysed. The integration table will be updated automatically.

Export

The result table for all traces can be exported into an ASCII-file. The file is stored in

C:\Users\USERNAME\AppData\Local\Azura\Knauer\Export

This file can be used to create a calibration curve for calculation the amount of an analyte in a sample.

In case of changing the integration parameters and re-analysing the chromatogram, the new resulting values will be added to this ASCII-file.

KNAUER Firmware Wizard

This chapter contains detailed information on how to perform an update of all possible firmware components for the various devices using either RS-232 or LAN interface.

Installing the Mobile Control

De-install a previous version of the Firmware Wizard, if applicable.

Process

1. Download the installer from the KNAUER website.
2. Execute the installation.



Fig. 34 Program window

Note: After the installation a re-start of the computer or tablet is recommended.

Starting the Firmware Wizard

Process

1. The Firmware Wizard is started by tapping the symbol



The Firmware Wizard supports the firmware update of the following AZURA devices:

1. AZURA RID 2.1L
2. AZURA ASM 2.1L
3. AZURA DAD 6.1L
4. AZURA DAD 2.1L
5. AZURA MWD 2.1L
6. AZURA UVD 2.1L
7. AZURA UVD 2.1S
8. AZURA P 6.1L
9. AZURA P 2.1L
10. AZURA P 2.1S/P 4.1S
11. CT 2.1

Note: The firmware update of other KNAUER devices (Smartline, BlueShadow) is possible but not fully supported. In case of issues please contact KNAUER.

RS-232

This chapter contains information on how to browse for the various devices using either RS-232 or LAN interface and how to establish connection with a particular device.

- Interface -> RS-232
- COM Port -> Select the COM-port to which the device is physically connected
- Baud rate -> Select the baud rate (has to correspond with the baud rate in the device)
- Connect -> Connects the device with the Firmware Wizard
- Disconnect -> Disconnect the device from the Firmware Wizard

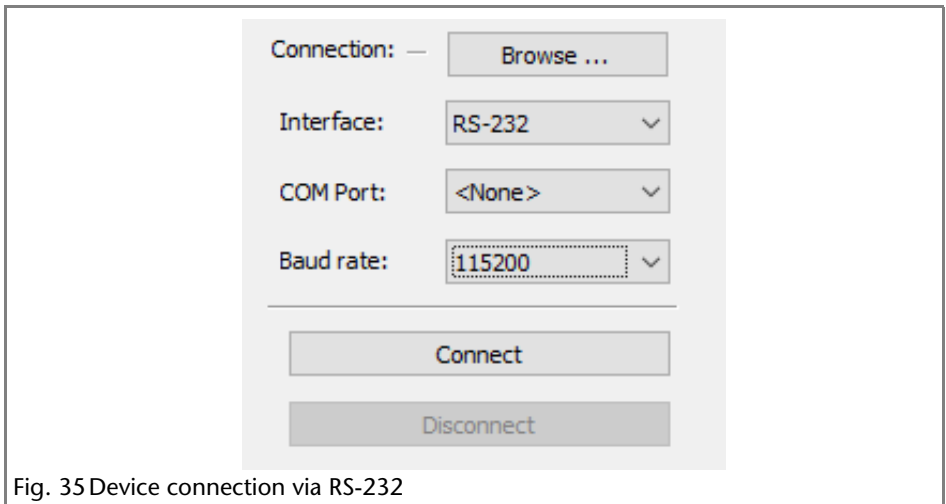


Fig. 35 Device connection via RS-232

LAN (WIFI)

- Interface -> LAN

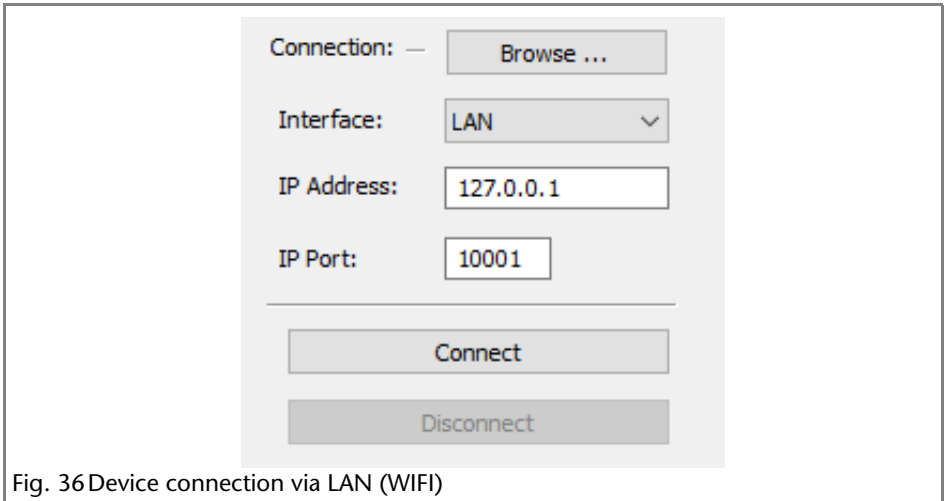


Fig. 36 Device connection via LAN (WIFI)

Process

1. The **Browse** button opens the device selection window.
2. Select the device and press the **SELECT** button.

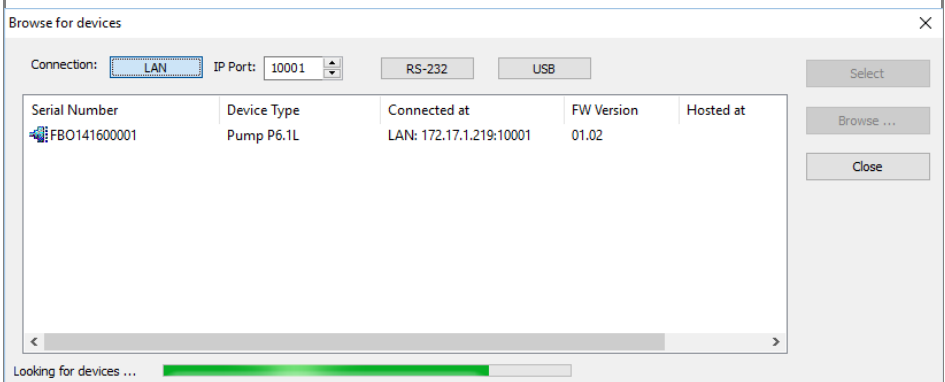


Fig. 37 Program window

3. Press the **CONNECT** button to establish a connection between the computer and the device.
4. Press the **BROWSE** button (**...**) next to the device line (e.g. HPLC Pump P 6.1L) to select the firmware package file. The file has the suffix .fwxxx.

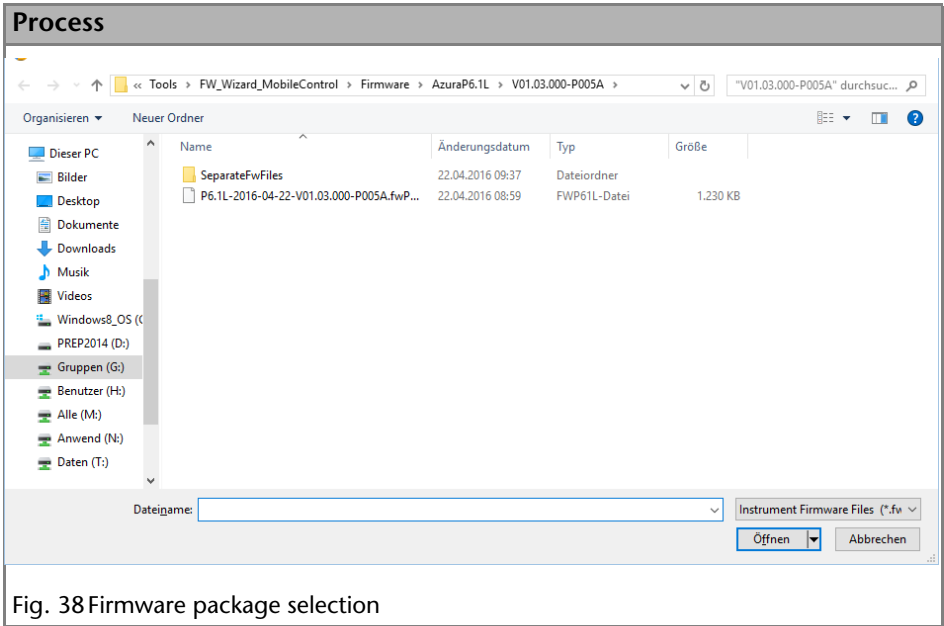


Fig. 38 Firmware package selection

Note: The The Service Tool automatically ticks the checkbox next to “Do not update this component” if the chosen firmware of the component is older than the firmware in the device. In this case the firmware of this component will not be updated.

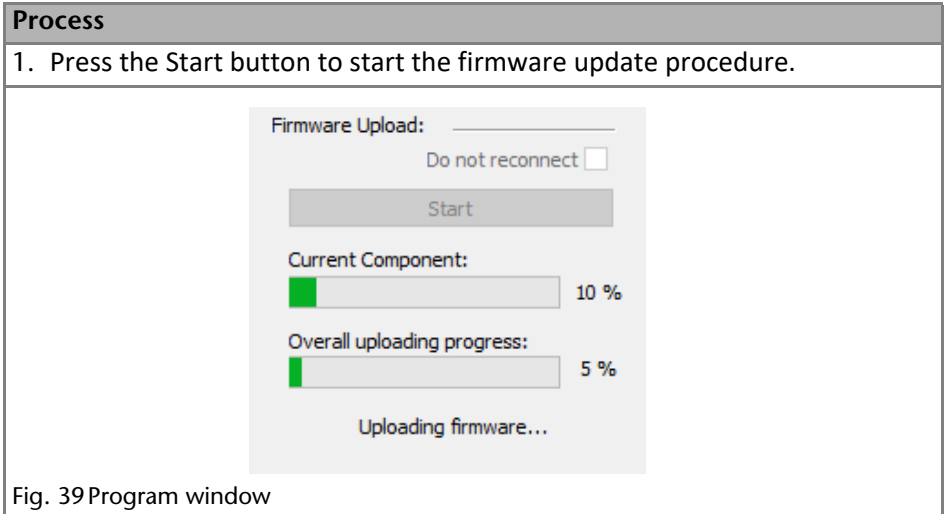


Fig. 39 Program window

When the upload is complete and successful, a green tick on the left side of the component line will be shown. The Firmware Wizard can be closed.


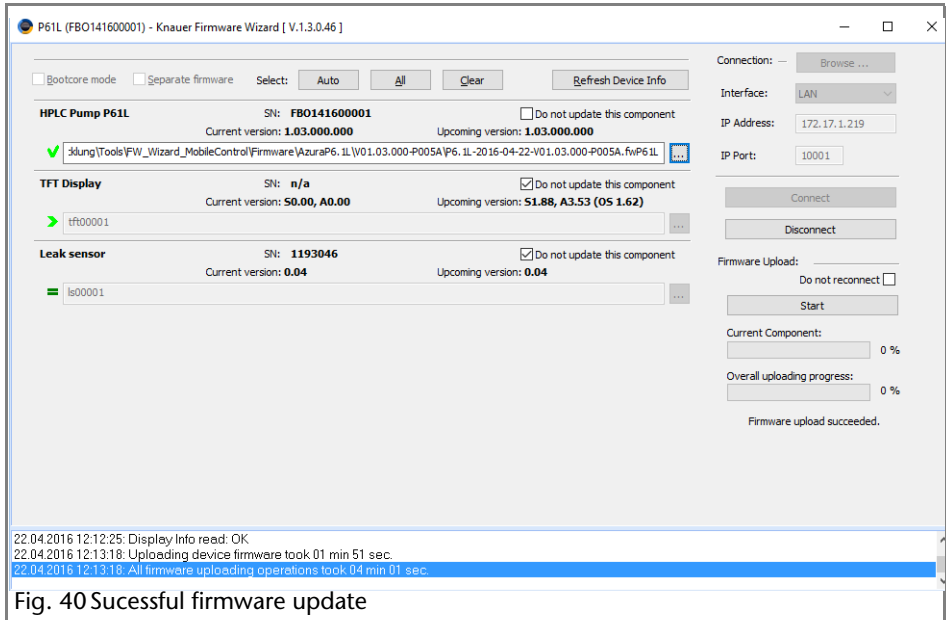


Fig. 40 Successful firmware update

Run Queue

The run queue is used to manage and schedule programs and sequences. Once a sequence or program is initiated, it is entered into the run queue automatically.

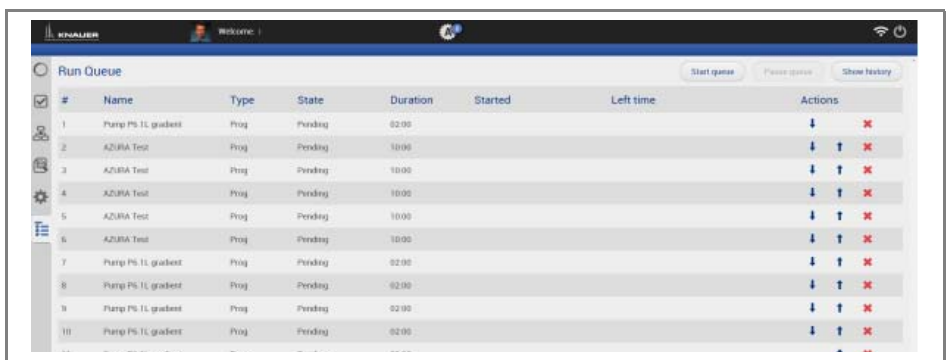

To view the current run queue, select the *RUN QUEUE* button. Each row in the run queue represents a program or sequence that is in process or waiting. From the run queue, you can view details about each run or sequence in the queue, including the following:

- Type - program run or sequence
- Name - name of resulting data file
- State - Pending, Running, Completed
- Duration
- Started
- Left time

If you are running a sequence, the Status shown in the sequence spreadsheet displays the status of an individual run in the sequence.

When there are items in the run queue, buttons at the top of the window will become available:

- *START* will start the run queue again after the *PAUSE* button has been pressed. This button is disabled when a sample is running or if the run queue is empty.
- *PAUSE* will pause the run queue as soon as one run is completed and wait for your action. You do not have to wait for the entire sequence to be completed.
- *ABORT* will immediately terminate the item currently running in the queue and pause the sequence.
- *SHOW HISTORY* allows to see all already processed programs or sequences.

Logs and Errors

Under *LOGS AND ERRORS*, all errors, system logs and communication logs are displayed.

SYSTEM LOGS are commands and signals which comprise the settings of an device. *COMMUNICATION LOGS* are control commands.

PERIOD: Changes the display period.

SYSTEM LOG, ERRORS, COMMUNICATION LOG: Activate the check box to display corresponding entries.

FILTER: Filters by user or source.

EXPORT TO TEXT, EXPORT TO XML: Saves all displayed messages to a text or XML file. The file is saved under the following path:

Note: In order to not overwrite the file, change the name of the existing file before you save the new file.

- Windows 8: Users/YourUserName/AppData/Local/Azura/tmp

Saving the file overwrites the existing file.

Energy Options

In the energy options, single devices or a system can be put into standby mode and woken up from standby mode.

Putting Devices into Standby Mode Manually

To put the device into standby mode manually, go to *DETAIL VIEW* and tap the button *STANDBY*.

Putting Devices into Standby Mode Automatically

To put the system or single devices into standby mode automatically, tap *SYSTEM SETUP · ENERGY OPTIONS · SYSTEM STANDBY*.

1. Tap the date and time buttons to enter the respective data.
2. To activate the standby mode, tick the check box.
3. Activate one of the options under *INSTRUMENT TO STANDBY*. *ALL* puts all devices which are part of the configuration into standby mode. Individual devices can be selected with *SELECT INSTRUMENT*. *NOT ONE* deactivates the standby mode for all devices.

Waking Up Devices from Standby Mode Manually

To wake up the device from standby mode, tap *DETAIL VIEW · POWER UP*. Note the waiting period which the lamp of the detector needs to be ready for use. Find the necessary data in the user manual of the device.

Waking Up Devices from Standby Mode Automatically

To put the system or single devices into standby mode automatically, tap *SYSTEM SETUP · ENERGY OPTIONS · SYSTEM WAKE-UP*.

1. Tap the date and time buttons to enter the respective data.
2. Under *SYSTEM WAKE-UP*, tick the check box.

Activate one of the options under *INSTRUMENT TO WAKE-UP*. Activating *ALL* wakes up all devices which are part of the configuration into standby mode. Individual devices can be selected with *SELECT INSTRUMENT*. *NOT ONE* deactivates the wake-up functionality for all devices.

User Management

The user account of the user that is logged-in is displayed under *USER MANAGEMENT*. The administrator can create new user accounts or assign rights to any user by activating the check boxes.



Fig. 41 User account settings *SYSTEM SETUP · USER MANAGEMENT*

CREATE NEW USER (for administrator only): Creates a new user.

CHANGE PASSWORD: Changes the password of the user. The user must have the respective permissions to do so (*SETTINGS > READ AND MODIFY*).

SETTINGS: *SETTINGS* refers to device and system settings. Select between *READ* and *READ AND MODIFY*.

GLP: Assigns reading rights for GLP data to the user.

PROGRAMS: The user may start programs/*RUN* or create programs, change or start them/*EDIT, ADD NEW AND ETC.*

LOGS: If checked, allows the reader to read the messages under *LOGS AND ERRORS*.

Creating User

If you execute the Mobile Control for the first time, you are prompted to enter a user name and a password. This user (administrator) has full access to the Mobile Control and can create additional users, edit them or assign rights.

1. Log in as administrator.
2. Select administrator under *SYSTEM SETUP*.
3. Tap *CREATE NEW USER*.
4. Enter a user name under *LOGIN*.
5. Tap *PASSWORD* and enter a password.
6. Tap *CONFIRM PASSWORD* and confirm password.

Changing the Password

Each user can change their own password, if they got the respective permissions from the administrator. The administrator can change the password of every user, without knowing their current password.

Changing Own Password

1. Select user under *SYSTEM SETUP*.
2. Tap *CHANGE PASSWORD*.
3. Tap *CURRENT PASSWORD* and enter the current password.
4. Tap *NEW PASSWORD* and enter the new password.

5. Tap *CONFIRM NEW PASSWORD* and confirm the new password.
6. Confirm with *OK* and set with *APPLY*.

Changing the Password of Another User

1. Log in as administrator.
2. Under *SYSTEM SETUP*, select the user whose password shall be changed.
3. Tap *CHANGE PASSWORD*.
4. Tap *NEW PASSWORD* and enter the new password.
5. Tap *CONFIRM NEW PASSWORD* and confirm the new password.
6. Confirm with *OK* and set with *APPLY*.

GLP Data

Under GLP, you find a list of all devices for which GLP data can be displayed. Tap the respective device to display GLP data.

General

All Devices	Serial number
	Firmware version
	Operation time
	Installation date
	Last service
Leak sensor	Serial number
	Firmware version

Detector

Optical Properties	Optical bandwidth at 656 nm FWHM	
	Stray light (AU)	
	Light intensity I-Sig at UV-maximum (nW)	Light intensity of the signal channel at UV-maximum (nW)
	Light intensity I-Ref at UV-maximum (nW)	Light intensity of the reference channel at UV-maximum (nW)
Lamps - Power Supply	Number of filter wheel switches	
	Serial number	
	Operation time	
	Firmware version	

	Supply number	
Lamps - Deuterium Lamp	Deuterium serial number	
	Operation time	
	Starts	Displays how often the lamp has been switched on.
Light Source (RID 2.1L)	Operating time-Installation date-Light source number	
Optical Cell (RID 2.1L)	Cell batch number	
Validation Data (RID 2.1L)	Last measured span	
	Last measured span date	
Pump		
Head left/right	Serial number	
	Operation time	
	Cycles	Piston cycles
	Volume	Total volume conveyed to date
	P-index	Unit for the current strain of the pump
	Head type	Pump head type
Pump	Operation time	
Assistant		
Left, Middle, Right device	Serial number	
Pump	Firmware version	
	Operation time	
	Head type	Pump head type
Valves	Configuration	
	Switching cycles	Displays how often the valve has been switched.
Detector	Firmware version	

Operation time

Starts

Displays how often the lamp has been switched on.

The view of the assistant depends on the installed devices.

Index

Symbols

- (Disconnected) 16
- + (Connected) 16
- + Add 31

A

- Activate 7, 8
- Activation code 7, 8
- Add new device 16
- Admin
 - administrator settings 51
- Administrator account 8
- Administrator settings
 - administrator 51
- Apply 11, 25
- Assistant
 - device settings 19
 - Device status 30
 - GLP data 54
 - serial number 54
- Autosampler 20, 24, 30
- Autozero 26

B

- Browse 16
- Buttons 11

C

- Cancel 11, 25
- Certificate 6
- Change password 52
- Chromatography software 14, 17
- Column thermostat 24, 30
 - settings 29
- Communication Logs 50
- Configuration 20
- Connected 16
- Create new user 52

D

- D2 lamp 19
- Delay (Fraction Collection)
 - device settings 18
- Demo version 7
- Detail View 13, 14, 24, 25
 - Pump 25
- Detector
 - device settings 18
 - Device status 29
 - GLP data 53
 - initiate autozero 26
 - insert wavelength 26
 - Optical bandwidth 53
 - Stray light 53
 - switch on lamp 26
 - System Overview 23
 - System Setup 18
 - wavelength 29
- Deuterium lamp
 - Serial number 54
- Device
 - Add new device 16
 - change selection 16
 - firmware version 53
 - handheld 7
 - Operation time 53
 - Run time 53
 - select 16
 - Serial number 53
 - serial number 17, 29
 - standby 51
- Device code 7, 8
- Device settings
 - Assistant 19
 - Detector 18
 - DHCP 17
 - I-Sig Abs 18
 - Instrument Lookup 17

- Integration 18
- IP Address 17
- Leak Sensor 18
- Mixing chamber 19
 - name 17
 - offset 18
- Port (Fraction Collection) 18
- Pump 19
- Pump head 19
- S/N 17
- Scale 18
- Sensitivity 18
- serial number 17
- Static 17
- System Setup 17
- Device Status 24, 25, 29
 - Assistant 30
 - Detector 29
 - Pump 29
 - valve 30
- DHCP 17
 - device settings 17
- Direct signal 26
- Disconnected 9, 16
- E**
- Emergency mode 15
- Energy Options 12, 51
- Errors 50
 - messages 12, 13, 14
- Events 25
 - set 25
- Export to text 50
- Export to XML 50
- F**
- Filter 50
- Firmware version
 - device 53
 - Lamp and Power Supply 53
 - Leak sensor 53
 - Flow rate settings 27
 - Full version 7
 - activate 7
- G**
- Gateway 14, 18
 - device settings 18
 - System Setup 18
- GLP 12, 14, 52
- GLP data 53
 - Assistant 54
 - Detector 53
 - firmware version 53
 - general 53
 - installation date 53
 - Operation time 53
 - Pump 54
 - Pump head type 54
 - Run time 53
- Gradient Type
 - Isocratic 19
- Gradients
 - composition 27
 - high-pressure gradient (HPG) 27
 - low-pressure gradients (LPG) 27
- H**
- Handheld devices 7
- High Sensitivity 18
- I**
- IMax sensitivity 20
- IMin sensitivity 20
- Initiate an autozero 26
- Instrument Lookup 12, 16
 - Browse 16
 - device settings 17
- Instrument lookup 9
 - Broadcast Port 14
- Instruments to standby 51
- Instruments to wake-up 51

- Integration
 - device settings 18
- Inverted signal 26
- IP Address 14, 17
 - allocate automatically 17
 - allocate manually 17
 - device settings 17
 - System Setup 17
- Isobar Mode 28
- Isocratic 19
- L**
- Lamp
 - switch on 26
- Lamp and Deuterium Lamp
 - Operation time 54
 - Run time 54
- Lamp and Power Supply
 - Firmware version 53
 - Operation time 53
 - run time 53
 - Serial number 53
 - Supply number 54
- LAN 6, 9, 14, 17
- Leak Sensor 18
 - adjust sensitivity 18
 - device settings 18
 - set 18
 - switch on 18
- Leak sensor
 - Firmware version 53
 - Serial number 53
 - switch off 18
- License 7
 - activate 7
 - demo version 7
 - full version 7
 - validity 9
- License validity 9
- Log lifetime 15
- Logout 13
- Logs 52
- Logs and Errors 12, 14, 50
- Low Sensitivity 18
- M**
- MAC address 7
- Medium
 - Sensitivity 18
- Mixing chamber
 - device settings 19
 - set 19
- Mobile Control app
 - exit 10
 - install 7
 - install on handheld device 7
 - integrate into a network 9
 - operate 11
 - start 9
 - update 10
- Monitor Mode 10
- N**
- Name 17
 - device settings 17
 - system setup 17
- Network adapter 7
- Network Settings 12, 14
- O**
- Offset 18
 - device settings 18
 - System Setup 18
- Operation 11
- Operation time
 - Lamp and Deuterium Lamp 54
 - Lamp and Power Supply 53
 - Pump 54
 - Pump head 54
- P**
- Password 9, 52
 - change 52

- Period 50
- PMax Mode 28
- Preferences 12, 15
- Prerequisites 11
- Pressure settings 27
- Pressure Units 15
- Program
 - see Programs
- Programs 12, 14, 30, 52
 - + Add 31
 - add device 31
 - Copy 31
 - create 31
 - Delete 31
 - Edit 31
 - Load 31
 - Save 31
 - Start 32
 - Stop 32
- Pump
 - Detail View 25
 - device settings 19
 - Device status 29
 - GLP data 54
 - gradients 27
 - Isobar Mode 28
 - Operation time 54
 - PMax Mode 28
 - Pump head type 54
 - purge 28
 - run time 54
 - Serial number 54
 - set flow rate 27
 - set pressure 27
 - start 28
 - System Overview 23
- Pump head
 - device settings 19
 - Operation time 54
 - run time 54
- Pump head type 54
- Purge 28
- R**
- Router 16
- Run time
 - Lamp and Deuterium Lamp 54
 - Lamp and Power Supply 53
 - Pump 54
 - Pump head 54
- S**
- S/N
 - serial number 29, 53
- Scale
 - device settings 18
 - System Setup 18
- Sensitivity 18
 - device settings 18
 - High 18
 - Low 18
 - Medium 18
 - System Setup 18
- Serial number
 - Assistant 54
 - Deuterium lamp 54
 - device 17, 29, 53
 - general 29, 53
 - Lamp and Power Supply 53
 - leak sensor 53
 - Pump 54
 - S/N 29, 53
- Set Date 15
- Set Time 15
- Setting temperature safety limits 29
- Setting the target temperature 29
- Settings 52
- Short-Circuit Bridge
 - see Events
- Signal Mode
 - signal settings 26

- Signal source 19
 - Software 6, 14, 17
 - Standby 25, 51
 - Start 9
 - Start Input 19
 - Static 17
 - Device settings 17
 - Status bar 14
 - Stop 13
 - Subnet Mask 14, 17
 - allocate manually 17
 - device settings 17
 - Supply number
 - Lamp and Power Supply 54
 - System detector units 15
 - System Logs 15, 50
 - System Overview 12, 13, 14, 23, 25
 - Detail View 24
 - System Setup 9, 12, 14
 - Detector 18
 - energy options 12
 - gateway 18
 - General 17
 - Instrument Lookup 12
 - IP Address 17
 - Network Settings 12
 - offset 18
 - Preferences 12
 - Scale 18
 - Sensitivity 18
 - Static 17
 - Subnet Mask 17
 - User management 12
 - System setup DHCP 17
 - Leak Sensor 18
 - name 17
 - S/N 17
 - System standby 51
 - System wake-up 51
- T**
- Time constant 18, 19
- U**
- Update 10
 - User
 - change 10
 - create 52
 - log-out 10, 14
 - User account 6
 - create 52
 - User management 12, 51
 - User name 8, 9, 52
 - User settings 51
- V**
- Valve
 - change position 29
 - choose 28
 - Device status 30
 - System Overview 23
 - virtual keyboard 15
- W**
- Wait for temperature 15
 - wake-up 51
 - Wavelength 29
 - WLAN 9
 - router 9

© KNAUER Wissenschaftliche Geräte GmbH
All rights reserved.
Technical data are subject to change without notice.
Please check our website for latest updates and changes.
Version 3.1
Last manual update: 2016-08-22
Printed in Germany on environmentally friendly paper
from sustainable forests.

® AZURA are registered trademarks of:
KNAUER Wissenschaftliche Geräte GmbH

▶ See up-to-date manuals online:
www.knauer.net/downloads

www.knauer.net

HPLC · SMB · Osmometry

KNAUER Wissenschaftliche Geräte GmbH
Hegauer Weg 38
14163 Berlin, Germany

Phone: +49 30 809727-0
Telefax: +49 30 8015010
E-Mail: info@knauer.net
Internet: www.knauer.net



© KNAUER 2015 V6851/0.0/05.14/---